



alpha laboratories Fastrak Tip Refill System Instructions

[Home](#) » [alpha laboratories](#) » alpha laboratories Fastrak Tip Refill System Instructions 

Contents

- 1 alpha laboratories Fastrak Tip Refill System
- 2 Product Information
- 3 Product Usage Instructions
- 4 INSTRUCTIONS FOR USE
- 5 Documents / Resources
 - 5.1 References

A L P H A

alpha laboratories Fastrak Tip Refill System



Product Information

The product is a rack system designed for efficient storage and dispensing of tips. It is manufactured by Alpha Labs and the model number is JBN370A.

Product Usage Instructions


1. Remove the plastic wrapping from the box.
2. Before opening, tap the base of the box on the bench to ensure all tips are fully nested.
3. Undo the paper tab and open the box by lowering both sides simultaneously.
4. Grasp the transfer top and lift the tips straight up and clear of the stack.
5. Align the transfer top over the rack and place it onto the rack.
6. Push down on the transfer top to release the refill plate and click it onto the rack.
7. Once in use, do not invert the pack.
8. Store the boxes upright to ensure good tip nesting.

INSTRUCTIONS FOR USE


1. Remove plastic wrapping. Before opening, tap the base of the box on the bench to ensure all tips are fully nested.



2. Undo the paper tab and open the box by lowering both sides simultaneously. Grasp transfer top. Lift tips straight up and clear of stack.

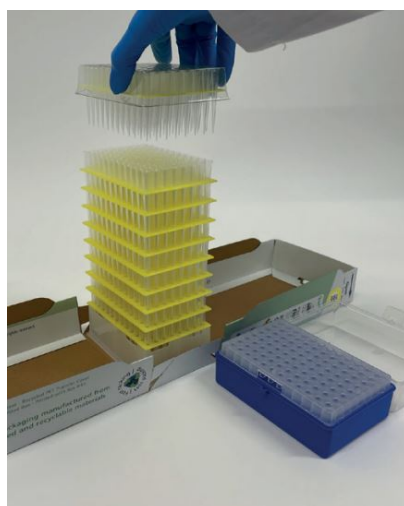

3. Align transfer top over rack and place onto rack.

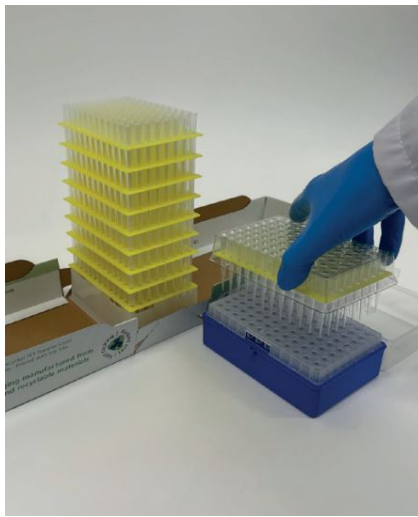

4. Push down on transfer top to release refill plate and click onto rack. Return transfer top to Fastrak® refill stack.



5. Close the Fastrak® carton, lifting both sides simultaneously and re-seal the paper tab. Once in use, do not invert the pack.



6. Store boxes upright to ensure good tip nesting.



NOTE: For more information and to watch a demonstration video, please visit www.alphalabs.co.uk/fastrak.

JBN370A 01.23

Alpha Laboratories Ltd 40 Parham Drive Hampshire SO50 4NU UK

- **Tel:** +44 23 8048 3000
- **Email:** sales@alphalabs.co.uk
- **Web:** www.alphalabs.co.uk

Documents / Resources

	<p>alpha laboratories Fastrak Tip Refill System [pdf] Instructions Fastrak Tip Refill System, Fastrak, Tip Refill System, Refill System</p>
--	---

References

- [Fastrak® Pipette Tip Refill System | Eco-friendly Pipette Tips](#)

Manuals+