


AVEN
26700-220-479
Mighty Scope
Clear Vue



AVEN 26700-220-479 Mighty Scope Clear Vue Owner's Manual

[Home](#) » [AVEN](#) » AVEN 26700-220-479 Mighty Scope Clear Vue Owner's Manual 

Contents

- [1 AVEN 26700-220-479 Mighty Scope Clear Vue](#)
- [2 Product Specifications](#)
- [3 Product Usage Instructions:](#)
- [4 Frequently Asked Questions \(FAQ\)](#)
- [5 INTRODUCTION](#)
- [6 Features](#)
- [7 SPECIFICATION](#)
- [8 SETTING INSTRUCTION](#)
- [9 Documents / Resources](#)
 - [9.1 References](#)
- [10 Related Posts](#)



AVEN 26700-220-479 Mighty Scope Clear Vue



Product Specifications

- **Item #:** 26700-220-479
- **Magnification:** 8x-25x | 40x-120x (on 24" Monitor)
- **Effective Pixels:** 1920(H) x 1080(V)
- **Image Sensor:** 2M CMOS Sensor
- **Working Distance:** 45mm-95mm
- **Interface:** Mini USB – Type A / HDMI
- **PC Software:** EZImage X3
- **Illumination:** Integrated Ring Light

Product Usage Instructions:

Setup:

1. Connect the Mighty Scope ClearVue digital microscope to a power supply using the USB cable.
2. Adjust the magnification level using the focus wheel to achieve the desired zoom level.

Illumination Adjustment:

Use the adjustable LEDs to control the lighting for optimal image clarity.

Image Capture:

1. Press the snap shot remote control to capture images.
2. Save captured images directly to the included 32GB MicroSD card for easy storage.

Viewing and Sharing:

1. View captured images on the crystal-clear 5" screen.
2. Process and share images using the ezImage X3 software included with the product.

Frequently Asked Questions (FAQ)

- **Can I adjust the magnification level of the Mighty Scope ClearVue?**

Yes, you can adjust the magnification from 8x to 25x or up to 120x when connected to a 24" monitor.

- **What is included in the package contents?**

The package includes the Mighty Scope ClearVue digital microscope, USB cable, power supply, snap shot remote control, 32GB MicroSD card, 2600mAh 18650 Li-ion battery, user guide, ClearVue post stand, gliding stage, and free ezImageX3 software download.

INTRODUCTION

Mighty Scope ClearVue has a built in 5" high resolution LED screen and offers a magnification range of 8x -25x. The included rechargeable battery allows you to operate the microscope without a power supply for hours thus making it suitable for field projects. You can also connect this microscope to an external monitor via an HDMI output for viewing images on a larger screen. Images can be saved on the included Micro SD card. Can also be mounted onto a variety of stands to suit your application. Includes the Mighty Scope ClearVue Gliding Stage, with horizontal and vertical movement, as well as a rotating base. Ideal for smooth scanning of objects.

Features

- Compact All-in-one system
- 1080p Full HD Imaging at 60 FPS
- HDMI Output for additional monitor viewing
- ezImageX3 software download included
- 5" built-in HD screen
- Save high resolution image/video files directly to microSD card
- Enhances productivity – eliminates eye strain and body fatigue
- 1080p Full HD Output



Crystal Clear 5" Screen



EzImage X3 Download Included



Aven's Dual Gooseneck LED Task Light



Recommended for
improved imaging
& color control

Sold Separately **#26537**


SPECIFICATION

Item #	26700-220-479
Magnification	8x-25x 40x-120x (on 24" Monitor)
Effective Pixels	1920(H) x 1080(V)
Image Sensor	2M CMOS Sensor
Working Dist.	45mm-95mm
Interface	Mini USB – Type A / HDMI
PC Software	EZImage X3
Illumination	Integrated Ring Light
Package Contents	Mighty Scope ClearVue digital microscope, USB cable, Power supply, Snap shot remote control, 32gb MicroSD Card, 2600mAh 18650 Li-ion battery, User guide, ClearVue Post Stand, Gliding Stage, Free ezImageX3 software download

SETTING INSTRUCTION

	<p>Mighty Scope ClearVue on Spring Balanced Arm w/ Table Clamp #26700-220-559</p> <p>This arm can be raised, lowered, and tilted as needed. The swivel motion makes it easy to move the ClearVue out of the way when not in use.</p> <ul style="list-style-type: none"> • Fully Covered 31" Spring Balanced Arm • Adjustable Tension Knob
	<p>Mighty Scope ClearVue on Flex Arm Stand w/ Table Clamp #26700-220-557</p> <p>This arm easily bends and stays put in any position. Tilt, rotate, and position as needed</p> <ul style="list-style-type: none"> • 18" Steady Mobile Flex Arm • 360° Swivel Head • 2" Table Clamp
	<p>Mighty Scope ClearVue on Flex Arm Stand w/ Magnetic Base #26700-220-558</p> <p>This arm easily bends and stays put in any position. Tilt, rotate, and position as needed</p> <ul style="list-style-type: none"> • 18" Steady Mobile Flex Arm • 360° Swivel Head • 3" Magnetic Base
	<p>Mighty Scope ClearVue on Post Stand #26700-220-MNT</p> <ul style="list-style-type: none"> • 7"x8.5" Metal Base • Vertical adjustment along 10" Post
	<p>Compact Gliding Stage XYR #26800B-479</p> <ul style="list-style-type: none"> • Smooth vertical and horizontal movement w/ Built-in rotating stage • Compact design (5.3 x 4.7 x 3.6 inch base) • Metal rack and pinion construction • Designed to be used with optical and digital microscope systems
	<p>Dual Gooseneck LED Task Light With Magnetic Base And 4" Table Clamp #26537</p> <ul style="list-style-type: none"> • Smooth vertical and horizontal movement w/ Built-in rotating stage • Powerful, long lasting ultra-bright LEDs provide 500lm of flicker and ghost-free illumination • On/Off button toggles between the lamps for single-lamp or dual-lamp lighting • Can be powered 3 ways: with included power supply, 3 AA batteries, or with a rechargeable lithium ion battery

Documents / Resources

	<p>AVEN 26700-220-479 Mighty Scope Clear Vue [pdf] Owner's Manual 26700-220-479, 26700-220-559, 26700-220-557, 26700-220-558, 26700-220-MNT, 26700-220-479 Mighty Scope Clear Vue, 26700-220-479, Mighty Scope Clear Vue, Scope Clear Vue, Clear Vue</p>
-------------------------------------------------------------------------------------	------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------

References

- [User Manual](#)

[Manuals+](#), [Privacy Policy](#)

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.