



azure biosystems AzureSpot Western Blot Image Analysis Software Installation Guide

[Home](#) » [azure biosystems](#) » azure biosystems AzureSpot Western Blot Image Analysis Software Installation Guide 



AzureSpot Western Blot Image Analysis Software Installation Guide

Contents

- [1 Installing AzureSpot Pro on Windows](#)
- [2 Installing AzureSpot Pro on macOS](#)
- [3 Licensing](#)
- [4 Delegated Licensing](#)
- [5 Support](#)
- [6 Documents / Resources](#)
 - [6.1 References](#)
- [7 Related Posts](#)

Installing AzureSpot Pro on Windows

System Requirements

Item	Required	Recommended
Operating System	Windows 8.1 64bit	Windows 10 Professional 64bit
CPU	Intel Core i3	Intel Core i5 / AMD Ryzen 5
GPU	WebGL 1.0 Compatible	WebGL 2.0 Compatible
RAM	2GB	16GB
Storage	2GB	4GB

Installation Process

1. Download and install Microsoft .NET 5 Desktop Runtime: <https://dotnet.microsoft.com/download/dotnet/5.0>
2. Install Microsoft Visual C+ + x64 re-distributable: <https://www.microsoft.com/en-us/download/confirmation.aspx?id=52685&6B49FDFB-8E5B-4B07-BC31-15695CSA2143=1>
3. Run the MSI installer included in this distribution
4. Launch the app from the desktop shortcut or the icon in the start menu

Installing AzureSpot Pro on macOS

Installing AzureSpot Pro on macOS

Item	Required	Recommended
Operating System	macOS 10.15 Catalina	macOS 11 Big Sur
Architecture	Intel	Intel

Any mac supported by 10.15 Catalina is supported by Azure Spot Pro.

Installation Process


1. Download and install Microsoft .NET 5 Runtime: <https://dotnet.microsoft.com/download/dotnet/5.0>
2. Open the DMG included in this distribution
3. Drag the AzureSpot Pro icon into your Applications folder
4. Launch the app from the Applications folder or LaunchPad

Licensing

Applies to both macOS and Windows.

To run the app, you must provide your activation key and allow the system to perform a one-time activation from our license server.

- Launch AzureSpot Pro
- From the Home screen, click ConfigureLicense (pictured below)
- Enter your activation key into the AcquireLicense field and click Acquire
- Close the configure license window

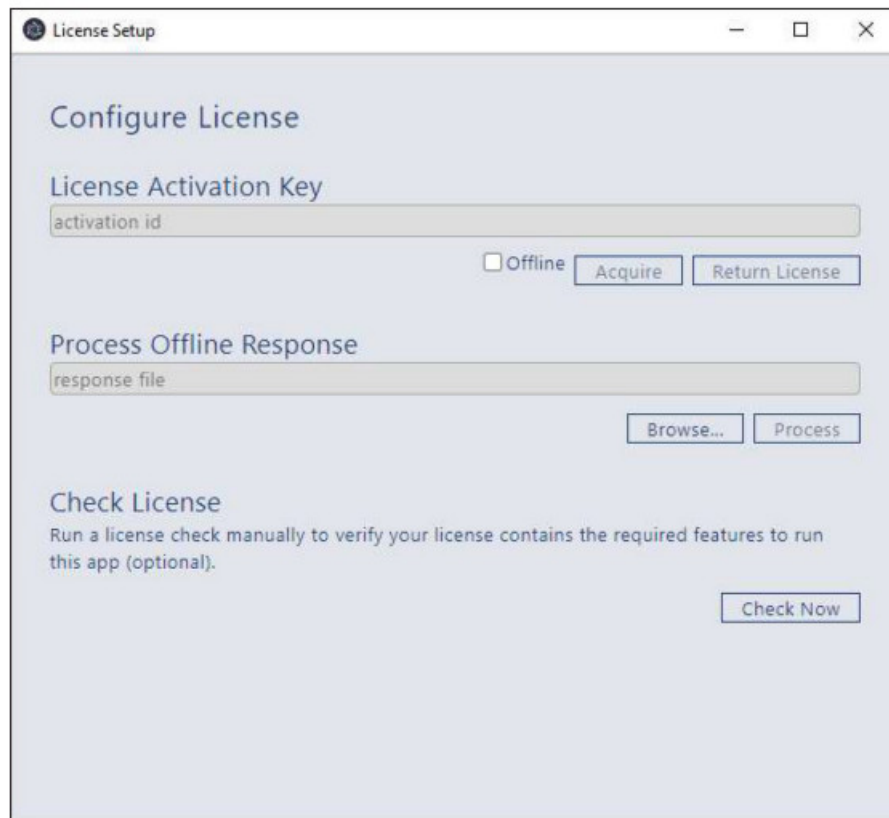

Delegated Licensing

Applies to both macOS and Windows.

License activation requires an active internet connection. If this is not available we can use a fall back where another machine that does have internet access can process the activation request as a proxy. Let's call the target, offline machine, PC1 and the support machine that has internet access PC2.

Start on PC1

- Launch AzureSpot Pro
- From the Home screen, click ConfigureLicense
- Enter your activation key into the activation id field
- Check the Delegate box
- Click Acquire
- Save the request file to your desktop

Move to PC2

- Install AzureSpot Pro
Note: AzureSpot Pro can be uninstalled from the 2nd machine after this process is complete and it will not require, or consume a license
- Launch AzureSpot Pro and open the Configure

License window

- Click Browse in the Delegated Activation section
- Select the activation request file from PC1

- Click Process
- Notice a new file, the response file, is generated alongside the request file
- Send the response file to PC7
- Uninstall AzureSpot Pro (if necessary) Back on PC1
- Launch AzureSpot Pro and open the Configure License window
- Click Browse in the Delegated Activation section
- Locate and open the processed response file
- Click Integrate
- Close the configure license window

Support

If you have any trouble, or need any further help, please contact us: support@azurebiosystems.com



www.azurebiosystems.com
info@azurebiosystems.com

Copyright © 2021 Azure Biosystems. All rights reserved. The Azure Biosystems logo, and Azure Biosystems® are trademarks of the
 Company. All other trademarks, service marks and trade names appearing in this brochure are the property of their respective owners.

Documents / Resources

	<p>azure biosystems AzureSpot Western Blot Image Analysis Software [pdf] Installation Guide</p> <p>AzureSpot, Western Blot Image Analysis Software, AzureSpot Western Blot Image Analysis Software, Blot Image Analysis Software, Image Analysis Software, Analysis Software, Software</p>
--	--

References

- www.azurebiosystems.com
- [Download .NET 5.0 \(Linux, macOS, and Windows\)](#)