



[Home](#) » [azure biosystems](#) » **azure biosystems Western Blot Image Analysis Software Installation Guide** 

Contents [[hide](#)]

- [1 azure biosystems Western Blot Image Analysis Software](#)
- [2 Installing on Windows OS](#)
- [3 Installing on Mac OS](#)
- [4 Licensing](#)
- [5 Delegated Licensing](#)
- [6 Support](#)
- [7 FAQ](#)
- [8 Documents / Resources](#)
 - [8.1 References](#)



azure biosystems Western Blot Image Analysis Software



Installing on Windows OS

System Requirements:

Item	Required	Recommended
OS	Windows 10 Professional 64-bit	Windows 11 Professional 64-bit
CPU	Intel Core i3	Intel Core i5 / AMD Ryzen 5
GPU	WebGL 1.0 Compatible	WebGL 2.0 Compatible
RAM	2 GB	16 GB
Storage	2 GB	4 GB

Installation Process:

1. Download and install Microsoft .NET 8 Desktop Runtime 8.x.x:
<https://dotnet.microsoft.com/en-us/download/dotnet/8.0>
 Important: From the .NET Desktop Runtime 8.x.x section choose specifically the x64 variant
2. Download and Install Microsoft Visual C++ x64 re-distributable:
https://aka.ms/vs/17/release/vc_redist.x64.exe
3. Run the installer included in this distribution

4. Launch the app from the desktop shortcut or the icon in the start menu

Installing on Mac OS

System Requirements:

Item	Required	Recommended
OS	macOS 10.15 Catalina	Windows 11 Professional 64-bit

Note: Any Mac supported by macOS 10.15 Catalina is compatible with AzureSpot Pro.

Installation Process:

1. Download and install Microsoft .NET 8 Runtime 8.x.x:

<https://dotnet.microsoft.com/en-us/download/dotnet/8.0>

Important: From the .NET Runtime 8.x.x section under the Installers column, choose the Arm64 (if you have an M1-M3 processor in Mac) or choose x64 (if you had Intel processor)

2. Drag AzureSpot Pro icon into your Applications folder
3. Launch the app from the Applications folder or LaunchPad

Licensing

This following instruction applies to both macOS and Windows. In the following guides, when we'll instruct to "launch the Configure License tool", please follow these two steps:

- Launch AzureSpot Pro
- From the Home screen, click on Configure License

To run the app, you must provide your activation key and allow the system to perform a one-time activation from our license server:

- Launch the Configure License tool
- Enter your activation key into the Acquire License field and click Acquire
- Close the Configure License window

Delegated Licensing

License activation requires an active internet connection. If this was not available on the target PC, then you may use a fallback where another machine that does have active internet access can process the activation request as a proxy. Let's call the target offline computer as PC1 and the support online computer PC2.

START ON PC1:

- Launch the Configure License tool
- Enter your activation key into the Activation ID field
- Check the Delegate box
- Click Acquire
- Save the request file to your desktop
- Move the request file somehow to PC2 (e.g. USB drive)

MOVE TO PC2:

- Install AzureSpot Pro (note: AzureSpot Pro can be uninstalled from the 2nd machine after this process is complete and it will not require nor consume a license)
- Launch the Configure License tool
- Click Browse in the Delegated Activation section
- Select the activation request file from PC1
- Click Process
- Notice a new file, the response file, is generated alongside the request file
- Move the response file somehow back to PC1 (e.g. USB drive)
- You may uninstall the AzureSpot Pro from PC2 (if necessary)

BACK ON PC1:

- Launch the Configure License tool
- Click Browse in the Delegated Activation section
- Locate and open the processed response file
- Click Integrate
- Close the Configure License window

Support

If you needed any further assistance, please contact our support team at:

support@azurebiosystems.com

Headquarters: 6747 Sierra Court, Suites A-B • Dublin, CA 94568 • Tel +1.925.307.7127 •

Fax +1.925.905.1816

4810 Jean Talon 0., Suite 328 • Montreal, QC, Canada H4P 2N5 • Tel +1.4382902000


info@azurebiosystems.com • vomu.azurebiosystems.com

FAQ

- **Q: What should I do if I encounter an error during installation?**

A: If you encounter errors during installation, make sure to check that your system meets the minimum requirements specified in the manual. Additionally, ensure you have a stable internet connection for licensing activation.

Documents / Resources

	azure biosystems Western Blot Image Analysis Software [pdf] Installation Guide Western Blot Image Analysis Software, Image Analysis Software, Analysis Software, Software
---	--

References

- [User Manual](#)

■ azure biosystems

◆ Analysis Software, azure biosystems, Image Analysis Software, Software, Western Blot Image Analysis Software

Leave a comment

Your email address will not be published. Required fields are marked *

Comment *

Name

Email

Website

☐ Save my name, email, and website in this browser for the next time I comment.

Post Comment

Search:

e.g. whirlpool wrf535swhz

Search

[Manuals+](#) | [Upload](#) | [Deep Search](#) | [Privacy Policy](#) | [@manuals.plus](#) | [YouTube](#)

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.