Home » BROOKFIELD AMETEK » BROOKFIELD AMETEK AMETEK DVPlus Viscometer User Guide



BROOKFIELD AMETEK AMETEK DVPlus Viscometer User Guide

Contents

1 BROOKFIELD AMETEK AMETEK DVPlus

Viscometer

- **2 Product Information**
- **3 Product Usage Instructions**
- 4 Steps to update firmware
- 5 FAQ
- 6 Documents / Resources
 - **6.1 References**

BROOKFIELD AMETEK AMETEK DVPlus Viscometer

Product Information

Specifications:

• Model: DVPlus

• Manufacturer: Brookfield Engineering • Contact: 800-628-8139 or 508-946-6200 • Website: brookfieldengineering.com

Product Usage Instructions

Steps to Update Firmware:

- 1. Insert a USB Flash Drive (FAT32 Format) with the Firmware Update loaded in the root folder into the USB A Port on the back ofthe DVPlus.
- 2. Do not remove the USB Flash Drive during the upgrade process.
- 3. Turn on the DVPlus.
- 4. Level and AutoZero the device accordingly.
- 5. Go to the Settings Menu and select Firmware Update.
- 6. Select the desired Firmware version from the USB Drive.
- 7. Confirm your choice when prompted.
- 8. Press Update to start the firmware update process. The DVPlus will automatically update the firmware to the selected version.
- 9. Wait for the update to complete. The DVPlus will automatically reboot. Do not remove the USB Flash Drive during this time.
- 10. Verify the installed firmware version by viewing the startup Splash Screen.
- 11. Once the DVPlus finishes booting, you can safely remove the USB Flash Drive.

Steps to update firmware

WARNING: It is critical to only install Firmware intended for the DVPlus on your instrument. Please ensure that no other .fwu files exist on your USB Flash Drive and that the only firmware files that exist are those you are trying to install.

- 1. Insert a USB Flash Drive (FAT32 Format) with the Firmware Update loaded in the root folder into the USB A Port on the back of the DVPlus. DO NOT Remove the USB Flash Drive at any point during the upgrade.
- 2. Turn the DVPlus On.
- 3. Level and AutoZero accordingly.
- 4. Go to Settings Menu and select Firmware Update.
- 5. Select the desired Firmware version from the USB Drive
- 6. DVPlus will have you confirm your choice
- 7. Press Update and the DVPlus will automatically update the firmware to the desired selection. When the update is complete, the DVPlus will automatically reboot. DO NOT Remove the USB Flash Drive at any point during the update.
- 8. Verify that the desired Firmware version has been installed by viewing the startup Splash Screen.
- 9. After the DVPlus finishes booting, the USB Flash Drive can be removed.
- 800-628-8139 or 508-946-6200
- brookfieldengineering.com

FAQ

Q: Can I use any USB Flash Drive for firmware updates?

A: It is recommended to use a USB Flash Drive formatted in FAT32 for firmware updates to ensure compatibility.

Q: What should I do if I encounter an error during the firmware update process?

A: If you encounter an error, contact customer support at 800-628-8139 or 508-946-6200 for assistance.

Documents / Resources



BROOKFIELD AMETEK AMETEK DVPlus Viscometer [pdf] User Guide AMETEK DVPlus Viscometer, DVPlus Viscometer

References

User Manual

Manuals+, Privacy Policy

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.