



CONTROL SOLUTIONS VFC400 Vaccine Temperature Data Logger User Guide

[Home](#) » [CONTROL SOLUTIONS](#) » CONTROL SOLUTIONS VFC400 Vaccine Temperature Data Logger User Guide 



VFC400 Download of Data Downloading data from your the VFC400

Contents

- [1 Downloading data](#)
- [2 Overview](#)
- [3 Documents / Resources](#)
 - [3.1 References](#)
- [4 Related Posts](#)

Downloading data

Buttons Actions

Overview



Review/Mark button

Press to enter REVIEW mode and then press again to advance through the min/max temperatures. This action also places an inspection mark in the log if RECORDING is active. This validates your twice daily temperature logs.

Start/Clear/Stop button

Press to start recording
 Press to Stop recording
 Press to exit day summary


Downloading data

- The first thing we want to do before we download data from the VFC400 is stop the data logger from logging temperatures.



Downloading data


Stopping the data logger



When it is time to download the data from the data logger we suggest stopping the logger. To accomplish this you hold down the Start/Clear/Stop button for 1-2 seconds. The word STOPPED will appear on the screen as shown on the top left picture. The REC symbol will disappear after 1-2 seconds and you take your finger off the button quickly.

The logger will stop and the screen will only say STOPPED and show the number of days collected.

Downloading data

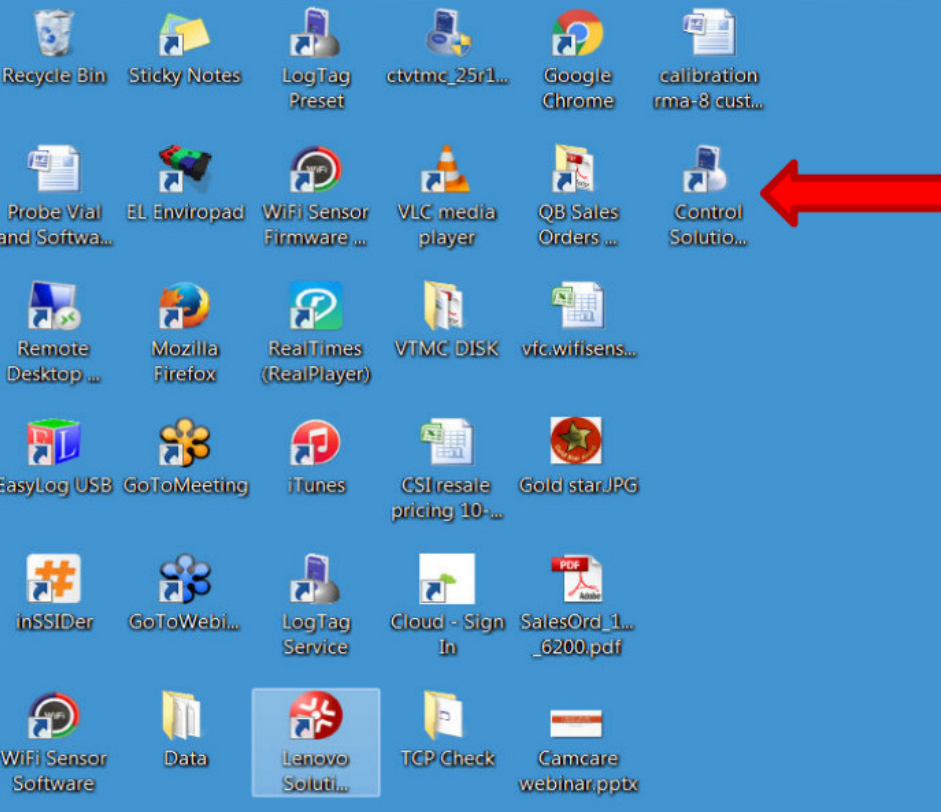

At your PC open up the Control Solutions VTMC software by double clicking on the Icon.

VFC400 Data Logger Configuration

- Plug your Docking Station into the USB port of your PC
- Insert data logger into the Docking Station firmly




Note: Make sure to download your data from your data logger prior to configuring/re-configuring as this will erase any and all data stored on your data logger.


Docking Station

Downloading data

- After you insert the logger into the docking station you will see the following message:

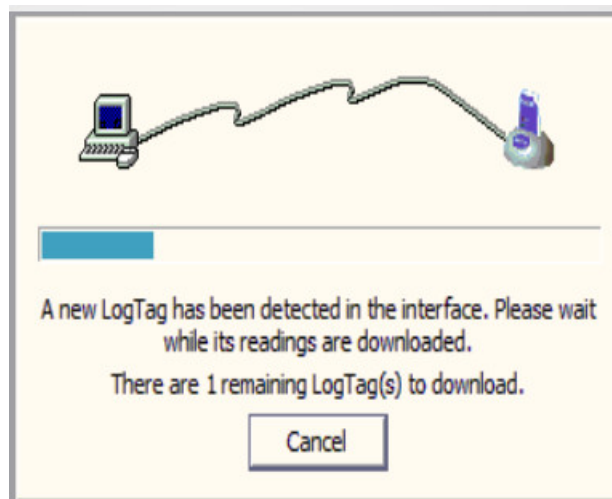


A few seconds later the chart/data will pop-up on the screen and then the following message: Click OK in the box and remove the data logger from the docking station.

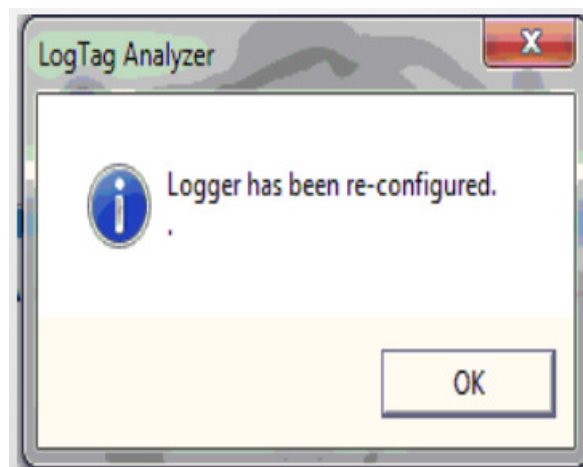


Downloading data

Starting The Logger

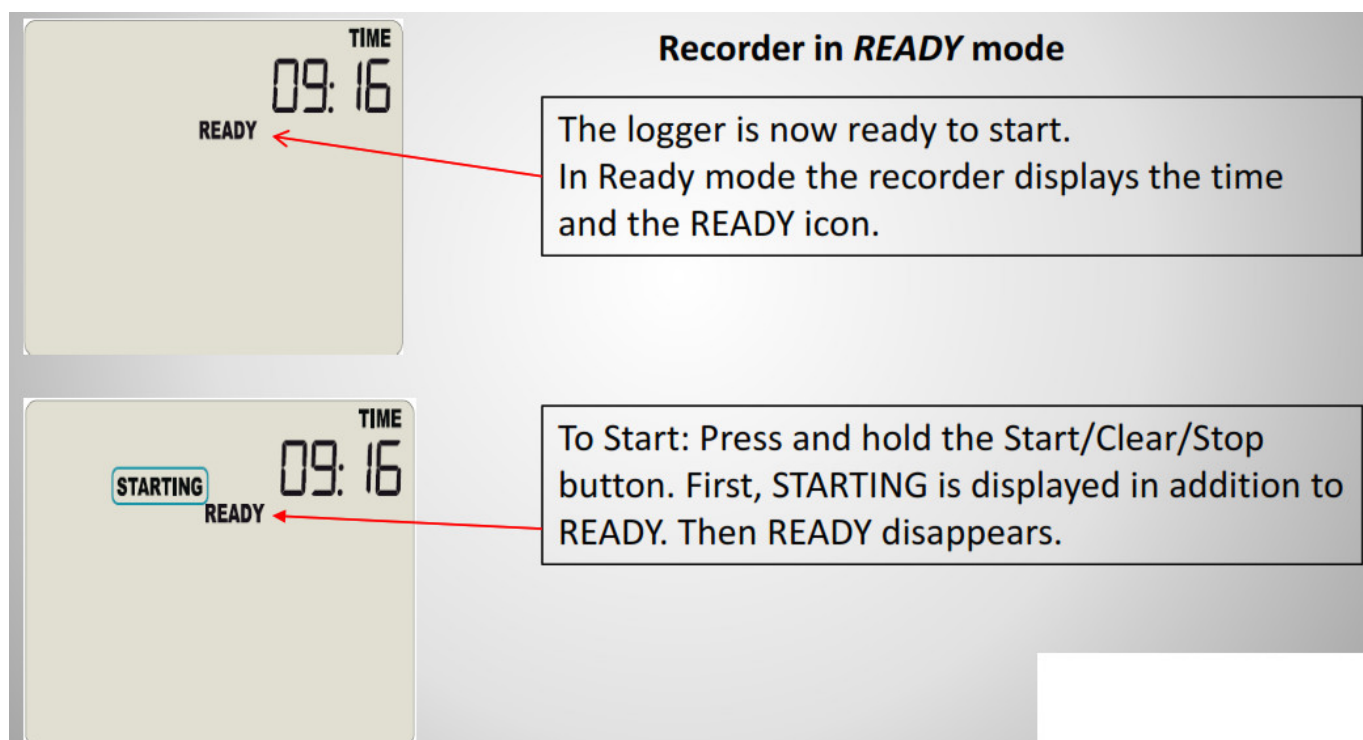


Downlaoding data Starting The Logger

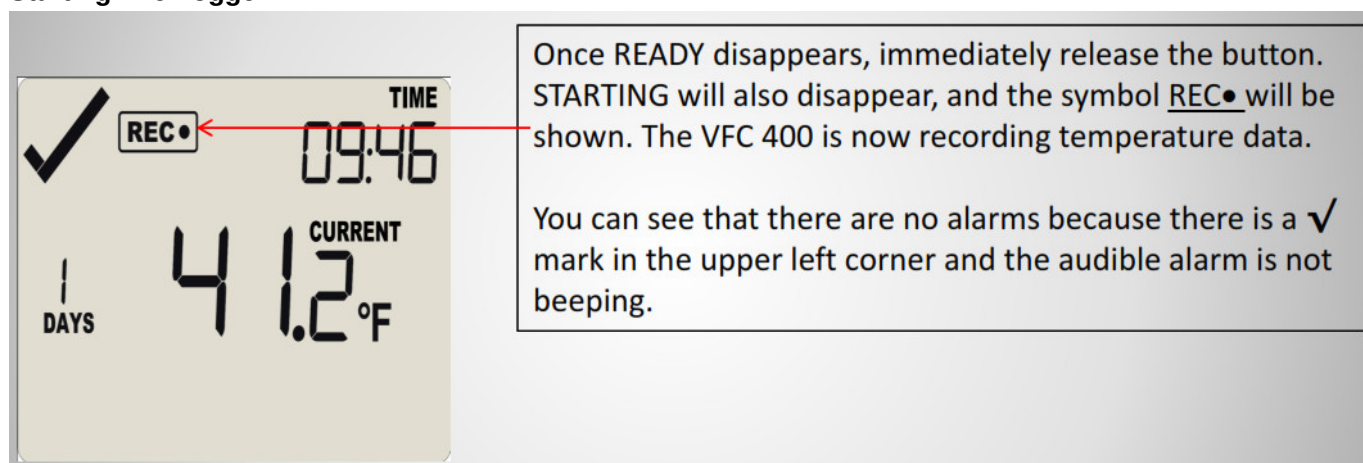

Control Solutions, Inc.
www.vfcdataloggers.com
 503-410-5996
 Thank you for your business
 Control Solutions, Inc.
 35851 Industrial Way, Suite D
 St. Helens, OR 97051
 503-410-5996
www.vfcdataloggers.com

Documents / Resources

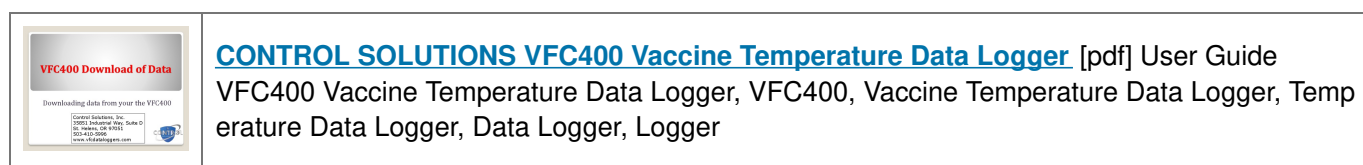

References

-  [Home](#) | [Control Solutions](#)

Manuals+.