

COREMORROW H63 6-Axis Motion Piezo Stage User Manual

Home » COREMORROW » COREMORROW H63 6-Axis Motion Piezo Stage User Manual








This document describes the following products:

• H63.XYZTR1S/K

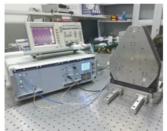





Contents

- 1 DECLARATION
- 2 Introduction
- 3 Safety Guide
- **4 Product Overview**
- 5 Operating environment
- 6 Installation and operation
- 7 Transportation and

inspection

- **8 Maintenance and Treatment**
- 9 Connector
- 10 Customer Service
- 11 Contact us
- 12 Documents / Resources
 - 12.1 References
- 13 Related Posts

DECLARATION

This user manual is applicable to H63 Six-Axis Piezo Stage, for your smooth and correct use of this product, please read this manual carefully before using it. Please follow the instructions in the installation and use process.

Improper operation may cause injury to the operator and/or damage to the product, so it is required that the entire installation and operation process should be completed by professionals with a certain foundation and a certain understanding of product principles, or under the guidance of a professional.

If you disassemble or transform the product without permission, our company will not be responsible for any consequences arising therefrom.

To change the product model and other false sales of our company's products are illegal, users should be vigilant. Once found, you should actively report or contact our company, jointly crack down on illegal acts, and prevent fraud and economic losses. CoreMorrow will investigate for the legal responsibility of the illegal subject and related party.

With the continuous development and innovation of technology,

CoreMorrow will timely add the latest information to the manual as required. If you need, please contact us, or to our website www.coremorrow.com for download, we apologize for any inconvenience caused to you.

Introduction

Purpose and content

- The manual is mainly for smooth and correct use of H63 piezo stage.
- The manual contains all the necessary information for H63 piezo stage.
- Introduction of the precautions during installation and use.

Symbol and meaning



DANGER: Improper handling may could cause injury to the operator.



Attention: Improper handling can cause damage to the equipment.

1. Please operating in order.

Users of This Manual

The default user of this manual has knowledge of piezoelectric products and controllers, and has a certain basis for their installation and use.

Manual Lost and Reissued

If this manual is lost, please contact CoreMorrow. We will send you the latest version of the user manual.

Safety Guide

H63 Six-Axis Piezo Stage is designed with the most advanced technology and safety standards for your own safety and the correct use of the product, please observe the following points when in use.

A change in temperature and pressure will charge the piezoelectric actuator, and the actuator will remain energized for a period of time when it is disconnected from the controller, So the operator should pay attention to when in use:

- Do not disassemble H63 Six-Axis Piezo Stage.
- Discharge H63 before installation, which can be done by connecting to the controller.
- Do not unplug the wire connecting with controller during operation.
- Before use, check whether the connection wire of the product is in good condition and whether the product and
 the controller are effectively grounded. During operation, strictly standardize the operation. Do not touch the
 product (up to 150V voltage) with your hand when power on, in case of danger.

In order to prevent damage to the product caused by improper operation, the following points should be paid attention to when in use.

- H63 Six-Axis Piezo Stage is used in a dust-free, oil-free and lubricant-free environment.
- As the product is designed with flexible structure, it is recommended that the load should not exceed load
 capacity of product. When loading, DO NOT apply torque force and DO NOT pry or twist the load surface, so
 as not to damage the structure.
- Recommended voltage range is 0~120V for long-term use and lifetime.
- Do not disassemble the product to avoid damage.
- Avoid stretching and bending the cable interface to prevent damage to the cable.

- Use the dedicated cable provided by CoreMorrow to connect H63 Six-Axis Piezo Stage to controller.
- Do not connect via an extension cable casually, if you need a longer cable, please contact CoreMorrow.

Product Overview

Product Introduction

H63 Six-Axis Piezo Stage has the following advantages:

High performance piezo stack:

H63 Six-Axis Piezo Stageis driven by a highly reliable ceramic actuator with a fully insulated structure, so its performance and service life are far better than the traditional actuator. The insulating layer can effectively prevent the piezoelectric ceramic from being damaged when exposed to moisture. Therefore, the stability of piezoelectric ceramics can be guaranteed under extreme circumstances. Compared with traditional driving mechanism, there are no transmission parts and friction. Therefore, ceramic actuator has the characteristics of no rebound, no maintenance, and no wear, etc.

Flexible hinge structure:

H63 Six-Axis Piezo Stage adopts flexible hinge mechanism has advantages of no friction, high positioning accuracy. Flexible hinge is an element not affected by static and dynamic friction, based on the elastic deformation of metal material (e.g., steel) , there is no rolling and sliding parts. Flexible unit has very high load capacity and stiffness. Hinge guide needs no maintenance and no wear and tear. Suitable for a wide range of temperatures, no lubricating oil is required

Strain Sensor

Strain sensor detects the position information by its own expansion. They are fixed in the proper position of the transmission part and measure the displacement of the moving part. This type of position measurement is made by contact and indirect measurement. So the displacement of the moving platform is measured by measuring lever, guide hinge, or piezoelectric ceramic stack.


The H63 can run up to 1mrad in $\theta X \theta Y \theta Z$ and $\theta \mu m$ in XYZ, and features fast response speed, high repeatable positioning accuracy, high flexibility, easy to operate, so it can meet different applications. This product can also be customized according to the needs.

Product Features

- Deflection Angle to 1mrad/axis, Travel rage to 6µm/axis.
- High precision sensor, high positioning accuracy.
- Sub-radian resolution.
- · Fast response.

Motion direction

FIG. 1 H63 Motion Direction

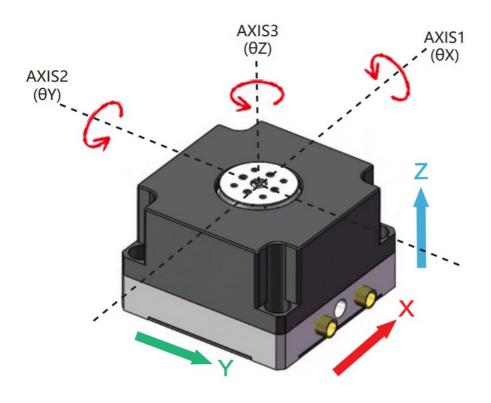

Model	Motion direction	Function	Label
H63	θΧ	PZT Drive	CH1
		Senser output	1
	θΥ	PZT Drive	CH2
		Senser output	2
	θΖ	PZT Drive	СНЗ
		Senser output	3
	Constant voltage	PZT Drive	CH4
	X	PZT Drive	CH5
		Senser output	5
	Υ	PZT Drive	CH6
		Senser output	6
	Z	PZT Drive	CH7
		Senser output	7

Operating environment

- A clean, dustfree and non-corrosive environment
- Operating temperature : -20~80°C
- Relative humidity: <60%
- Recommended voltage range is 0~120V for long-term use and lifetime


Installation and operation

Safety check should be carried out before installation, such as whether the connection wire is damaged or broken, whether the drive power supply is grounded, etc.

Installation and Precautions

Installation diagram of H63 is as follows

FIG. 2 H63 installation diagram

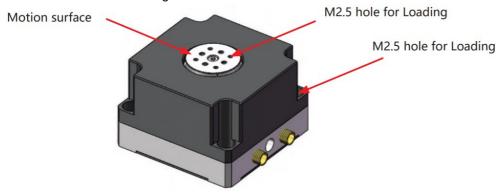


- 1. Ensure that the installation plane is relatively flat.
- 2. Use M4 screws to mount the H63 and tighten.
- 3. Mounting the load and tighten it properly to ensure the positioning accuracy.
- 4. Connect H63 connector with the controller.

Load

When mounting the load, ensure that the load quality is less than the maximum load capacity of H63, and try to mount the load in the center of the moving surface.

The correct position to mount the load is shown below.


FIG.3 The load is applied to the center of the moving surface

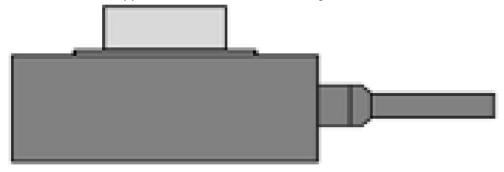

For the protection of the product, we do not recommend that the load is too high. If a high load must be used, please appropriately low the position of the center of gravity of the load.


The wrong mounting way is shown in the figure below:

FIG.4 Try not to let the position of the center of gravity far exceed the active platform

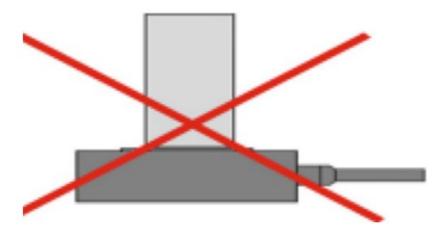

It is forbidden to mount the load on one side of the moving platform, which will cause serious damage to the product, as shown in the figure below.


FIG.5 Do not mount long loads on one side of the platform

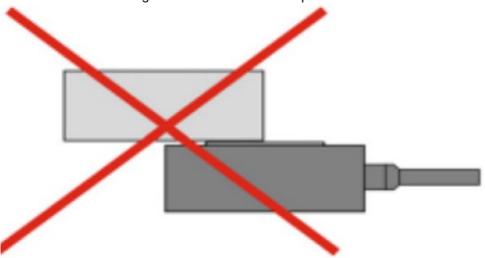

Excessive length of screws selected in the fastening and connecting process will lead to failure in installation or leakage of bolts, which will expose the product and scratch the product, affecting the motion accuracy of the product. Short bolts will lead to insufficient compactness of the connection and easy looseness, which will damage the internal structure of the product during operation.


We recommend using the connection as shown below:

FIG.6 The correct bolt connection

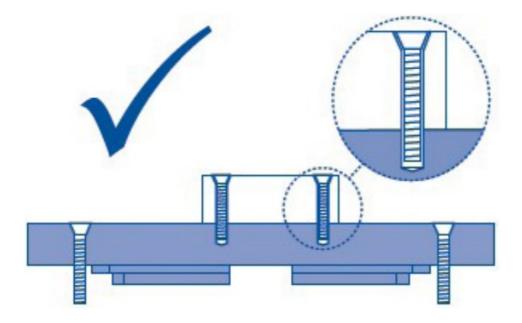

Please do not connect in this way as shown below:

FIG. 7 Inappropriate bolt selection

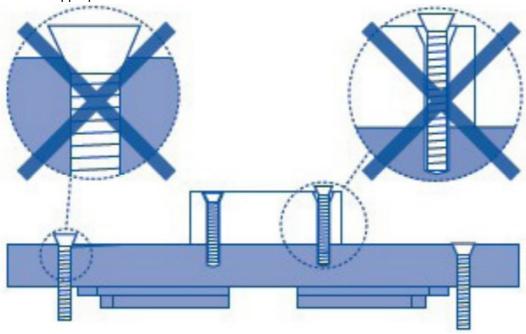

Notice:

- When mounting the load, tighten the load slowly. The torque should not be too large to prevent damage to the mechanism.
- · Avoid stretching and bending cables.
- Use the dedicated cable provided by CoreMorrow to connect H63 and the controller.
- Do not connect via an extension cable casually, if you need a longer cable, please contact CoreMorrow.

Operating

Please reconfirm the following information before turning on the power:

• The H63 has been grounded.

- The operating environment meets the operating environment requirements of H63.
- · Fixed screws must be tightened.

After the above information is confirmed, relevant controller operations can be carried out:

- 1. Please read the user manual of the controller carefully before operation, and then start the controller.
- 2. Set the controller' output voltage to zero first.
- 3. Check no abnormal phenomenon, then input voltage value (via keyboard oranalog control) according to specified input voltage range.
- 4. Operating normally within the rated voltage range of piezoelectric ceramics (0~120V, maximum 150V).
- 5. After use, please carry out voltage reset operation

Please take light steps when H63 in operation, do not knock the product, so as not to affect the use of the product or even damage the product.

In order to achieve better motion precision of H63, it is recommended to use our company's supporting controller. Please refer to the user manual of the controller for the specific operation to ensure correct operation before relevant operation.

Notes

- The driving voltage shall not exceed the voltage range of H63.
- If any abnormal noise or oscillation occurs during use, turn off the power immediately and check the parameter setting.

Transportation and inspection

Transportation

This product is packed in cartons. Transportation is carried out under the packaging conditions. This product can be transported by various means of transport under normal circumstances. In the process of transportation, direct rain and snow, corrosive substances, strong collision, extrusion, irregular placement and other improper behaviors should be avoided as far as possible.

Open-Box Inspection

Inspection content:

- Before opening the package, please check the outer packing carefully to see if there is any breakage, wetting, dampness, deformation, etc.
- Please check the completeness of package contents one by one according to the supply contract and packing list.
- Check whether the surface of the product has obvious damage
- Check whether the product identification is clear and complete.
- Check whether the connector is loose and the connector wire is broken or damaged.

In case of any of the above, please make a detailed record and take a photo of it.

Meanwhile, please contact us for processing immediately

Maintenance and Treatment

Problem Handling

FAQ are shown in the table below.

Problem description	Possible reason	Solution
	The connection wire is not connect ed properly or poor contact	Check the connection wire
No displacement or less displacem ent	Excessive load	Load according to the product's loa d capacity
	Sensor zero d	Check t
	Mountingsurface unevenness	Mount the platform on a flat surface
Low accuracy	Loose connection	Fasten connection
	Wrong connection	Check if the wiring is correct
Vibration or inaccurate positioning	Incorrect input control param	Stop immediately and check the pa rameter Settings
at startup	Resonance occurs	Make drive frequency less than res onance frequency

If the user cannot solve the problem by himself in the process of using, please record the problem in detail and contact CoreMorrow, and professional technical personnel will help to solve.

Problems caused by improper operation are not covered by the warranty.

CoreMorrow is not responsible for any problems caused by dismantling products without permission.

Daily maintenance

If the product is not used for a long time, please store it in a dust-free, well-ventilated, clean and non-corrosive indoor environment after packaging.

In the daily maintenance, please do not loosen the bolts on the product structure, so as not to affect the movement accuracy of the product.

Daily cleaning:

Please disconnect the product from the controller before cleaning the surface of the product.

When cleaning, gently wipe the surface of the product with a towel dipped in a small amount of detergent, and then clean it with alcohol. Please do not use ultrasonic cleaner.

When cleaning the surface, try not to pick up the product. If you have to pick up the product, it is better to put the

sponge under the product for protection. The product should not be too high from the sponge, in case the product falls off

Disposal

Waste products should be disposed according to national and local rules and regulations.

In order to fulfill our responsibility as a product manufacturer, we will dispose all old equipments on the market in an environmentally friendly manner. If you have equipment that cannot be disposed, you can ship it to CoreMorrow.


Address: 1F, Building I2, No.191 Xuefu Road, Nangang District, Harbin, Heilongjiang Tel: +86-451-86268790

Connector

Driving and Sensor Connector

H63 can be controlled by E51.D7S piezo controller that adopts DB15-HD (female) connector. H63's connector is DB15-HD (male), and the definition is as follows:

FIG.8 db15-HD (male) connector

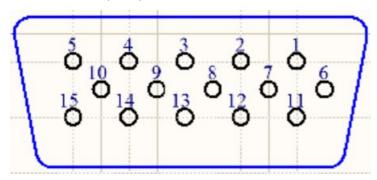

The pin definition of DB15-HD (male) connector(in left) is as follows:

No	Pin Definition
1	CH2 drive outpu
2	CH3 drive output
3	CH3 sensor input signal –
4	CH3 sensor input signal +
5	Sensing power supply:DC10V
6	CH1 drive output
7	CH2 CH3drive output GND
8	Constant voltage drive output
9	CH2 CH3 sensing power supply GND
10	CH1 sensing power supply GND
11	CH1 drive output GND Constant voltage output GND
12	CH2 sensor input signal
13	CH2 sensor input signal+
14	CH1 sensor input signal+
15	CH1 sensor input signal-

Table 3 Connector pin definition

The pin definition of DB15-HD (pin) connector(in right) is as follows:

No	Pin Definition
1	CH5 drive output
2	CH6 drive output
3	CH6 sensor input signal –
4	CH6 sensor input signal +
5	Sensing power supply:DC10V
6	CH4 drive output
7	CH5 and CH6 drive output GND
8	Do not connect any cables
9	CH5 and CH6 sensing power supply GND
10	CH4 sensing power supply GND
11	CH4 drive output GND
12	CH5 sensor input signal
13	CH5 sensor input signal+
14	CH4 sensor input signal+
15	CH4 sensor input signal

Customer Service

If you have questions about the products you are currently using, please let us know the following information:

- Product model and relevant number
- The controller model
- Software driver version
- Computer operating system that installs supporting software

Contact us

Harbin Core Tomorrow Science & Technology Co., Ltd.

Tel: +86-451-86268790

Email: info@coremorrow.com
Website: www.coremorrow.com

Address: Building I2, No.191 Xuefu Road, Nangang District, Harbin, HLJ, China

CoreMorrow Official and CTO WeChat are below:







Documents / Resources



<u>COREMORROW H63 6-Axis Motion Piezo Stage</u> [pdf] User Manual H63 6-Axis Motion Piezo Stage, H63 6-Axis, Motion Piezo Stage, Piezo Stage

References

· (0) - -

Manuals+,