



dynamic BIOSENSORS 10X BUFFER HE140 PH 7.4 Running Buffer User Manual

[Home](#) » [dynamic BIOSENSORS](#) » dynamic BIOSENSORS 10X BUFFER HE140 PH 7.4 Running Buffer User Manual 

dynamic
BIOSENSORS
heliX+
User Manual
10X BUFFER HE140 PH 7.4
running buffer
Dynamic Biosensors GmbH & Inc.
BU-HE-140-10 v2.1



Contents

- [1 Product Description](#)
- [2 Preparation](#)
- [3 Contact](#)
- [4 Documents / Resources](#)
- [4.1 References](#)

Product Description

Order Number: BU-HE-140-10

Table 1. Contents and Storage Information

Material	Composition	Amount	Storage
10x Buffer HE140 pH 7.4	100 mM HEPES, 1.4 M NaCl, 500 μ M EDTA, 500 μ M EGTA and 0.5 % Tween20; 0.2 μ m sterile filtered	50 mL	2-8°C

For research use only.

This product has a limited shelf life, please see expiry date on label.

Preparation

Dilute the complete solution 10x Buffer HE140 pH 7.4 (50 mL) by mixing with 450 mL ultrapure water.

After dilution HE140 Buffer is ready for use (10 mM HEPES, 140 mM NaCl, 50 μ M EDTA, 50 μ M EGTA and 0.05 % Tween20).

The diluted buffer should be stored at 2-8°C.

Contact

Dynamic Biosensors GmbH

Perchtinger Str. 8/10
81379 Munich
Germany

Dynamic Biosensors, Inc.

300 Trade Center, Suite 1400
Woburn, MA 01801
USA

Order Information order@dynamic-biosensors.com

Technical Support support@dynamic-biosensors.com

www.dynamic-biosensors.com

Instruments and chips are engineered and manufactured in Germany.

©2024 Dynamic Biosensors GmbH | Dynamic Biosensors, Inc. All rights reserved.

dynamic
BIOSENSORS

BU-HE-140-10 v2.1

www.dynamic-biosensors.com

Documents / Resources

	dynamic BIOSENSORS 10X BUFFER HE140 PH 7.4 Running Buffer [pdf] User Manual BU-HE-140-10, 10X BUFFER HE140 PH 7.4 Running Buffer, 10X BUFFER HE140 PH 7.4, Running Buffer, Buffer
---	--

References

- [HomePage | Biosensors International Ltd](#)
- [Home - Dynamic Biosensors](#)
- [Home - Dynamic Biosensors](#)
- [User Manual](#)

Manuals+ Privacy Policy

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.