

illumina v4.0.5 Convert Standalone

illumina SBS (HiSeq GAllx)	Total
Step 1 » Library Normalization (Illumina SBS)	0
Step 2 » Library Pooling (Illumina SBS)	0
Step 3 » qPCR QC	0
Step 4 » Cluster Generation (Illumina SBS)	0
Step 5 » Illumina Sequencing (Illumina SBS)	0
Step 6 » Bcl Conversion & Demultiplexing (Illumina SBS)	0



illumina v4.0.5 Convert Standalone User Guide

[Home](#) » [illumina](#) » illumina v4.0.5 Convert Standalone User Guide 

Contents

- 1 illumina v4.0.5 Convert Standalone
- 2 Product Information
- 3 FAQ
- 4 INTRODUCTION
- 5 FEATURES
- 6 RESOLVED ISSUES
- 7 Documents / Resources
 - 7.1 References



illumina v4.0.5 Convert Standalone

illumina SBS (HiSeq GAllx)	Total
Step 1 » Library Normalization (Illumina SBS)	0
Step 2 » Library Pooling (Illumina SBS)	0
Step 3 » qPCR QC	0
Step 4 » Cluster Generation (Illumina SBS)	0
Step 5 » Illumina Sequencing (Illumina SBS)	0
Step 6 » Bcl Conversion & Demultiplexing (Illumina SBS)	0

Product Information

Specifications

- Product Name: BCL Convert

- Version: 4.0.5
- Release Date: May 2, 2023
- Function: Converts per cycle binary data output by Illumina sequencers containing basecall files and quality scores to per read FASTQ files

Product Usage Instructions

Introduction

BCL Convert is a software tool designed to convert binary data output by Illumina sequencers into FASTQ files.

New Features

The latest release of BCL Convert includes new features that enhance the conversion process.

Resolved Issues in v4.0.5

This version addresses specific issues that were present in the previous release to improve performance and stability.

Known Issues

There may be some known issues listed in the release notes that users should be aware of while using BCL Convert.

FAQ

- **Q: What is the primary function of BCL Convert?**
 - A: BCL Convert is primarily used to convert binary data from Illumina sequencers into FASTQ files.
- **Q: How often are updates released for BCL Convert?**
 - A: Updates for BCL Convert are released periodically to address issues and introduce new features.

INTRODUCTION

These Release Notes detail the latest release of BCL Convert, including known issues. BCL Convert converts per cycle binary data output by Illumina sequencers containing basecall files and quality scores to per read FASTQ files.

FEATURES

NEW FEATURES

- Per tile and per cycle primary statistics output to report files
 - Adapter_Cycle_Metrics: provides, for each cycle and sample, the number of reads mapping to the sample where the adapter was detected beginning at that cycle.
 - Demultiplex_Tile_Stats.csv: includes all columns that exist in the Demultiplex_Stats.csv output file but at the per sample and per tile level

- Quality_Tile_Stats.csv: includes all columns that exist in the Quality_Stats.csv output file but at the per sample and per tile level
- Support Sample_Name column in the Sample Sheet Data section
 - Only allowed when --sample-name-column-enabled is enabled
 - Must be specified for every sample when enabled
 - Fastq filename will be named according to Sample_Name in sample sheet.
 - Fastq files will be output to subdirectory name by Sample_ID when --sample-name-column-enabled is enabled and --bcl-sampleproject-subdirectories is enabled
 - Reports will include Sample_Name and Sample_Project values when these features are enabled on the command line.
- Allow legacy FindAdaptersWithIndels sample sheet setting (default off)
 - Provides identical output to bcl2fastq2 2.20 for default adapter trim settings.
- Allow for no sample sheet to be provided using the --no-sample-sheet command
 - Default is disabled (sample sheet required)
 - When enabled, all sequences will go to the 'Undetermined' fastq files
- Cannot be enabled with any of the following options
 - bcl-sampleproject-subdirectories
 - sample-name-column-enabled
 - bcl-only-matched-reads
 - num-unknown-barcodes-reported
 - bcl-validate-sample-sheet-only
 - sample-sheet
- Can specify fastq gzip compression level
 - Command line option is --fastq-gzip-compression-level and default is 1
 - 0 through 9 are allowed to be specified
- Corruption detection for BCL input files
 - In strict mode, will abort conversion with an error
 - In robust mode, will skip processing lane

RESOLVED ISSUES

RESOLVED ISSUES IN V4.0.5

- Fix a crash that can occur when using per-sample-settings with higher sample counts in a lane, due to a hash-table pre-size using a signed integer as input that is overflowed
- Fix BCL FASTQ file paths in the fastq_list.csv when ORA-interleaved compression format is used. Previously the fastq_list.csv file contained two files (instead of one single interleaved file) under the "Read1File" and "Read2File" columns and the files were not named correctly.
- Some versions of RTA3 outputs CBCL with 0 qual + nonzero base for 0/"N#". In other words, masked nibbles (0 qual) do not zero out the base. bcl2fastq2 has a masking step, but DRAGEN™ and bcl-convert did not. This fix adds masking so that DRAGEN™ and bcl-convert matches bcl2fastq2 for those RTA3 outputs.
- Fix for no-sample-sheet setting that omits index sequences from fastq headers.
- Fix to remove BCL conversion thread settings limit of 64, a regression in v3.10, to allow runs on high core count systems.
- Fix a crash due to threading error on a 150k sample dataset.

- Fix for Index_Hopping_Counts.csv containing incorrect Same_{Name,Project}.

RESOLVED ISSUES IN V4.0.3

- tiles command line option did not work correctly with bgzf inputs, resulting in fastq files with incorrect tile header. This will now be correct.
- Barcode Collision Error and Solution: Previous versions of BCL Convert allowed a conversion to continue if either index (i7 or i5) had sufficient hamming distance from all other samples in the lane. BCL Convert 3.10.5 uses stricter barcode collision logic to support increased high-throughput and complex sample pooling.
- Each index in a dual setup must individually meet the hamming distance requirements set by the BarcodeMismatchesIndex# value. If either i7 or i5 does not meet the hamming distance requirements the program will error. For default BarcodeMismatchesIndex1 and BarcodeMismatchesIndex2:
 - Barcodes must differ by at least three bases.
 - If any two samples in i7 differ by fewer than 3 bases, an error is produced and the run will not proceed, regardless of their i5 values.
 - If any two samples in i5 differ by fewer than 3 bases, an error is produced regardless of their i7 values.
- If you receive errors with current versions of DRAGEN or BCL Convert, lower the mismatch tolerance for the index producing the error by using the BarcodeMismatchesIndex1 or BarcodeMismatchesIndex2 sample sheet settings.

KNOWN ISSUES

- If a directory is specified as input to ‘–sample-sheet’, BCL Convert will hang at the beginning of a run while trying to copy that path as a file to <outdir>/Reports/SampleSheet.csv
- BCL Convert does not validate when “Logs” or “Reports” is provided for a Sample_Project, and the software will be unable to create the subdirectories if these string are provided.
- BCL Convert does not support the –first-tile-only option being specified for SP flow cells, but the new –tiles option can be used as a substitute.
- Does not error when no tile list exists in the RunInfo.xml file and –tile or –exclude-tiles is specified in the command line
- ORA Compression cannot be specified on the command line when the –no-sample-sheet option is used. Therefore gzip compression is the only type of compression supported when no sample sheet is provided

RELEASE HISTORY

Revision	Release Reference	Originator	Description of Change
00	1085822	Daniel Tracy	Initial release

Template, Software Customer Release Notes, Doc. No.: 15048849, Rev. 01 Effective Date: 23-Jan-2023
Proprietary & Confidential

- BCL Convert v4.0.5
- Customer Release Notes

- Document ID: 200038045_v00
- Release Date: May 2, 2023 Page 4 of 4

Documents / Resources

	illumina v4.0.5 Convert Standalone [pdf] User Guide v4.0.5 Convert Standalone, v4.0.5, Convert Standalone, Standalone
---	--

References

- [User Manual](#)

[Manuals+](#), [Privacy Policy](#)

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.