



JOYALENS JL246MS 3 Lens LCD HDMI Digital Microscope Instruction Manual

[Home](#) » [JOYALENS](#) » JOYALENS JL246MS 3 Lens LCD HDMI Digital Microscope Instruction Manual 

JOYALENS JL246MS 3 Lens LCD HDMI Digital Microscope Instruction Manual



Contents

- [1 Components list](#)
- [2 Assembly steps](#)
- [3 Documents / Resources](#)
 - [3.1 References](#)
- [4 Related Posts](#)

Components list

Thank you for choosing our products. Please read the assembly steps and user manual carefully before use. If you have any problems with the use and assembly, please feel free to contact us. We will reply within 24H (there may be a delay on holidays). We appreciate the opportunity to serve you, your satisfaction is our utmost priority!


1. @ Microscope Monitor



2. @ Base



3. Beam

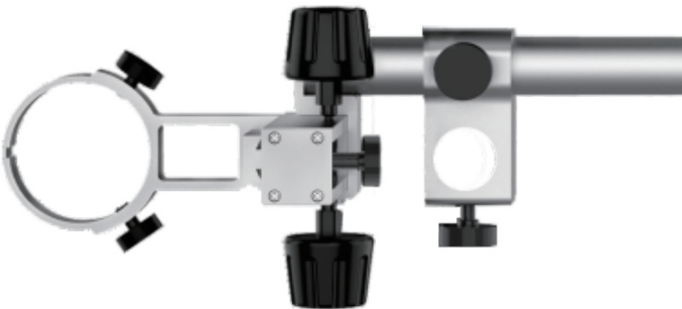


4. Column



5. Screw x2



6. Spring x2

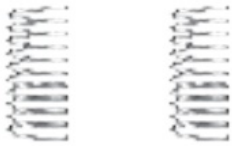



7. Metal clip x2



Assembly steps

Step 1

Assemble 4 to 2.

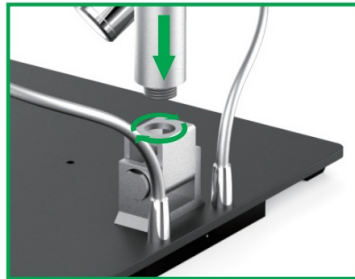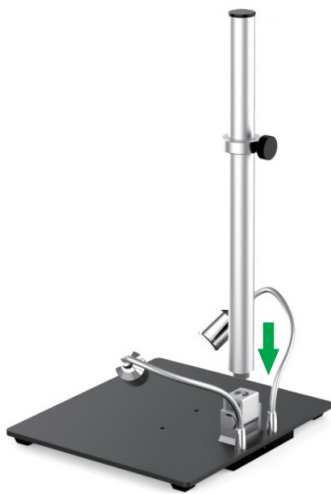


Step 2



(2.1)

Assemble 3 to 4.

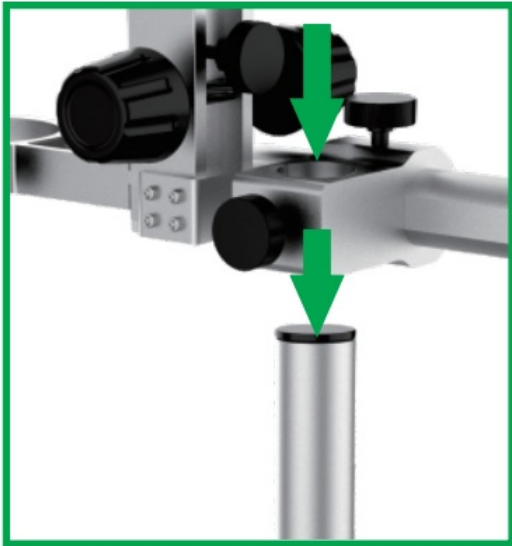

(2.2)
Tighten the screws.



Step 3



Unscrew the screws and remove the lens.) tighten the screws. *Be careful not to drop the lens.




Step 4



(4.1)
Insert the Protrusion into the slot.




(4.2)
Assemble the lens and tighten the screws.

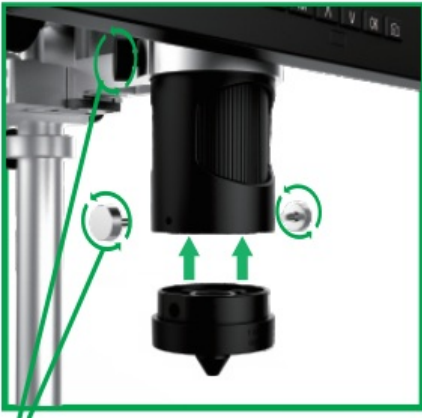


Step 5



Metal clips assembly step.

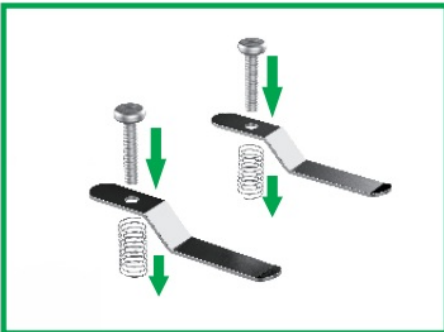

Assemble the clips to the base if necessary.

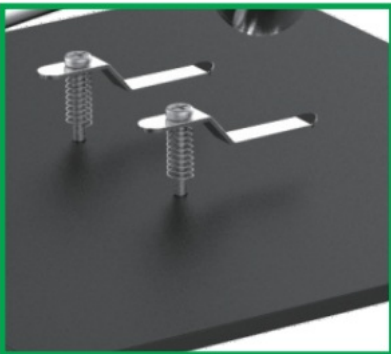


Step 6



When observing with lens D, slot in the DC power supply.



Documents / Resources

	<p>JOYALENS JL246MS 3 Lens LCD HDMI Digital Microscope [pdf] Instruction Manual JL246MS, JL246MS 3 Lens LCD HDMI Digital Microscope, 3 Lens LCD HDMI Digital Microscop e, LCD HDMI Digital Microscope, HDMI Digital Microscope, Digital Microscope, Microscope</p>
---	--

References

- [User Manual](#)