



# Kingfar BioSensor ErgoLAB Wireless System Instructions

[Home](#) » [Kingfar](#) » Kingfar BioSensor ErgoLAB Wireless System Instructions 

## Contents

- [1 Kingfar BioSensor ErgoLAB Wireless System](#)
- [2 ErgoLAB wireless physiological recoder](#)
- [3 Core advantages](#)
- [4 Parts list](#)
- [5 Manner of wearing](#)
- [6 BioSensor Specific operation](#)
- [7 Installation](#)
- [8 BioSensor Specification param-](#)
- [9 EXGSensor Specific operation](#)
- [10 EXGSensor Specification parameter](#)
- [11 Data collection](#)
- [12 FCC STATEMENT](#)
- [13 Documents / Resources](#)
- [14 Related Posts](#)


# Kingfar

**Kingfar BioSensor ErgoLAB Wireless System**



## ErgoLAB wireless physiological recorder

The new generation of ErgoLAB wireless human factors recorder includes Biosensor sensor and EXGSensor sensor types. All hardware adopts modular design, built-in Type-C interface, and unified collector and different signal source acquisition lines are configured to realize synchronous lead of multiple systems. At the same time, the equipment built-in acceleration, gyroscope magnetometer, in the collection of EMG, RMS, ECG, EDA, RESP SKT, SpO2, HR and other multidimensional physiological signals, but also can obtain high-precision human motion posture data. The product has the characteristics of intelligence, high precision and portability, which can meet the diverse needs of laboratory research, real field research and various simulated cabin environment research.

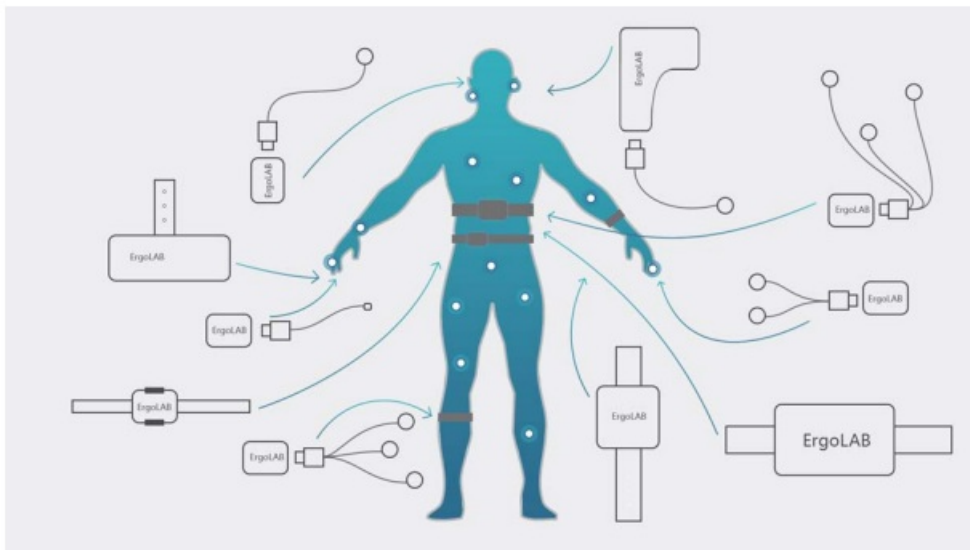

BioSensor wireless human physiological recorder adopts signal self-identification technology and is equipped with a unified collector. According to research requirements, the collection line of corresponding physiological signals can be arbitrarily connected. The equipment includes PPG blood volume pulse, EDA skin electricity, RESP respiration, SKT skin temperature, FSR fingertip pressure and other indicators, each sensor can collect multiple signal indicators.

You can complete the measurement mode of any index by adding sensors and electrode wires at any time.



EXGSensor wireless human factor sensor device adopts signal self-identification technology and is equipped with a unified EXG collector and acquisition line. According to research requirements, it can be equipped with EMG/ECG/EOG/EEG signal types, supporting patch mode and external acquisition line mode. The sampling frequency is up to 4096Hz, which can realize synchronous lead of multiple systems. At the same time, built-in acceleration, gyroscope, magnetometer, can obtain the human body's motion posture data.




## Core advantages

- Adopting multi-signal and multi-sensor fusion design, each single sensor can collect multiple data indicators at the same time.
- The equipment selects high-precision components, the system resolution can reach 16-24bit, single-channel sampling rate is the highest 4096Hz
- One to many data transmission protocol, using Bluetooth 5.0 ultra wideband transmission, can collect multiple devices at the same time;
- All products adopt modular design, support intelligent insertion detection signal Type, support type-c external sensor,
- The new sensor is designed to be portable and wearable. Including a band, buckle, bracelet, ear clip, finger, chest strap six wearing mode
- All sensors are uniformly calibrated for NTP network time before delivery to ensure time synchronization of all data collection and eliminate clock differences between different sensors and different computers. At the same time, the sensor and the system will automatically real-time synchronous calibration;
- New appearance design. Material and process upgrade to ensure product comfort and beauty
- Datalogger APP PAD, PC multi-terminal device collection,

- With marking function. t can be controlled by the subjects Treely and mark the time at a certain test stage to help better data analysis

## Parts list

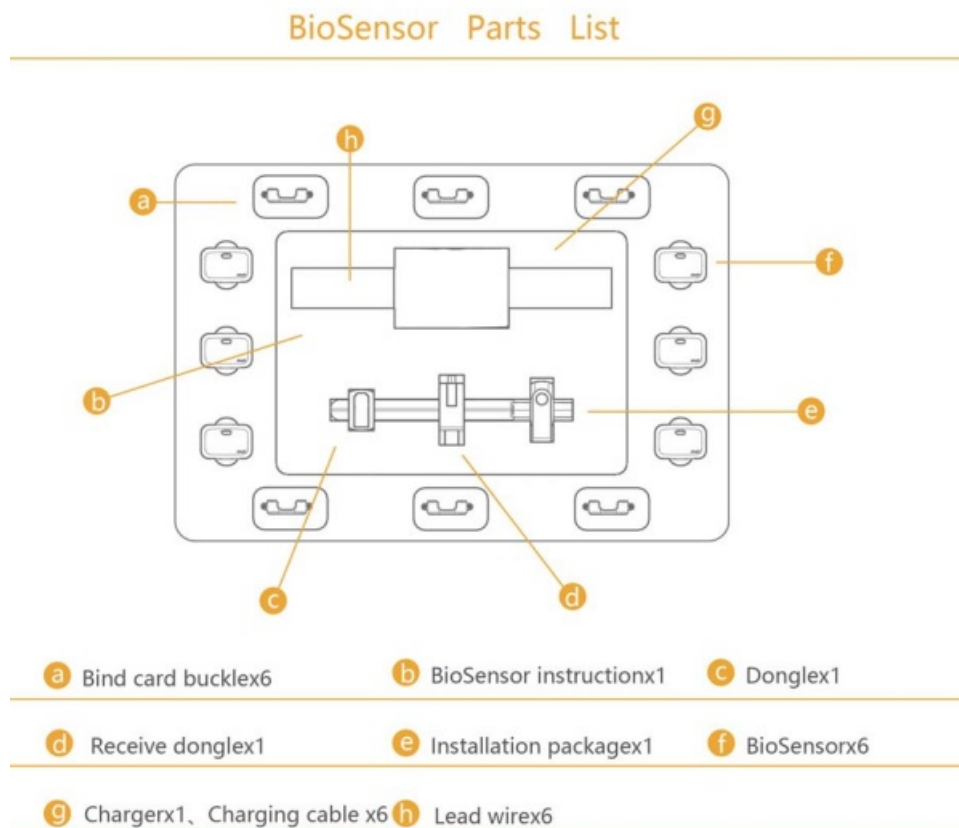

## Manner of wearing

1. Exg-patch acquisition mode: three silver/silver chloride electrodes are embedded in the sensor, which only need to be fixed on the skin surface of human muscles with double-sided adhesive to remove the influence of cables, especially suitable for fine research of small muscle groups; New appearance design. Material and process upgrade to ensure product comfort and beauty:
2. BIO, EXG-Type-C external sensor acquisition: the system can automatically insert detection, customize the electrode measurement position, support the long-distance test of large muscle groups, better adapt to the requirements of different muscle group position data acquisition.





## BioSensor Specific operation

BioSensor process includes four main steps: installing sensor, wearing sensor, starting sensor and connecting software for data collection.


Such as:

### Wear way

### Installation

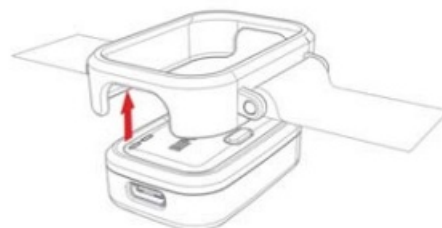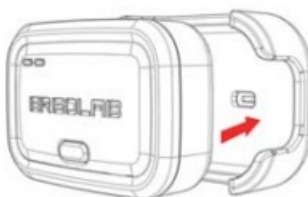


The sensor is fixed in the card slot, and the card slot is connected with the strap by means of penetration

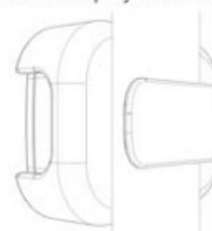

The sensor is fixed in the card slot, and the card slot and the strap are connected by a clamp

### Wear (wireless wear)

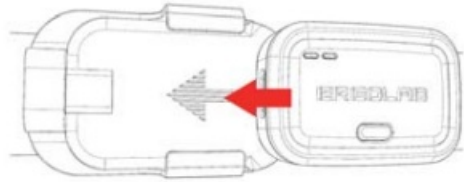

### Wear (wired wear)

Connect the sensor to the lead wire




The heart rate ear clip is inserted into the sensor, and the sensor is fixed in the card slot, and the card slot is connected with the strap by means of insertion



Push the sensor into the card slot, and the card slot is directly connected with the strap

Worn in the appropriate position of the body

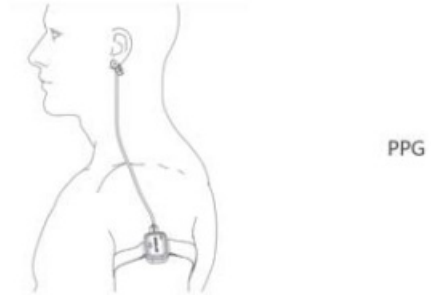

Sensors are worn on the upper arm and heart rate ear clips are clipped to the ears


### Wear (wired wear)

Connect the sensor to the lead wireThe heart rate ear clip is inserted into the sensor, and the sensor is fixed in the card slot, and the card slot is connected with the strap by means of insertion

### BioSensor Specific operation



Insert the lead wire with electrode buckle into the sensor,  
and the sensor is fixed in the clamp.  
The clamp is connected with the strap.

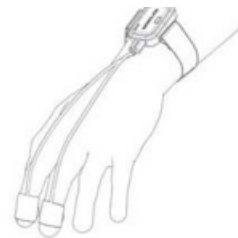


EDA

The sensor is worn on the wrist and the  
electrode buckle is fixed on the finger




Insert the lead wire with electrode buckle into the sensor,  
and the sensor is fixed in the clamp.  
The clamp is connected with the strap.

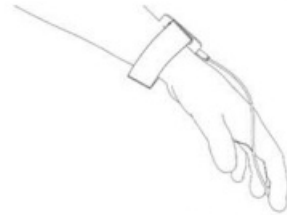


SKT

The sensor is worn on the wrist. Attach the  
collecting end to the finger

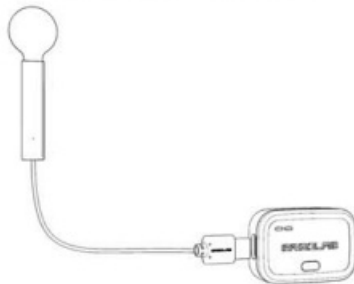

Insert the lead wire with electrode buckle into the sensor,  
and the sensor is fixed in the clamp.  
The clamp is connected with the strap.

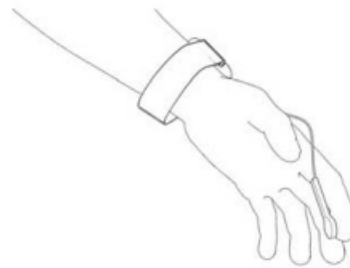

FSR


The sensor is worn on the wrist. Press or  
pinch the sensor

## Start Connection


## BioSensor Specification param-

## BioSensor Specification parameter

BIO Sensor Acceleration and magnetometer axis orientation. As shown in the figure below (Take Type-c to the left as the front):

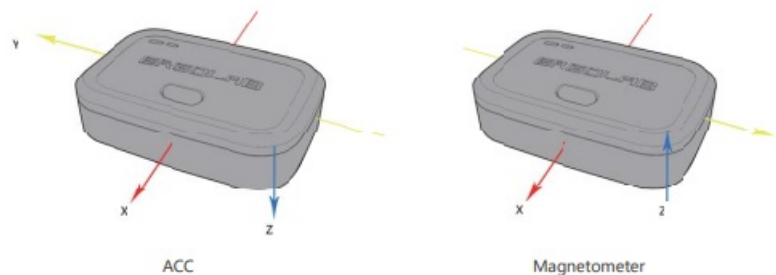

EXG Sensor Acceleration and magnetometer axis orientation. As shown in the figure below (Take TvDe-c to the left as the front):

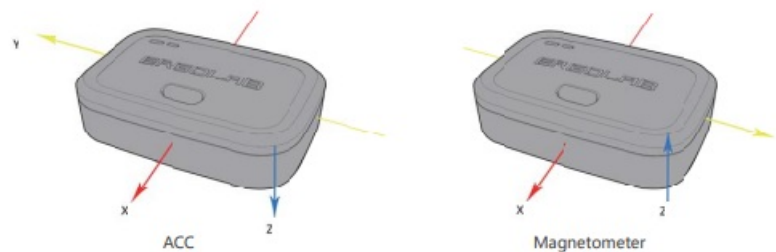


## EXGSensor Specific operation

EXG Sensor process consists of four main steps: installing the Sensor, wearing the Sensor, starting the Sensor and connecting the software for data collection.

AS shown below:

### Wear way (Wireless wear)

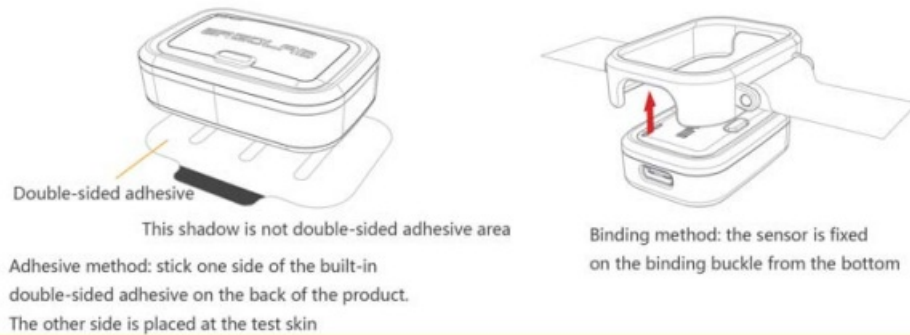


## Installation

### Fixed



### Connection Start

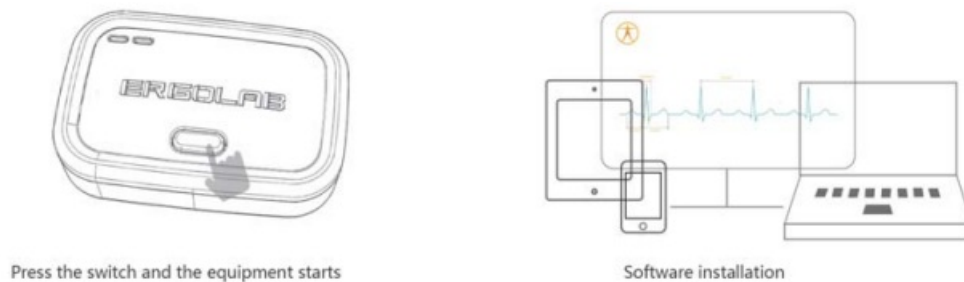

### Other way (Cable wear)



## EXGSensor Specification parameter

### Indicator st status










The following describes the status of sensor LED lights in use, and the corresponding meaning:

#### Battery indicator light:

 >60%    59%~40%    59%~40%    59%~40%    >60%

When charging, the battery light flashes at low frequency and the color is determined by the electric quantity. After charging, the green light is steady on, and the Bluetooth light is off. After startup, the battery light will be on every 5s and off for the rest of the time:

Bluetooth light color	Bluetooth light status	Meaning
	Put out	When the device is in charge mode, the Bluetooth light is off
	Normally on	(Wired) The device is disconnected after being turned on
	High frequency blinking	(Bluetooth) After the device is turned on, it is in broadcast state
	Normally on	It is connected to the receiving end
	Low frequency blinking	After the software is connected, data is reported
	Low frequency blinking	The collection end is abnormal
	Power off after steady on	When the power is less than 3%

#### PS:

1. Press on key equipment to start
2. Long press the key device to shut down
3. After startup, press the open key to mark the event
4. Bluetooth indicator on the left and battery indicator on the right
5. If no data is transmitted after the device is started, the indicator is in blue
6. After the device is started and data is being transferred, the indicator is green
7. When the Bluetooth indicator turns orange and flashes at low frequency, it indicates that data collection is abnormal. Check whether the electrodes are properly connected
8. If the device is not connected 3 minutes after startup, it will be in low-power mode. At this time, press the on

key again to turn on

## Data collection


### ErgoLAB synchronization test cloud platform acquisition system

ErgoLAB is independently developed by Jifa Technology, which can be applied to scientific research, personnel status identification, product design evaluation, and intelligent measurement of the human-information systems. Professional and efficient cloud architecture comprehensive platform in the direction of rating. It includes subjective and objective experimental design, multi-equipment data acquisition and data analysis Statistical report in one and build a complete test process system.



### ErgoLAB Datalogger APP portable acquisition system.

Datalogger APP enables real-time collection of all physiological data without any cables or computers, making it the best choice for outdoor research. At the same time of physiological data collection, temporal and spatial trajectory data can be collected synchronously, and then imported into ErgoLAB system for physiological signal analysis and temporal and spatial analysis, supporting synchronous analysis and comprehensive statistics with other multimodal data.

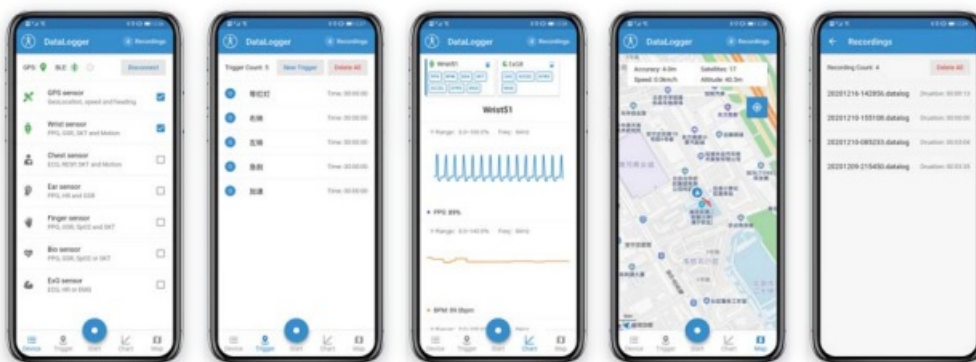

## FCC STATEMENT

This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) this device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation. Any changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.


**NOTE:** This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses, and can radiate radio frequency energy

and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- client or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

The device has been evaluated to meet general RF exposure requirements. The device can be used in portable exposure conditions without restriction FCC ID: 2ADKA-SENSOR

## Documents / Resources

	<a href="#">Kingfar BioSensor ErgoLAB Wireless System</a> [pdf] Instructions SENSOR, 2ADKA-SENSOR, 2ADKASENSOR, BioSensor ErgoLAB Wireless System, BioSens or ErgoLAB, Wireless System
---	--

[Manuals+](#).