




KoPa HW1200 Binocular Stereo Microscope Camera Solution Smart Embedded Camera User Manual

[Home](#) » [kopa](#) » KoPa HW1200 Binocular Stereo Microscope Camera Solution Smart Embedded Camera User Manual 



KS096500-Ver1.0
Binocular Stereo Microscope Camera
Solution (Smart Embedded
Camera User Manual

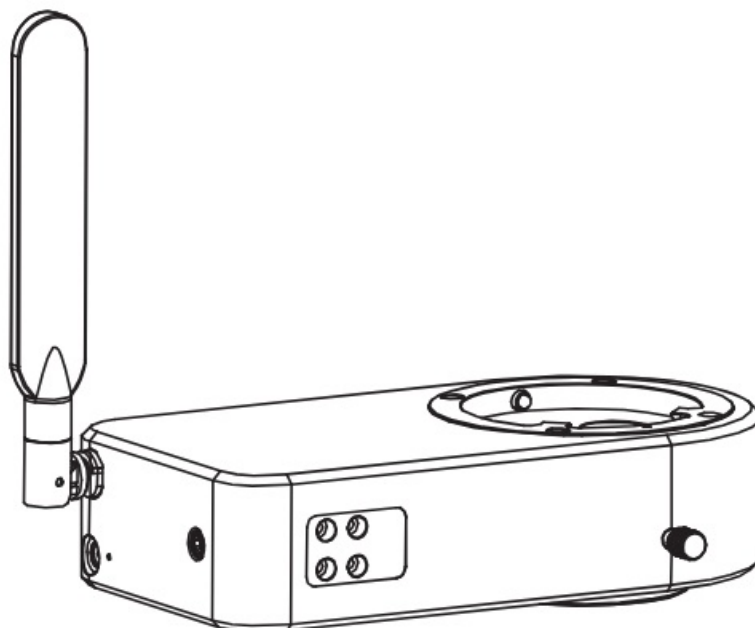

Contents

- 1 HW1200 Binocular Stereo Microscope Camera Solution Smart Embedded Camera
- 2 Notes And Safety Requirements
- 3 Package List
- 4 Hardware Introduction
- 5 Installation Methods
- 6 Connection Methods
- 7 Troubleshooting And Frequently Asked Questions
- 8 Disclaimer
- 9 Documents / Resources
- 9.1 References

HW1200 Binocular Stereo Microscope Camera Solution Smart Embedded Camera

Applicable brand	Olympus	Nikon	Leica	Olympus	Nikon	Leica
Model	HW1200	HW1200-M	FS12		FS20	
Category		N/A	L0	A0	N0	L0
Physical resolution		12.0MP			20.0MP	

Notes And Safety Requirements

Assemble

1. To avoid danger or damage incurred to the lens, do not touch the lens or sensor directly with your fingers.
2. To avoid failure or electric shock hazard and so on, do not disassemble or modify the internal structure of the device.
3. Do not plug in or unplug the USB port or the HDMI port when hands are wet.
4. Do not use alcohol and other organic solvents to clean.
5. If the lens or sensor is dirty or damp, you should better use dry and non-linen fabric or professional lens tissue to wipe them.
To avoid scratches on the surface, do not touch the lens with your fingers. Wipe the lens or sensor lightly.
6. The products are not specifically designed for an outdoor use. Do not expose it to outdoor environment without any protection.
Excessive temperature and humidity will damage the lens. Please avoid using the product under the following environment: high temperature or high humidity environment, places with direct sunlight, dirt or vibration and places near heat source.
7. Please use and store in the following environment:
Operating temperature 0 °C~ 40 °C;
Storage temperature -10 °C~ 60 °C;
Operating Humidity 30~60%RH;
Storage Humidity 10~80%RH.
8. If any foreign matter, water or liquid enter into the device by accident, disconnect the USB cable immediately.

Please send it to the maintenance center and do not use the hair dryer to dry it by yourself.

9. To prevent microscope from being tripped over or dropped, please put away the device's USB cables in use or standby.
10. To avoid electric shock by accident, please power off microscope before you move your Windows PC or laptop.
11. The cleanliness of the device lens will directly affect clarity degree of contents from the computer screen during preview.

Problems like various circles or spots on the screen may mostly be incurred by dirt on the lens. When cleaning, please use professional lens tissue or other professional detergent to clear the dirt on the lens.

12. Registered trademark and copyright information: this product is copyrighted by GUANGZHOU OSTECH ELECTRONIC TECHNOLOGY CO.,LTD. Without written authorization of the company, any organization or individual shall not copy, print or translate any part of this article into another language.

Package List

1. Smart embedded camera



2. HDMI cable(2m)



3. Wired mouse and keyboard



4. Allen key(3mm)



5. Power adapter and power cord (Optional Chinese, American, European, Australian, Korean, British standard, etc.)



6. One paper with QR code and link to download software

Software Manual Download

Windows download link:

<https://ostec.com.cn/downloads/>



<https://ostec.com.cn/downloads/>

Windows Download



For 5.0/8.0MP



For 12.0MP



KoPa WiFi Lab For Android/iOS Download KoPa WiFi Lab


<https://ostec.com.cn/downloads/kopa-wifi-lab>



Software Manual Download

<https://ostec.com.cn/downloads/software-manual>

Hardware Introduction

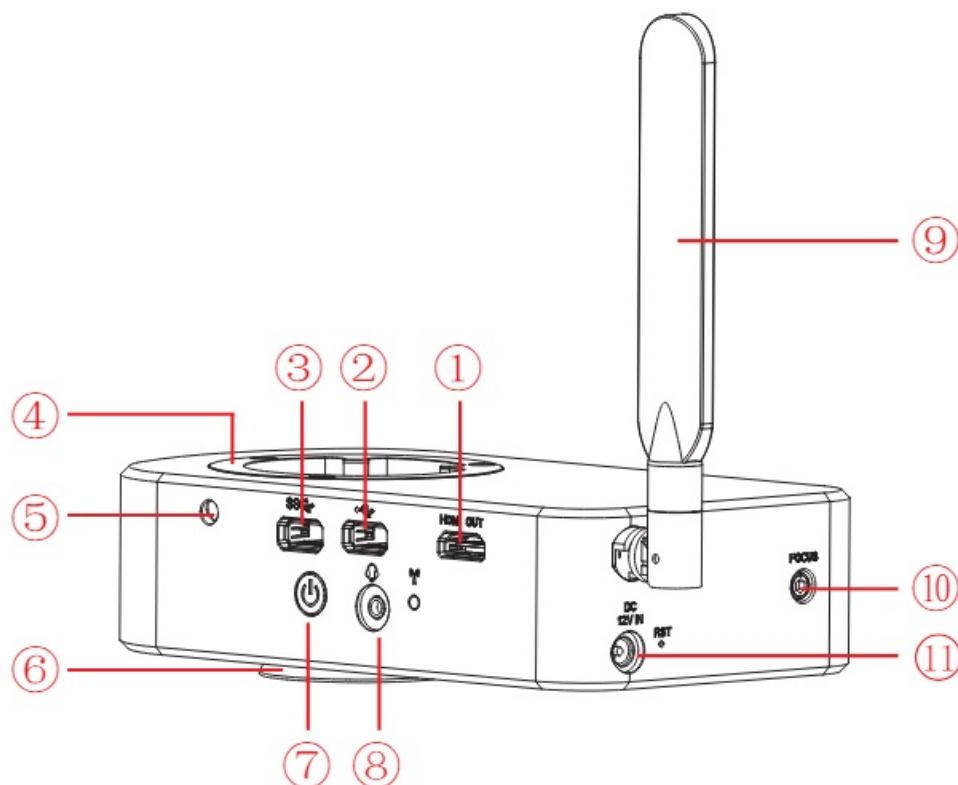

1	HDMI output interface	through HDMI cable, connect with external monitor with HDMI interface to realize video, signal transmission, etc.
2	USB 2.0 interface	can be connected to a mouse, keyboard, U disk, convenient for the use of software.
3	USB 3.0 interface	can be connected to a mouse, keyboard, U disk, convenient for the use of software.
4	Female interface	connects to the eyepiece end of the binocular microscope.
5	Tighten the screws	use the accessory Allen key to tighten the screws and fix the eyepiece of the microscope.
6	Male interface	connected to the objective end of the binocular microscope.
7	Power switch	switch on and off. Keep press for 4 seconds to power on / off.
8	Headphone and microphone interface	connect with headset cable to achieve transmission of audio signals.
9	5G WiFi antenna	enhance the 5G WiFi signal transmission, to achieve wireless connection to the camera for image acquisition.
10	Easy focus hole	internal filament focus structure is used to adjust the 0.43X field lens of the camera to achieve par-focal with electronic imaging and observation under eyepiece.
11	Power input interface (DC 12V 5A)	connect with the adapter in the accessories to achieve power supply to the camera.

Installation Methods

, A_k cp_ gl qr_jjrm_ gl masj_pk gapmqamne

1. Separate the objective end of the binocular microscope from the eyepiece end, as shown in Figure 1:
 - 1) Loosening the fixation screws of the binocular microscope;
 - 2) Remove the eyepiece end of the microscope.
2. Installation for smart embedded camera:
 - 1) Snap the smart embedded camera to the objective end of the microscope, oriented as in Figure 2;
 - 2) Tighten the fastening screws on the objective end of the microscope, fix the smart embedded camera on the microscope.
3. Installation for the objective end of the microscope, as shown in Figure 3:
 - 1) Snap the eyepiece end of the microscope to the female interface end of the smart embedded camera;
 - 2) Tighten the fastening screws of the smart embedded camera and fix the eyepiece end of the microscope on the smart embedded camera.

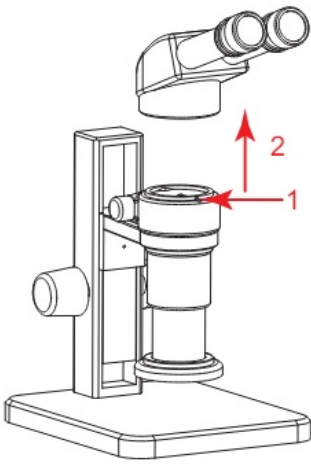


Figure 1

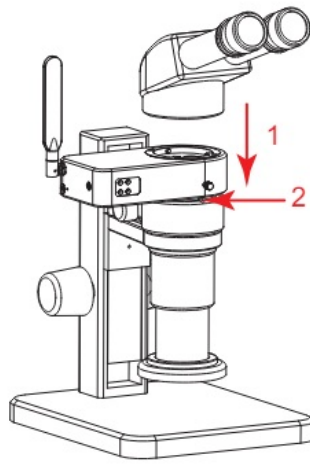


Figure 2

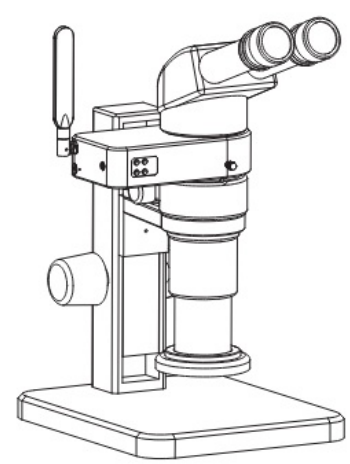


Figure 3

, Pck mtc a_k cp_dpmk k gapmgamnc


Reverse the order with the installation.

, N_ptdna _j

1. Place a flat specimen (such as a business cards, etc.) on the microscope's carrier table, adjust the diopter of the eyepiece to 0, adjust the microscope's magnification to the minimum, and adjust the microscope's objective distance to the clearest visualization;
2. Adjust the focal length in clockwise or counterclockwise direction with the allen key that we provided until the image from monitor is clear, you can realize parfocal in this way.

Note: as soon as parfocal is completed on the minimum magnification , the other magnification is basically free of adjustment and achieve parfocal as well.

Connection Methods

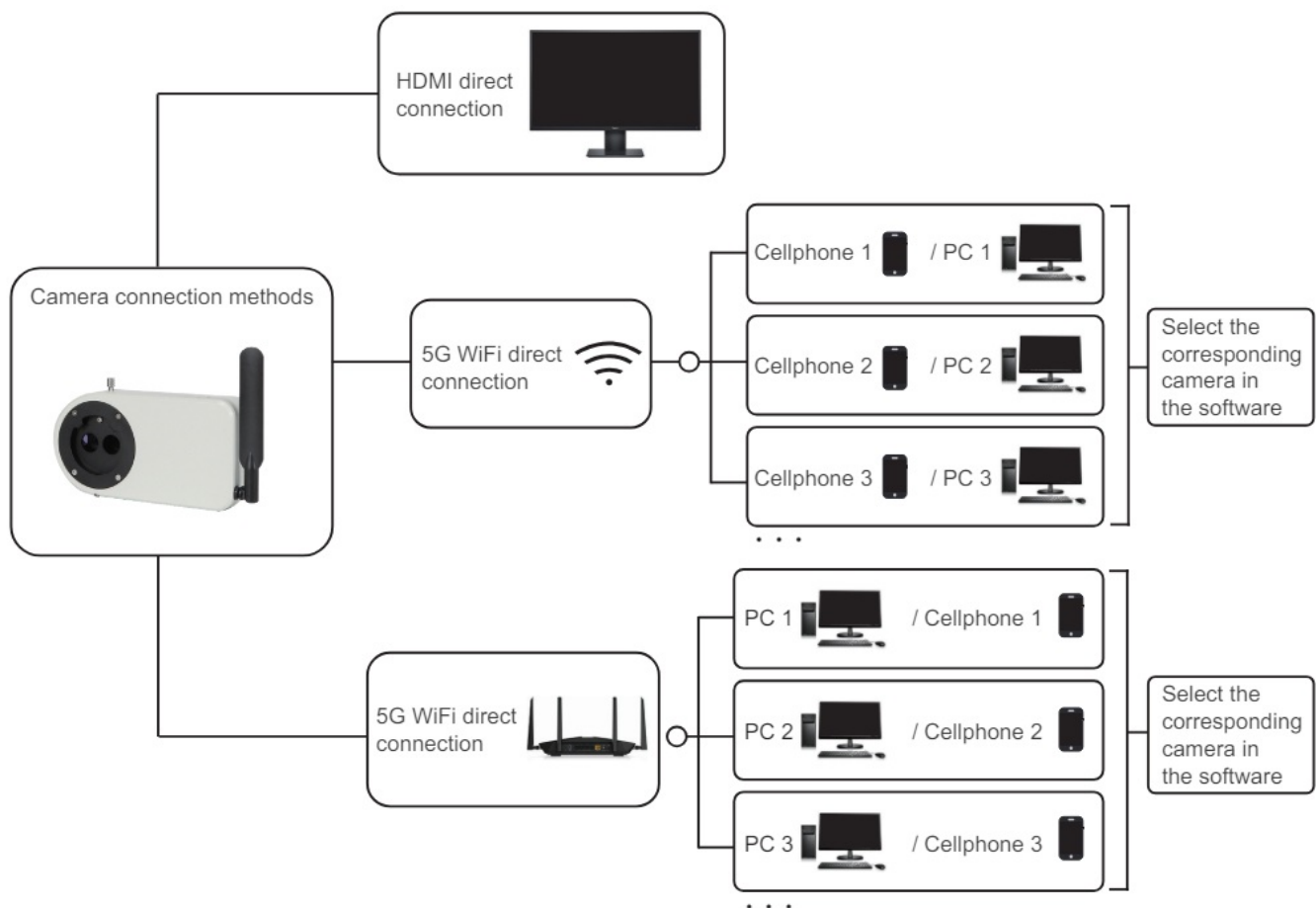

, FBK Gaml I cargml

(1) HDMI-connected display to be provided by user, recommended Dell 24°27-inch high color gamut display.
 (2) Since the smart camera has a built-in Android operating system, the software KoPa WiFi Lab AO will be displayed on the monitor to control the smart camera after the external monitor is connected to the camera via HDMI. Support HDMI and 5G WiFi output at the same time.

, EU gDganl I cargil

Desktop computers, cell phones, tablets, and laptops can be connected to the camera via 5G WiFi , supports multiple desktop computers, cell phones, tablets and laptops can realize microscopic view on the same screen.

, , Ugl bmuqNAqmdru_pe I mN_ A_aors pe Npm bau | jm_b jg i

<https://cn.ostec.com.cn/downloads/>;

Please refer to the “KoPa Capture Pro Software User Manual” for details.

Windows System Requirements

Microsoft Windows 10 (64 bit) or later version operating systems.

CPU: i5 8th generation or later version.

Hard drive: 512GB or more.

Memory: 16GB or more.

Graphics card: Core graphics.

Network card: 10/100/1000M adaptive.Wireless NIC supports 5G WiFi IEEE802.11ac.

, , Aml I cargnl k crf mb dmp acjj nf ml cq mpr_ jerq

Android System Requirements

Android System 7.0 or later version.


CPU dual-core 1.7Ghz or later version.

Memory RAM 3GB or more.
Storage ROM 32GB or more.
Support wireless protocol 5G WiFi IEEE802.11ac.

iOS System Requirements


IOS System 11.0 or later version.
CPU dual-core 1.8Ghz or later version.
Memory RAM 2GB or more.
Storage ROM 32GB or more.
Support wireless protocol 5G WiFi IEEE802.11ac.


1. Download and install the APP:

- 1)Click "KoPa WiFi Lab AO" software top left corner"  " settings icon, the bottom left corner will appear double QR codes.
- 2)The QR code on the left is used to download APP, WeChat code sweeping function does not support downloading APP, please use a browser or QR code of your cell phone or tablet to scan the QR to link to official website.
- 3)Official website link: <https://cn.ostec.com.cn/downloads/>.
- 4)Choose KoPa WiFi Lab for Android or KoPa WiFi Lab for iOS to download and install according to your phone and tablet system.
- 5)You can search and download KoPa WiFi Lab from Google Play app store.
- 6)You can search and download KoPa WiFi Lab from Apple App Store.

2. Scan QR code to live stream image automatically:

- 1)Run the built-in App KoPa WiFi Lab AO.
- 2)Click the QR code on the right to enlarge.

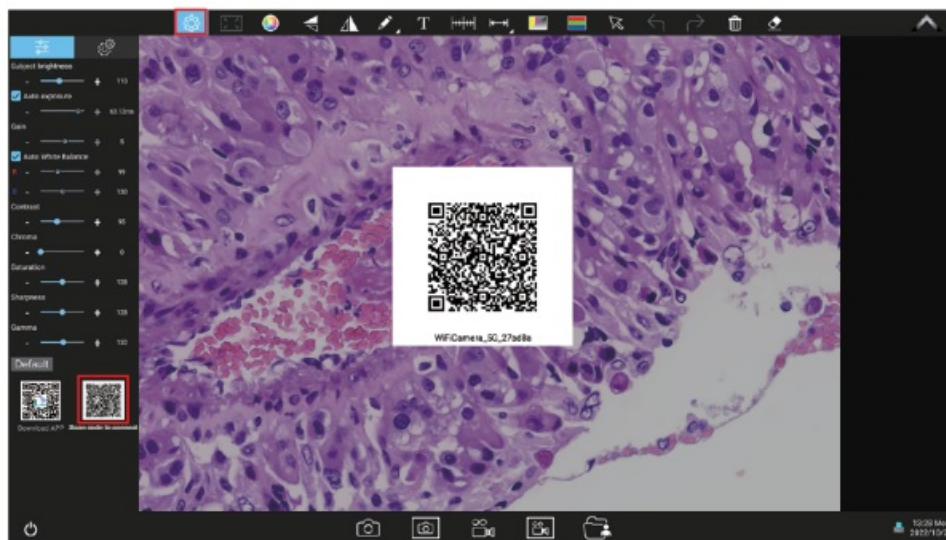


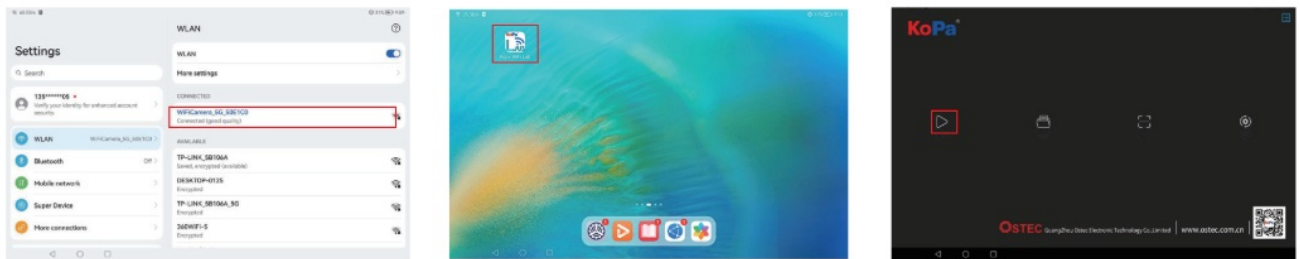
3. Click the "  " button on the navigation interface, scan the QR code on the camera body to connect to the microscope, no need to enter password, realize the same screen view.



(3)Manual connection method:

1)Search for the camera’s WiFi SSID: WiFi5Camera_xxxxxx in the WLAN of phone or tablet; default connection password: 12345678.




2)Run the App KoPa WiFi Lab, enter the App KoPa WiFi Lab navigation interface, and click  the button to connect to the microscope manually.



, , Jma_j_pe_l cru mpi ami 1 carginl

Once the camera connected to a local area network, it can be changed to a network camera. Multiple desktop computers, cell phones, tablets and laptops on the same local area network can realize synchronized real-time preview.

, , Rfck erf mb dmpantl | carrfca_k cp_rm_jma_j_pe_1 cru mpi

- (1) Click the “” setting icon in the upper left corner of KoPa WiFi Lab AO;
- (2) Click the icon “” ;
- (3) Click the icon “” at the bottom left, Open WLAN (as in Figure 1 and 2);
- (4) Select a local area network (Figure 3), enter password and click “Connect” button.

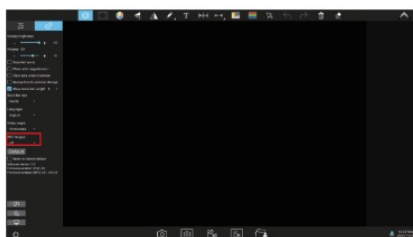


Figure 1

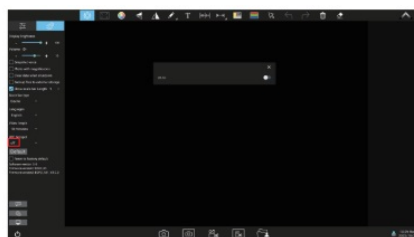


Figure 2

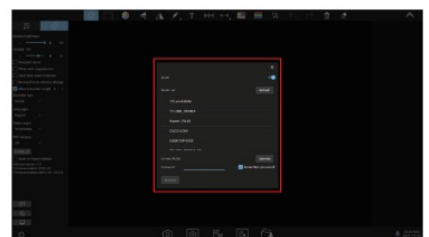



Figure 3

Notes:

1. The local area network must support 5G WiFi IEEE802.11ac protocol, otherwise it cannot be displayed in the list.
2. If the local area network opens the Internet function, the camera can also access the Internet.

, , Perpgete gk ecqtg rfc jma_j_pe_lcrumpi dpmk _Ug bmuqNAal rfeq_k c jma_j_pe_l cru mpi

1. Connect a Windows PC or laptop that supports 5G WiFi IEEE802.11ac protocol to the same local area network as the camera.
2. Install and run the software KoPa Capture Pro on the Windows PC or laptop.
3. When there is only one network camera in the local area network, the software can automatically get the live image from the camera.
4. When there are multiple network cameras in the same local area network, click the icon “” in the upper right corner of KoPa Capture Pro (Figure 1), the list automatically shows all cameras in the current local area network (Figure 2), select a camera as needed, support switching midway at any time.

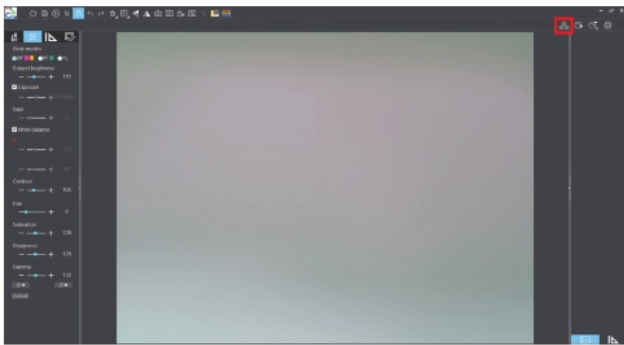


Figure 1

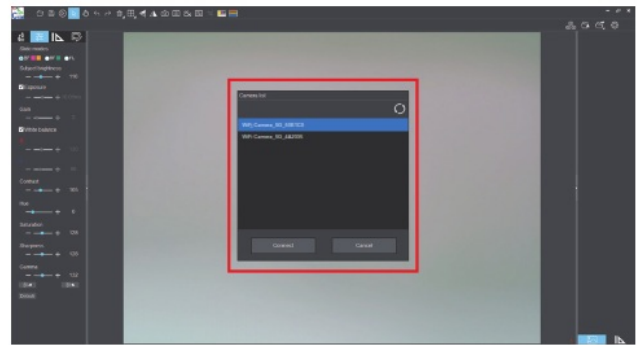




Figure 2

Notes:

When a Windows PC or laptop computer accesses the camera images, it does not affect the computer's ability to access the Internet, and if that local area network is open to the Internet, the Windows PC or laptop can also access the Internet.

, , Perpgete gk ecqtg rfc jma_j_pe_1l cru mpi dpmk acjjnf ml cq_1 br_— jerqgl rfc q_k c jma_j_pe_l cru mpi

1. Connect the cell phone or tablet that supports 5G WiFi IEEE802.11ac protocol to the same local area network as the camera.
2. Install and run the App KoPa WiFi Lab on the phone or tablet, enter the navigation interface and click the “” button.
3. When there is only one network camera in the local area network, the App can automatically get the live image from the camera.
4. When there are multiple network cameras in the same local area network, click the icon “” in the navigation interface (Figure 1), pop-up menu (Figure 2). The list automatically shows all cameras in the current local area network (Figure 3), select a camera as needed, support switching midway at any time.

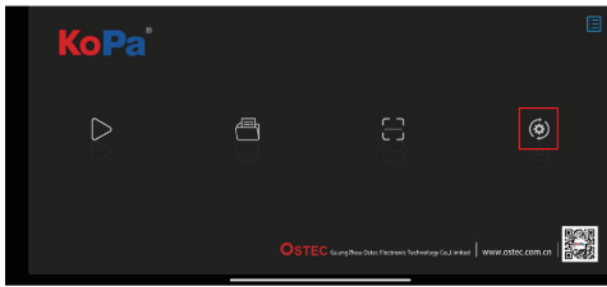


Figure 1



Figure 2

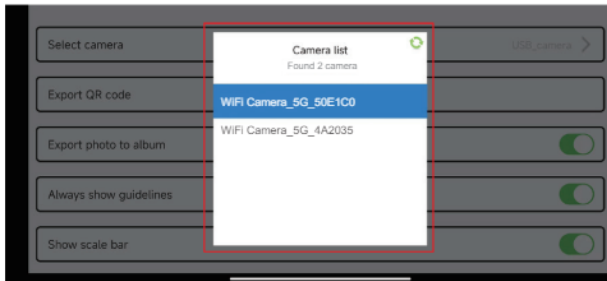


Figure 3

The Function Introduction And Connection Way For Video Set-top Box (Optional Accessory)

, N_a i g l e j g q r

1. Video set-top box (optional) Product model: NA001



2. HDMI cable (2m)



3. Wired mouse and keyboard




4. Power adapter and power cord (Optional Chinese, American, European, Australian, Korean, British standard, etc.)




, ?bt_lr_ecq_1b aml 1 cargml mdt gbcmqertrmn ' mvcq
 , , ?bt_lrecq

1. It is suitable for applications such as pathology seminars, demonstration teaching, and video sharing on the large display screens in research institutes.
2. It supports the whole series of KoPa smart embedded display camera, smart embedded camera and WiFi embedded cameras, and the factory has built-in high performance APP–KoPa WiFi Lab. which automatically connects with camera when it is turned on.
3. Built-in Android operating system, even if the large display screen has no operating system, it can wireless connect to the whole series of KoPa smart embedded display camera, smart embedded camera and WiFi embedded cameras.

, , Aml l c argml

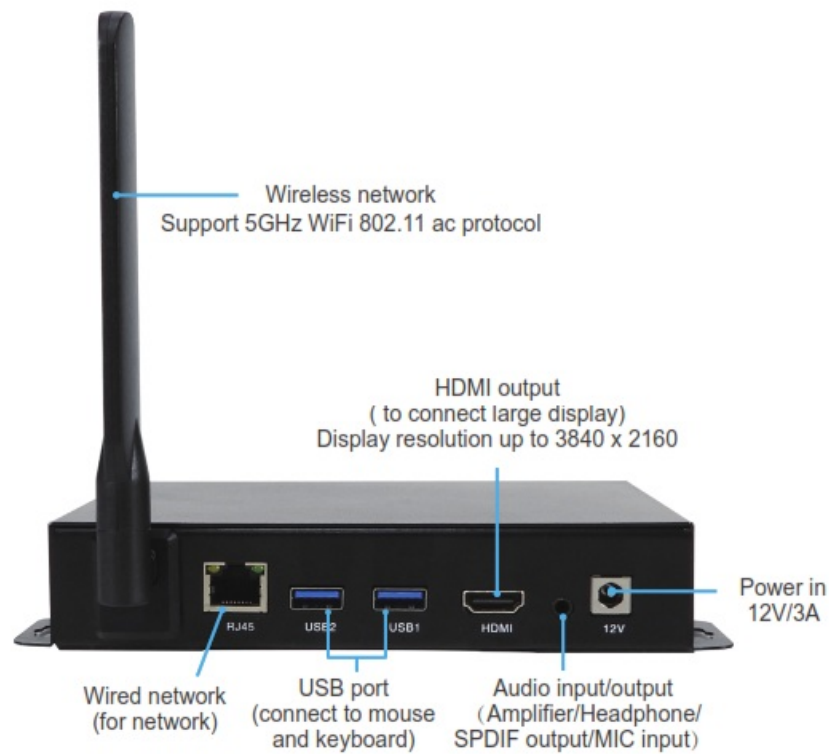

1. Connect the video set-top box to the oversized display with an HDMI cable, the display resolution supports up to 3840×2160.
2. The network cable is plugged into the video set-top box's wired network interface to give the video set-top box Internet access.
3. USB interface connects mouse and keyboard to facilitate the operation of the APP (KoPa WiFi Lab) on the oversized display.
4. Switching method for multiple network cameras in the same local area network:

Go to the APP (KoPa WiFi Lab) navigation interface, click the icon “” in the right (Figure 1), pop-up menu (Figure 2). The list automatically shows all cameras in the current local area network (Figure 3), select a camera as needed, support switching idway at any time.

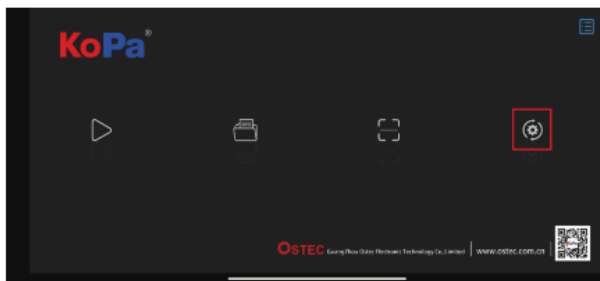


Figure 1

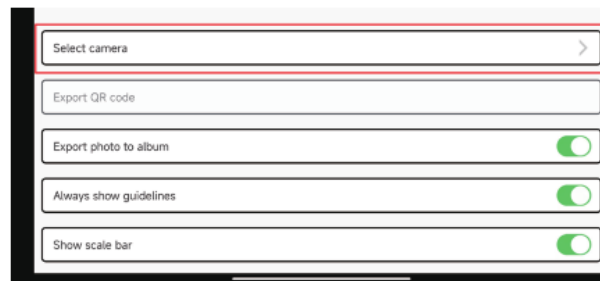


Figure 2

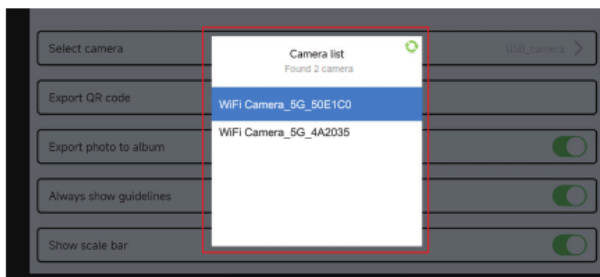


Figure 3

Troubleshooting And Frequently Asked Questions

1. Ensure that both ends of the power cord are plugged into the power socket and the power jack of the camera;

2. Ensure that the power on button has been long pressed for 24 seconds to turn on the camera. If the power indicator light still does not light up and the camera still cannot be turned on, please contact our customer service.

1. Ensure that both ends of the power cord are plugged into the power socket and the power jack of the camera;
2. Ensure that the power on button has been long pressed for 24 seconds to turn on the camera. If the power indicator light still does not light up and the camera still cannot be turned on, please contact our customer service.

1. Keep press the power switch for 4 seconds to turn off the camera, and keep press the power switch for 4 seconds to turn on the camera;

1. Keep press the power switch for 4 seconds to turn off the camera, and keep press the power switch for 4 seconds to turn on the camera;
2. If the power still cannot be turned on, please contact our customer service.

1. Keep press the power switch for 4 seconds to turn off the camera, and keep press the power switch for 4 seconds to turn on the camera;

1. Keep press the power switch for 4 seconds to turn off the camera, and keep press the power switch for 4 seconds to turn on the camera;
2. If the fault cannot be eliminated after restarting the camera, please contact our customer service

1. Ensure that the microscope has turned on the light and the image is visible to the eyepiece;

2. Keep press the power switch for 4 seconds to turn off the camera, and keep press the power switch for 4 seconds to turn on the camera;

- (1) Ensure that the microscope has turned on the light and the image is visible to the eyepiece;
- (2) Keep press the power switch for 4 seconds to turn off the camera, and keep press the power switch for 4 seconds to turn on the camera;
- (3) If the fault cannot be eliminated after restarting the camera, please contact our customer service.

, , rfc SQ@ qmai cr gq! mr u mpi gi e

(1) Ensure that the mouse and keyboard are working properly;

(2) Confirm that the USB socket is still not working after replacing mouse and keyboard, please contact our customer service.

, , Cvrepl _jF BK Gk ml grmp bmeq! mr bgqnj_w

(1) Ensure the HDMI cable can be used normally;

(2) Ensure that the external HDMI monitor can support 3840×2160 30Hz;

(3) If faulty cannot be eliminated, please contact our customer service.

, , Lmaqms! b dpmk fc_bnf aml c bai

(1) Ensure the headset currently in use can be used normally;

(2) Confirm that the headphone jack is still not working after replacing the headphones, please contact our customer service.

Disclaimer

1. In order to protect the legitimate rights and interests of users, please carefully read the instructions, disclaimers and safety instructions provided with this product before using it. The company reserves the right to update the above documents. Please operate the product according to the instructions and safety instructions.
2. Once you begin to use the product, you shall be deemed to have read, understood, recognized and accepted all terms and contents of the product's instructions, disclaimers and safety instructions. Users undertake to be responsible for their actions and all consequences. User undertakes to use the product only for legitimate purposes and agrees to these terms and any relevant policies or guidelines that the company may establish.
3. In the process of using this product, please strictly abide by and execute the requirements including but not limited to the instructions and safety instructions.

All personal injuries, accidents, property losses, legal disputes and other adverse events that cause conflicts of interest caused by violation of the safety instructions or irresistible factors shall be borne by users themselves, and the company shall not assume any responsibility.

4. Safety instructions:

- 1) Please do not use wet hands to plug and unplug the power supply of the equipment.
- 2) Please be sure to use a regular brand power socket, and make sure the grounding is well grounded to prevent electric shock.
- 3) Please be sure to make regular safety checks on sockets and plugs to avoid potential electrical safety hazards caused by aging and short circuit.
- 4) Please be sure not to use the product in a humid or hot environment to ensure the safety of the product.
- 5) Please be sure to loosen bundle cords of the power cable before using, to avoid electromagnetic induction and heating, thus increasing the heat dissipation speed.
- 6) Please be sure not to use the product equipment or socket or other places in high or easy to fall, so as to avoid damage.
- 7) Before opening the socket power supply, please be sure to turn off the load power switch of the product equipment.
- 8) The company reserves the rights to improve product, upgrade technology and change parameters without prior notice.

Documents / Resources



[KoPa HW1200 Binocular Stereo Microscope Camera Solution Smart Embedded Camera \[pdf\] User Manual](#)

HW1200 Binocular Stereo Microscope Camera Solution Smart Embedded Camera, HW1200, Binocular Stereo Microscope Camera Solution Smart Embedded Camera, Stereo Microscope Camera Solution Smart Embedded Camera, Microscope Camera Solution Smart Embedded Camera, Solution Smart Embedded Camera, Smart Embedded Camera, Embedded Camera

References

- [KoPa GUANGZHOU OSTEC ELECTRONIC TECHNOLOGY CO.,LTD](#)
- [KoPa GUANGZHOU OSTEC ELECTRONIC TECHNOLOGY CO.,LTD](#)
- [KoPa GUANGZHOU OSTEC ELECTRONIC TECHNOLOGY CO.,LTD](#)
- [User Manual](#)

[Manuals+](#), [Privacy Policy](#)

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.