

**LABOLYTIC
SMART 7 Inch
Colour Touch
Panel Advanced
Controller**



LABOLYTIC SMART 7 Inch Colour Touch Panel Advanced Controller Instructions

[Home](#) » [LABOLYTIC](#) » LABOLYTIC SMART 7 Inch Colour Touch Panel Advanced Controller Instructions 

Contents

- 1 LABOLYTIC SMART 7 Inch Colour Touch Panel Advanced Controller
- 2 Specifications
- 3 7" COLOUR TOUCH PANEL GUARANTEES INTUITIVE AND COMFORTABLE OPERATION
- 4 Advantages of SMART PRO Controller
- 5 FAQ
- 6 Documents / Resources
 - 6.1 References

LABOLYTIC

LABOLYTIC SMART 7 Inch Colour Touch Panel Advanced Controller




Specifications

- **Product Model:** SMART PRO
- **Controller Type:** Advanced 7 Colour Touch Panel
- Intuitive and Comfortable Operation
- **Temperature Sensor:** Pt 100 (option for extra sensor reading)
- **Control Parameters:** Air Flap, Fan, Humidity, CO2

Product Information

The SMART PRO is an advanced controller with a 7 colour touch panel that ensures intuitive and comfortable operation. It allows users to easily navigate through various functions and settings for precise control.

7" COLOUR TOUCH PANEL GUARANTEES INTUITIVE AND COMFORTABLE OPERATION

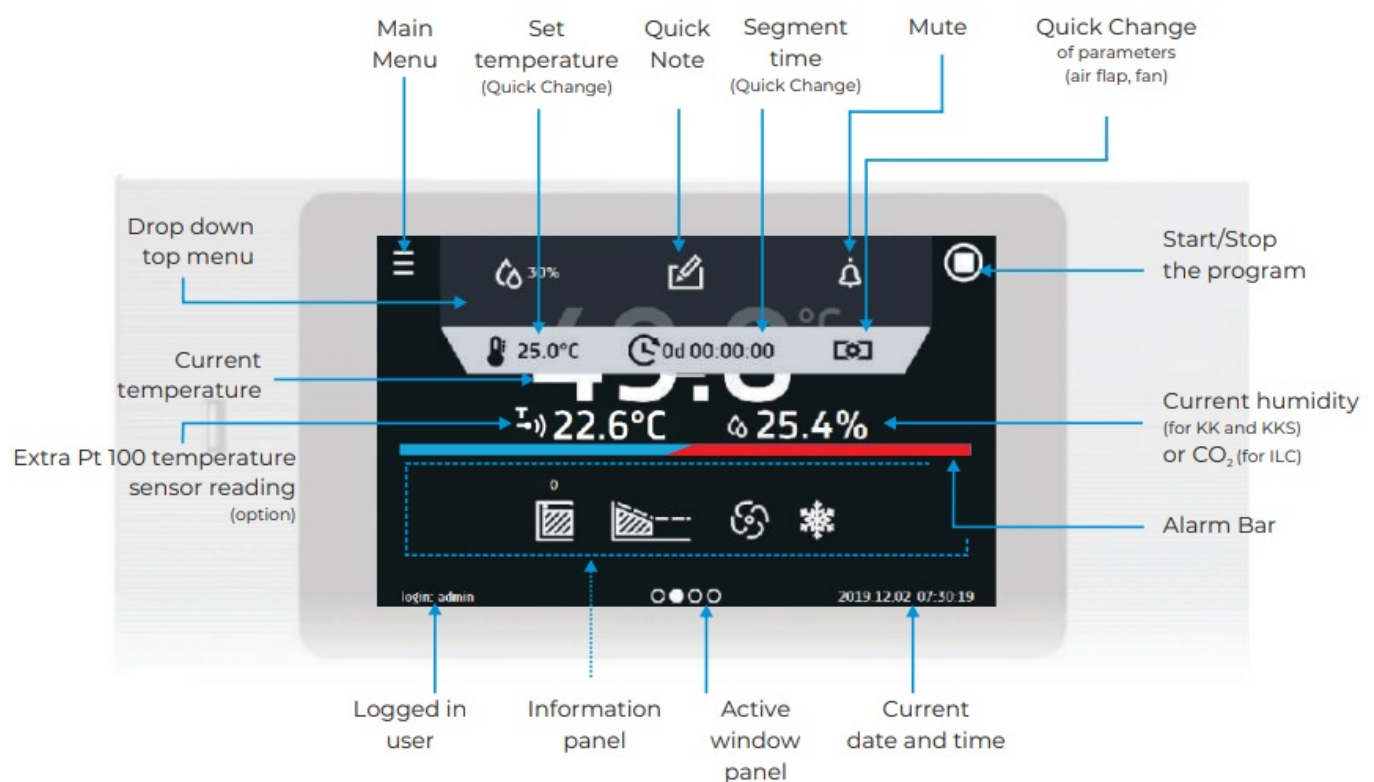

Usage Instructions

Main Menu:

- **Set Temperature:** Quickly change the set temperature.
- **Quick Segment Note Time:** Quickly change segment note time.
- **Mute:** Quick change settings like air flap and fan.
- **Drop Down Top Menu:** Access additional options.

Display Information:

- Current Temperature Displayed
- Extra Pt 100 Temperature Sensor Reading (if option installed)

- Current Humidity (for KK and KKS) or CO2 (for ILC)
- Alarm Bar for alerts
- Logged-in User Information
- Active Window Panel
- Current Date and Time

Advantages of SMART PRO Controller

- large (7"), clear, full-colour touch screen
- LAN, USB ports, Wi-Fi for communication and data transfer
- multi-segment time and temperature programs
- overview of data in tabular and graphic form
- visual and audible alarm
- administration functions for easy management
- password protected log-in
- internal memory for programs and data storage
- operating with gloves on
- event registry with user notifications

LabDesk software and instruction manual for direct download

- **Alarm Bar** – instant visual information about chamber status
- **Quick Note** – user can save text notes (50 characters) in Smart PRO controller memory
- **Quick Change of parameters:** temperature, humidity, time, air flap and fan (according to model)

FAQ


Q: How do I start/stop a program using the SMART PRO controller?

A: To start or stop a program, navigate to the main menu and select the option to start or stop the desired program.

Q: Can I adjust the set temperature quickly with the SMART PRO controller?

A: Yes, you can quickly change the set temperature by accessing the 'Set Temperature' option in the main menu.

Documents / Resources

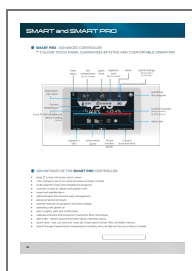

[LABOLYTIC SMART 7 Inch Colour Touch Panel Advanced Controller](#) [pdf] Instructions SMART, SMART 7 Inch Colour Touch Panel Advanced Controller, 7 Inch Colour Touch Panel Advanced Controller, Colour Touch Panel Advanced Controller, Touch Panel Advanced Controller, Advanced Controller, Controller

References

- [User Manual](#)

Manuals+. Privacy Policy

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.