

LIQUID INSTRUMENTS Moku:Go Frequency Response Analyzer User Manual

May 17,
2022

[Home](#) » [Liquid instruments](#) » **LIQUID INSTRUMENTS Moku:Go Frequency Response Analyzer User Manual** 

LIQUID INSTRUMENTS Moku:Go Frequency Response Analyzer User Manual

Moku:Go's Frequency Response Analyzer can be used to measure a system's frequency response from 10 mHz up to 30 MHz.

Frequency Response Analyzers are commonly used to measure the transfer functions of electrical, mechanical or optical systems by injecting a swept sinewave into the system and then comparing the output voltage to the input voltage. The resulting measurements of the system's magnitude and phase response can be used to optimise the closed-loop response of control systems, characterize resonant behaviour in non-linear systems, design filters, or measure the bandwidth of different electronic components. Frequency Response Analyzers are quite simply an indispensable tool in any electronics lab.

Contents [[hide](#)]

- 1 [User Interface](#)
- 2 [Main Menu](#)
- 3 [Export data](#)
- 4 [Signal Display Navigation](#)
- 5 [Settings](#)
- 6 [Normalization](#)
- 7 [Sweep modes](#)
- 8 [Cursors](#)
- 9 [Additional Tools](#)
- 10 [Power Supply](#)
- 11 [Documents / Resources](#)
 - 11.1 [References](#)
- 12 [Related Posts](#)

User Interface

User Interface

ID	Description	ID	Description
1	Main menu	6	Normalization*
2	Export data	7	Single/continuous mode switch*
3	Signal display navigation	8	Start / pause sweep*
4	Settings	9	Cursors

5	Control pane		
---	--------------	--	--

*Detailed information can be found in the Sweep mode section.

Main Menu

The main menu can be accessed by clicking the  icon on the top-left corner.

Main Menu

This menu provides the following options:

Options	Shortcuts	Description
Save/recall settings:		
Save instrument state	Ctrl+S	Save the current instrument settings.
Load instrument state	Ctrl+O	Load last saved instrument settings.
Show current state		Show the current instrument settings.
Reset instrument	Ctrl+R	Reset the instrument to its default state.
Power supply		Access power supply control window.*
File manager		Open file manager tool.
File converter		Open file converter tool.
Help		
Liquid Instruments website		Access Liquid Instruments website.
Shortcuts list	Ctrl+H	Show Moku:Go app shortcuts list.

Manual	F1	Access instrument manual.
Report an issue		Report bug to Liquid Instruments.
About		Show app version, check update, or license

Power supply is available on Moku:Go M1 and M2 models. Detailed Information about power supply can be found in Moku:Go power supply manual.

Export data

The export data options can be accessed by clicking the  icon, allowing you to:


Export data

Description

1. Select the type of data to export.
2. Select the file format (CSV or MAT).
3. Enter additional comments for the saved file.
4. Select the exporting location on your local computer.
5. Click to execute data export.
6. Click to close the export data window.


Signal Display Navigation




Signal display position

The displayed signal can be moved around the screen by clicking anywhere on the signal display window and dragging to the new position. The cursor will turn into a  icon once clicked, drag horizontally to shift along the frequency axis and drag vertically to shift along the amplitude/power axis.

The signal display can also be moved horizontally and vertically with arrow keys.

Display scale and zoom

Scrolling the mouse wheel zooms in and out along the primary axis. Access the scroll setting by hovering the cursor over the  icon.

Icons	Description
	Assign the horizontal axis as the primary axis.
	Assign the vertical axis as the primary axis
	Rubber band zoom: hold the primary mouse button to draw a region to zoom-in, release the button to execute.

Additional keyboard combinations are also available.

Actions	Description
Ctrl + Scroll Wheel	Zoom secondary axis
+/-	Zoom primary axis with keyboard
Ctrl +/-	Zoom secondary axis with keyboard.
Shift + Scroll Wheel	Zoom primary axis towards the center.


Ctrl + Shift + Scroll Wheel	Zoom secondary axis towards the center.
R	Rubber band zoom.


Auto scale

Double click anywhere on the signal display window to auto scale the traces.

Settings

The instrument control menu allows you to configure the Frequency Response Analyzer for your measurement, which will vary depending on the specific characteristics of the system under test.


Access the instrument Control menu by clicking the  icon.

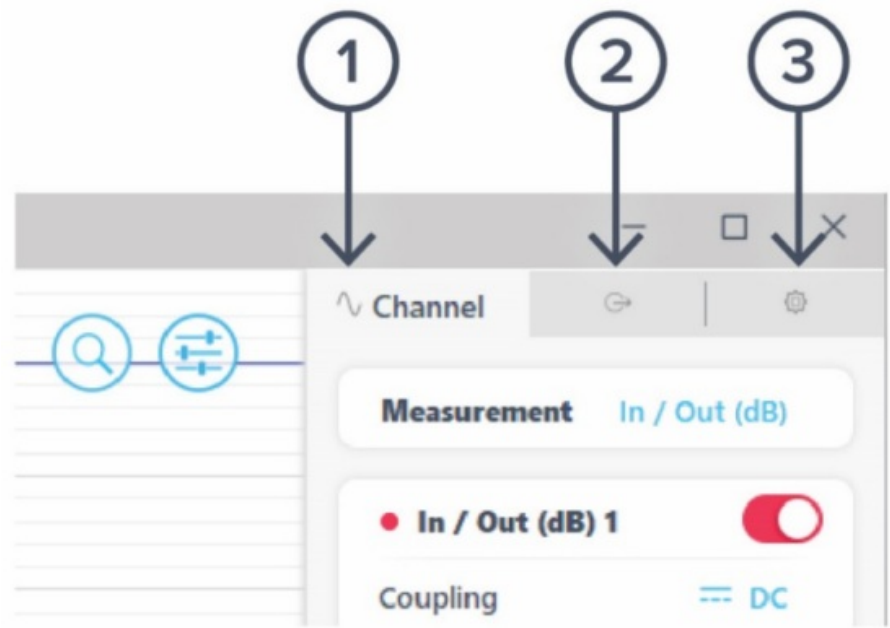

ID	Description
1	Channel

2	Swept sine
3	Advanced

Channels

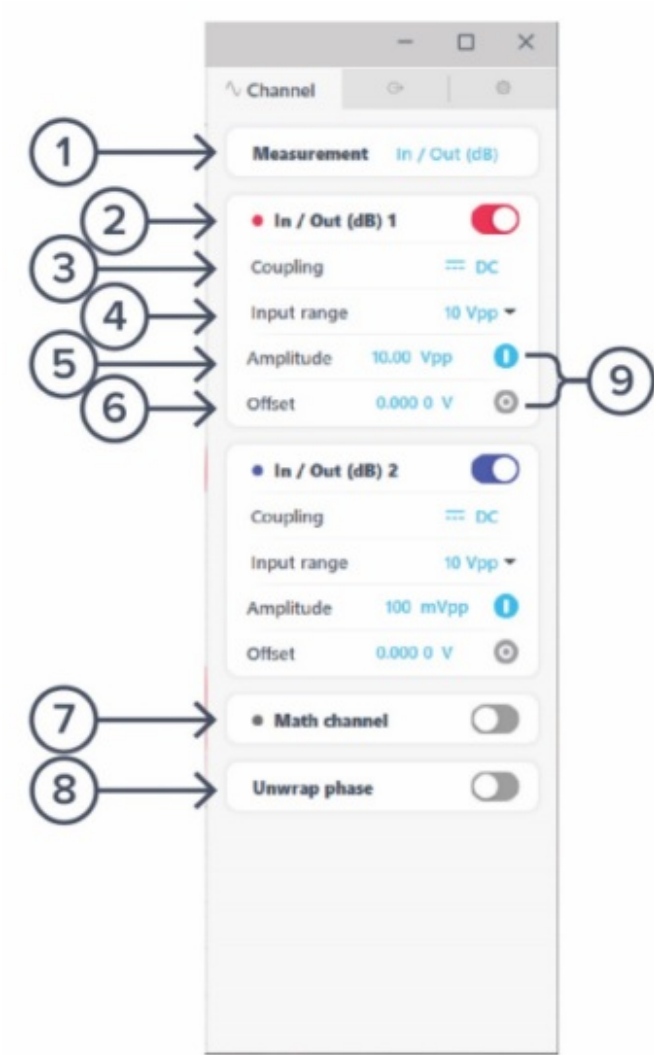

ID	Description	ID	Description
1	Select to display In (dBm) or In/Out (dB)	6	Swept sine (output) offset
2	Toggle channel on/off	7	Enable/disable Math channel

3	Select AC or DC coupling	8	Unwrap/wrap phase
4	Select input range 10 Vpp or 50 Vp p	9	Turn on/off amplitude and/or offset
5	Swept sine (output) amplitude		

Math channel

- Select between addition, subtraction, multiplication and division of the two channels.
- Compare transfer functions of channel 1 and 2 by configuring them identically.

Unwrap phase

- Phase is measured as a modulo of 2π . Enabling unwrapping will display an estimate of the total accumulated phase of the system.

Swept Sine



ID	Description	ID	Description
1	Configure the sweep start frequency	6	Configure the minimum averaging time
2	Configure the sweep stop frequency	7	Configure the minimum averaging cycles
3	Select the number of sweep point	8	Configure the minimum settling time
4	Select Linear or Log scale	9	Configure the minimum settling cycles
5	Reverse sweep direction	10	Total sweep time based on selected parameters

Sweep points

- Increasing the number of points in the sweep increases frequency resolution of the measurement allowing narrower features to be detected over a wider frequency range but will increase the total measurement duration.

Sweep scale

- The discrete points in the swept sine output can be spaced linearly or logarithmically. Logarithmic sweeps provide greater measurement resolution at lower frequencies.

Averaging

- Measurements at each point in the frequency sweep are averaged to improve accuracy and precision. You can configure the period over which each measurement is averaged to control the signal-to-noise ratio (SNR). Longer averaging times result in higher SNRs, allowing small features to be detected with greater precision. Shorter averaging times result in lower SNR measurements but the reduce total sweep time.
- The total averaging time is determined based on the minimum duration and minimum number of cycles over which each point in the sweep is averaged. Moku:Go's Frequency Reference Analyzer averages for the greater of the two values rounded upto the nearest number of integer cycles in order to avoid spectral leakage.

Settling time

- The settling time determines how long the Frequency Reference Analyzer waits before performing measurements at each frequency in the sweep. Settling time is important when characterizing resonant systems with high Q-factors in order to allow excitations to 'settle' between measurements. It can also be used to account for transmission delays in cables. When measuring a non-resonant system, the settling time should be set to equal the total propagation delay through the system.
- The total settling time is determined based on the minimum duration and minimum number of cycles over which the instrument will wait before beginning a measurement at each frequency in the sweep. The Frequency Response Analyzer


will wait for the greater effective duration of the two settings before beginning a measurement at each point in the sweep.

Advanced




ID	Description
1	Set the harmonic to demodulate for the frequency response
2	Set the phase difference between the output and local oscillator

Normalization


Moku:Go’s Frequency Reference Analyzer features a Normalization tool  that can be used to normalize subsequent measurements. Normalization is useful when

compensating for cable delays and comparing different devices under test.


Clicking the  icon will bring up the Normalization menu. Re-normalize will replace the current Normalization trace with a new one. Remove Normalization will erase all stored Normalization settings and cannot be undone.

Sweep modes


Single



Clicking the  icon will enable single sweep mode, which will pause the swept sine source at the end of the next full sweep. The swept sine signal will be turned off after the sweep completes and displayed data will not be updated.

Continuous


Clicking the  icon will enable continuous sweep mode, which will perform a new measurement as soon as the previous one has finished. This mode is commonly used to monitor systems with transfer functions that may change over time (e.g., control loops).

Pause / Restart


Clicking the  icon will immediately pause the current sweep. While paused, you can zoom in on features for more details, but no new data will be captured. Pressing the icon will also pause capture.

clicking the  or  icons will restart the sweep.

Cursors

Cursors can be accessed by clicking the  icon, allowing you to add power or frequency cursors, or remove all cursors. In addition, you can click and hold the cursors icon, and drag horizontally to add a frequency cursor, or drag vertically to add a magnitude or phase cursor.


User Interface

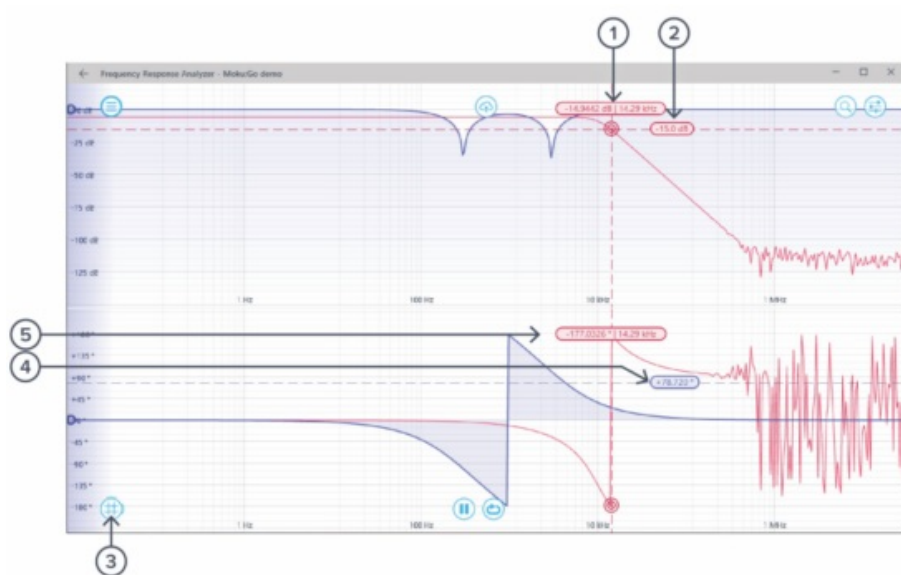

ID	Cursor item	Description
1	Frequency/Tracking cursor	Drag to reposition the cursor (Gray – Unattached, Red – channel 1, Blue – channel 2, Yellow – math).
2	Amplitude cursor	Drag to reposition, right click to set magnitude manually & other options.
3	Create cursor	Cursor options.
4	Phase cursor	Drag to adjust, right click to set phase manually & other options
5	Cursor label	Label depicting frequency, magnitude and phase of cursor. Drag to reposition.

Frequency Cursor


Right-click (secondary click) to reveal frequency cursor options:

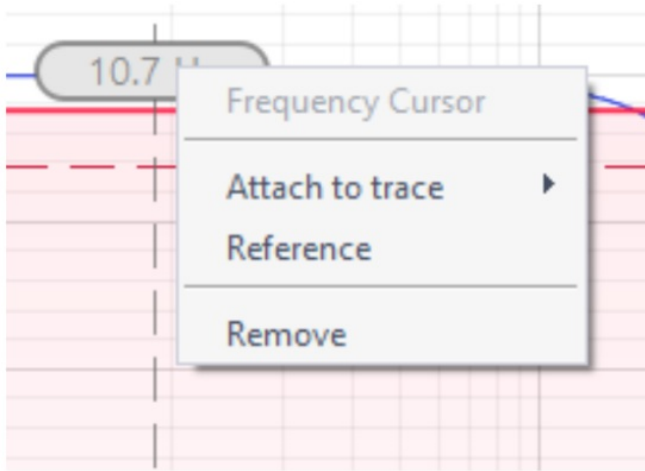

Options	Description
Frequency Cursor	Cursor type.
Attach to trace	Choose to attach the frequency cursor to channel A, channel B, or the math channel. Once the cursor is attached to a channel, it becomes a tracking cursor.
Reference	Set the cursor as the reference cursor. All other cursors in the same domain and channel measure the offset to the reference cursor.
Remove	Remove the frequency cursor

Tracking Cursor


Once a frequency cursor is attached to a channel, it becomes a tracking cursor. It displays the frequency and power level of the signal at the set frequency.

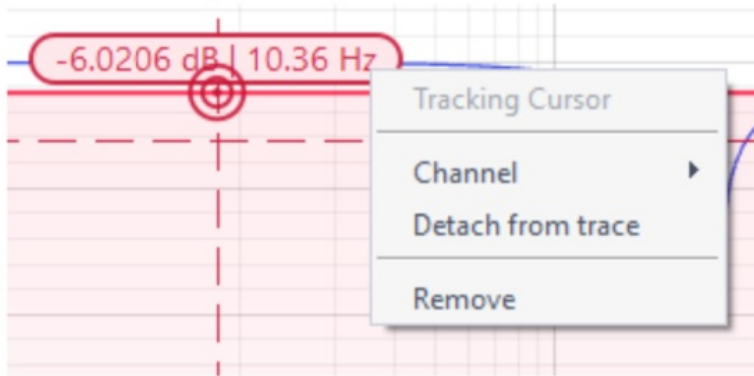

Options	Description
Tracking Cursor	Cursor type.
Channel	Assign tracking cursor to a specific channel
Detach from trace	Detach the tracking cursor from the channel to a frequency cursor.
Remove	Remove the tracking cursor

Magnitude/Phase Cursor

Right-click (secondary click) to reveal power cursor options:

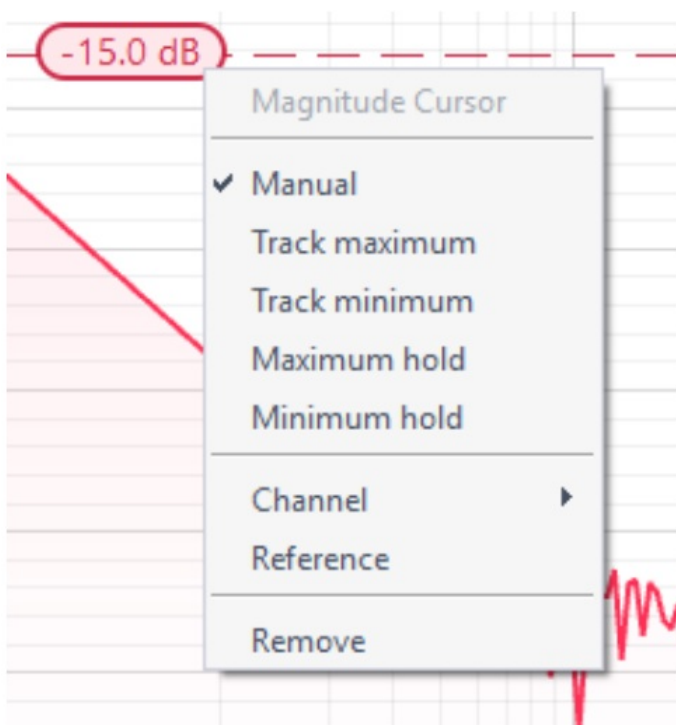


Options	Description
Manual	Manually set the vertical position of the cursor.
Track minimum	Track the maximum magnitude/phase.
Track minimum	Track the minimum magnitude/phase.
Maximum hold	Set the cursor to hold at the maximum magnitude/phase level.
Maximum hold	Set the cursor to hold at the minimum magnitude/phase level.
Channel	Assign the power cursor to a specific channel.
Reference	Set the cursor as the reference cursor.
Remove	Remove the magnitude/phase cursor.

Additional Tools

Moku:Go app has two built-in file management tools: file manager and file converter

File Manager


The file manager allows the user to download the saved data from Moku:Go to local computer, with optional file format conversion.

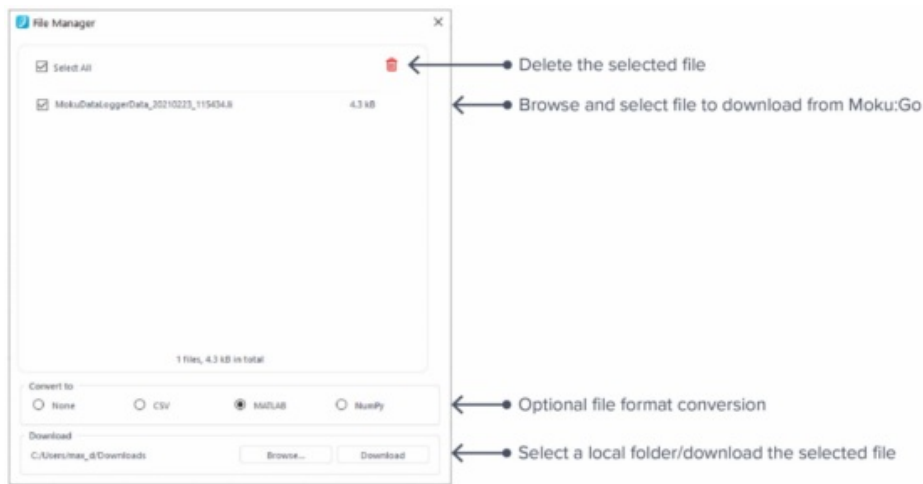

Once a file is transferred to the local computer, a  icon shows up next to the file.

File Converter

The file converter converts the Moku:Go's binary (.li) format on the local computer to either .csv, .mat, or .npz format.

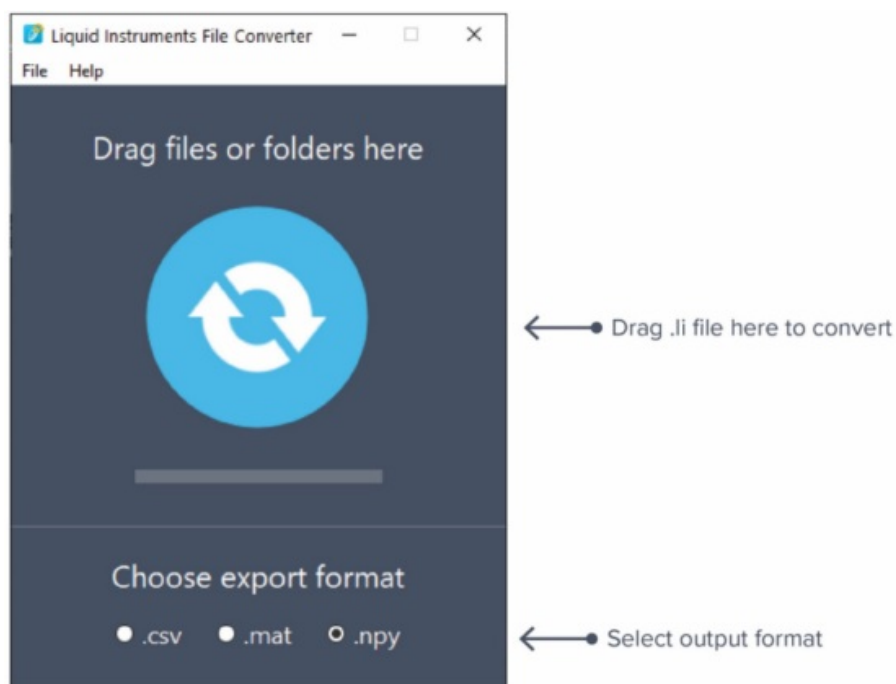

The converted file is saved in the same folder as the original file.

Liquid Instruments File Converter has the following menu options:


Options	Shortcut	Description
---------	----------	-------------

File			
·	Open file	Ctrl+O	Select a .li file to convert
·	Open folder	Ctrl+Shift +O	Select a folder to convert
·	Exit		Close the file converter window
Hel p			
·	Liquid Instruments w ebsite		Access Liquid Instruments website
·	Report an issue		Report bug to Liquid Instruments
·	About		Show app version, check update, or lice nse

Power Supply

Moku:Go Power supply is available on M1 and M2 models. M1 features a 2-channel power supply, while M2 features a 4-channel power supply. The power supply control window can be accessed in all instruments under the main menu.

The power supply operates in two modes: constant voltage (CV) or constant current (CC) mode. For each channel, the user can set a current and voltage limit for the output. Once a load is connected, the power supply operates either at the set current or set voltage, whichever comes first. If the power supply is voltage limited, it operates in the CV mode. If the power supply is current limited, it operates in the CC mode.

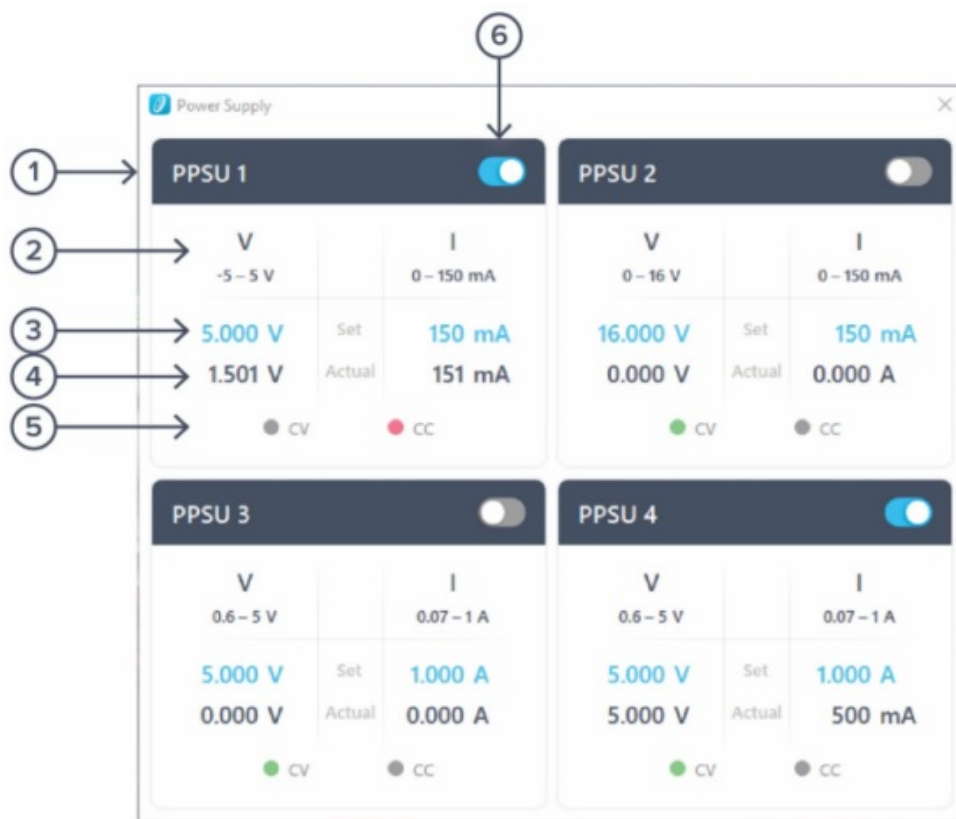


ID	Function	Description
1	Channel name	Identifies the power supply being controlled.
2	Channel range	Indicates the voltage/current range of the channel.
3	Set value	Click the blue numbers to set the voltage and current limit.
4	Readback numbers	Voltage and current readback from the power supply, the actual voltage and current being supplied to the external load.
5	Mode indicator	Indicates if the power supply is in CV (green) or CC (red) mode.
6	On/Off Toggle	Click to turn the power supply on and off.

Ensure Moku:Go is fully updated. For the latest information:

www.liquidinstruments.com

Documents / Resources

	<p>LIQUID INSTRUMENTS Moku:Go Frequency Response Analyzer [pdf]</p> <p>User Manual</p> <p>Moku Go, Frequency Response Analyzer</p>
---	--

References

- [Liquid Instruments](#)

Related Posts



[LIQUID INSTRUMENTS Moku:Go Portable Hardware Platform User Guide](#)

LIQUID INSTRUMENTS Moku:Go Portable Hardware Platform Moku:Go Layout Powering On & Off Connect the magnetic power supply to...



[Liquid instruments Moku Logic Analyzer /Patter Generator Liquid User Manual](#)

Liquid instruments Moku Logic Analyzer /Patter Generator Liquid User Interface Main menu Add channel Save data Output pattern...



[EHX Frequency Analyzer Ring Modulator User Guide](#)

EHX Frequency Analyzer Ring Modulator Congratulations! You have just purchased one of the most sophisticated and unusual musical...



[electro-harmonix FREQUENCY Analyzer Ring Modulator User Manual](#)

FREQUENCY ANALYZER RING MODULATOR

Congratulations! You have just purchased one of the most sophisticated and unusual musical accessories...

◆ Frequency Response Analyzer, Liquid instruments, Moku

■ Liquid instruments Go

—Previous Post

[LIQUID INSTRUMENTS Moku:Pro Device User Guide](#)

Next Post—

[LIQUID INSTRUMENTS Moku:Go Waveform Generator User Manual](#)

Leave a comment

Your email address will not be published. Required fields are marked *

Comment *

Name

Email

Website

☐ Save my name, email, and website in this browser for the next time I comment.

Post Comment



Search

Search

Manuals+,