

## **CLEANING SOLUTION 3**

for maintenance of a heliXcyto

Dynamic Biosensors GmbH & Inc.  
CS-50-3 v1.0



## Product Description

Order Number: **CS-50-3**

*Table 1. Contents and Storage Information*

Material	Amount	Storage
<b>Cleaning solution 3</b>	50 mL	RT

Handle with personal protection (gloves, goggles) and avoid direct skin contact. Rinse thoroughly with plenty of water in case of skin contact.

Store in a dark, cool, and well aired location. Avoid contact with acids.

For research use only.

This product has a limited shelf life, please see expiry date on label.

## Application Note

The Cleaning Solution 3 (CS 3) is ready-to-use for cleaning of a **helix<sup>cyto</sup>** device on a **maintenance chip**.  
**DO NOT** run over **helix<sup>cyto</sup>** trapping chips!

Follow the use of CS 3 immediately with RB 1 or another suitable buffer or plenty of H<sub>2</sub>O.

## Contact

**Dynamic Biosensors GmbH**  
Perchtinger Str. 8/10  
81379 Munich  
Germany

**Dynamic Biosensors, Inc.**  
300 Trade Center, Suite 1400  
Woburn, MA 01801  
USA

**Order Information**   [order@dynamic-biosensors.com](mailto:order@dynamic-biosensors.com)  
**Technical Support**   [support@dynamic-biosensors.com](mailto:support@dynamic-biosensors.com)

[www.dynamic-biosensors.com](http://www.dynamic-biosensors.com)

Instruments and chips are engineered and manufactured in Germany.  
©2024 Dynamic Biosensors GmbH | Dynamic Biosensors, Inc. All rights reserved.