

# USER MANUAL

Electronic Scale SF-400D-S



# **Specifications:**

Color: Black

Dimensions: (8.11 x 6.30 x 2.17) inch/(20.6 x 16.0 x 5.5) cm (L x W x H)

Weight: 17.4 oz/494 g

Material: Stainless Steel, Plastic

Maximum Capacity: 600 g Dividing Value: 0.01 g

Power supply: USB cord (to charge only) (Included), USB cord + AC/DC adapter (Included) or

2 x 1.5V AA battery (Not Included)
Weighing modes: g/oz/ct/dwt/T/gn/pcs

Package Dimensions: (8.07 x 9.45 x 2.95) inch/(20.5 x 24.0 x 7.5) cm (L x W x H)

Package Weight: 24.7 oz/700 g

## **Package Includes:**

1 x Electronic scale

1 x USB cord

1 x AC/DC adapter

1 x User manual

#### FIRST TIME USE

Remove the protective pads under the weighing pan and the protective film on it and on the display.

#### **NOTICE**

This is a precision instrument and must be handled with extreme care.

Avoid any exposure to extreme heat or cold. Scale works better when placed flat on the surface and started at normal room temperature. Always allow the scale to acclimate to normal room temperature for at least one hour before use.

Keep the scale clean. Dust, dirt, moisture, vibration, air currents, and a close proximity to other electronic devices can all cause an adverse effect on the reliability and accuracy of the scale. Handle with care. Gently apply all items to be weighed on the weighing platter. Although this scale is designed to be quite durable, try to avoid rough treatment as this may permanently damage the internal sensor. Avoid shaking, dropping or other shocking the scale.



# **Function Keys**

# Compact Electronic Scale





# Calibration

Long press CAL button. Press TARE button to switch between 200 g or 500 g calibration weight (not included). Press CAL button to confirm. Put the corresponding calibration weight. Display will show PASS – the calibration is finished.

#### On/Off

Switching on and off

#### Tare

Press one time to weigh to zero

### Units

Real-time convertible weighing units: Ounce (oz), Gram (g), Carat (ct), Pennyweight (dwt), Tael troy (T), Grain (gn)

# Pcs

Long press PSC button. Press MODE button to switch between 25/50/75/100 sample pool. Place the preset sample pool on the scale. Press PSC button to confirm. Place the unknown number of items and LCD will show the quantity calculated.

#### **Auto off function**

The auto off function only works when the scale is powered by batteries.

If it the scale is powered by USB or by using AC/DC adapter, the scale will not auto off. The auto off function can be configured or disabled by following the instructions below.

- 1. Start the scale and long press TARE key.
- 2. Then press Tare key to choose the mode you like.
  - 5 means auto power off in 5 minutes without any operation.
  - 3 means auto power off in 3 minutes.
  - 1 means auto off in 1 minute.
  - means the scale will not auto off.
- 3. Press MODE key to save setting.