# PLW 8683 Safety

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|        |        |
|--------|--------|
| Breite | 670 mm |
|--------|--------|

|       |        |
|-------|--------|
| Tiefe | 740 mm |
|-------|--------|

|                                                   |           |
|---------------------------------------------------|-----------|
| <b>Information on the installation plan .....</b> | <b>20</b> |
| Definition of terms .....                         | 20        |
| <b>Appliance dimensions.....</b>                  | <b>21</b> |
| Freestanding machines .....                       | 21        |
| Built-under machines .....                        | 21        |
| <b>Connections .....</b>                          | <b>22</b> |
| <b>Installation notes .....</b>                   | <b>26</b> |
| Environmental requirements.....                   | 26        |
| Vapour barrier for built-under machines .....     | 26        |
| Electrical connection .....                       | 26        |
| Equipotential bonding .....                       | 26        |
| Electromagnetic compatibility (EMC) .....         | 27        |
| Network connection .....                          | 28        |
| Water connection .....                            | 29        |
| Water inlet .....                                 | 29        |
| Drainage .....                                    | 30        |
| <b>Technical data .....</b>                       | <b>31</b> |
| Electrical connection .....                       | 31        |
| Water inlet .....                                 | 32        |
| Drainage .....                                    | 32        |
| Dimensions .....                                  | 33        |
| Weights .....                                     | 33        |
| Emission values .....                             | 34        |
| Heat dissipation .....                            | 34        |
| Sound pressure level .....                        | 34        |
| Operating conditions .....                        | 34        |
| Storage and transportation conditions .....       | 34        |

## en - Information on the installation plan

---

For safe installation and commissioning of the washer-disinfector please read the installation plan, the service documentation and the operating instructions.

This installation plan includes the dimensions of the washer-disinfector, the technical data and the installation requirements to be met on the customer's site.

The machine's equipment will vary by version and can be expanded with additional options. The installation plan describes the maximum equipment configuration.  
Certain connections are not required for every equipment version.

### **Installation requirements**

This machine must only be installed and commissioned by the Miele Customer Service Department or a suitably authorised technician.

Installation should only be performed in accordance with valid regulations, relevant standards and health and safety codes.

### **Definition of terms**


**Cleaning machine** In this installation plan, the washer-disinfector is referred to as "the machine".

**Wash water** The term "wash water" refers to water or to a mixture of water and process chemicals.


**Process chemicals** All media dispensed during a programme sequence are generally referred to as process chemicals, e.g. cleaning agents.

## Freestanding machines

Front view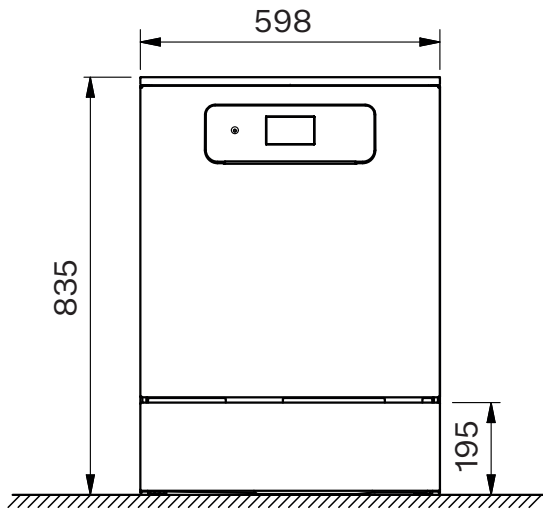


Side view

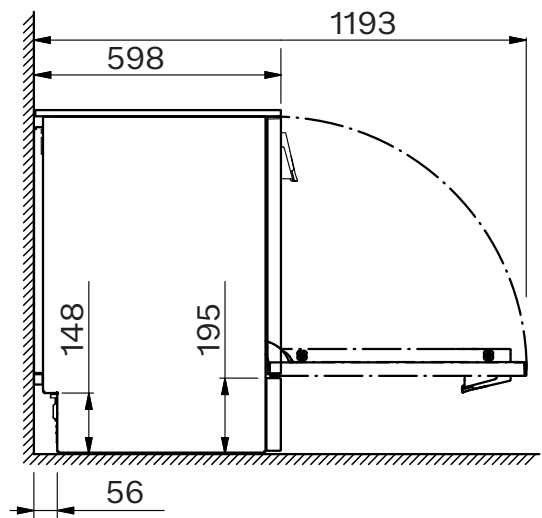
## Built-under machines


Front view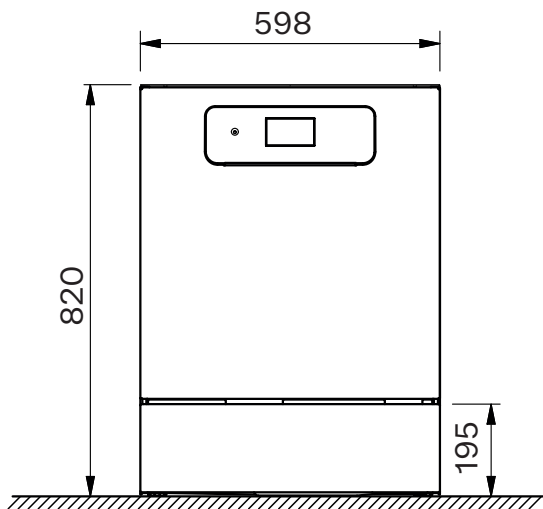

Side view

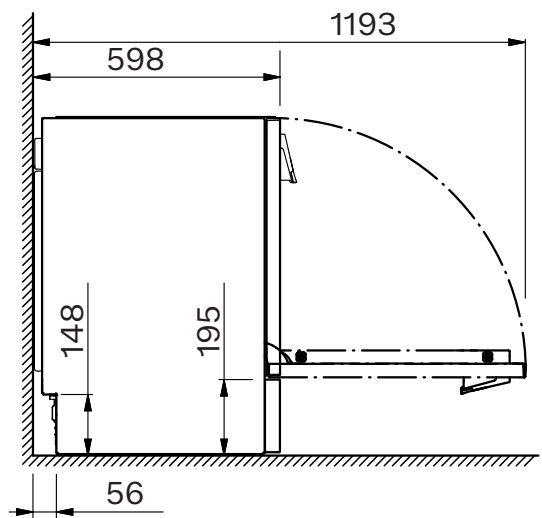
### Rear view

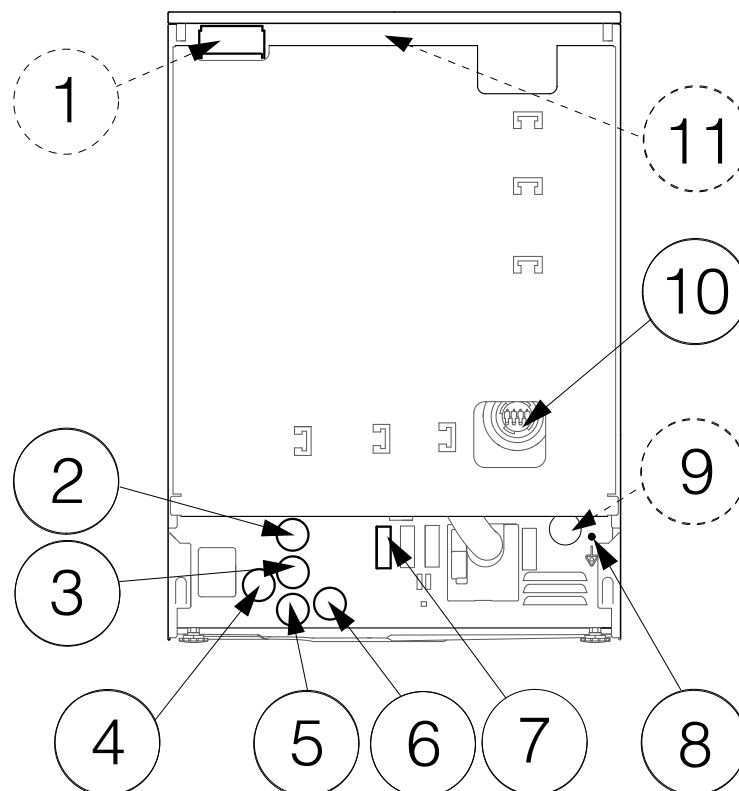
- ① Network connection (optional)
- ② Mains connection cable
- ③ Hot water
- ④ Demineralised water
- ⑤ Cold water
- ⑥ Waste water
- ⑦ External dispensing, power supply connection
- ⑧ Equipotential bonding
- ⑨ Water inlet with feed pump (optional)
- ⑩ External dispensing, connection for dispensing hose
- ⑪ Potential-free contact (optional)


## On-site connections

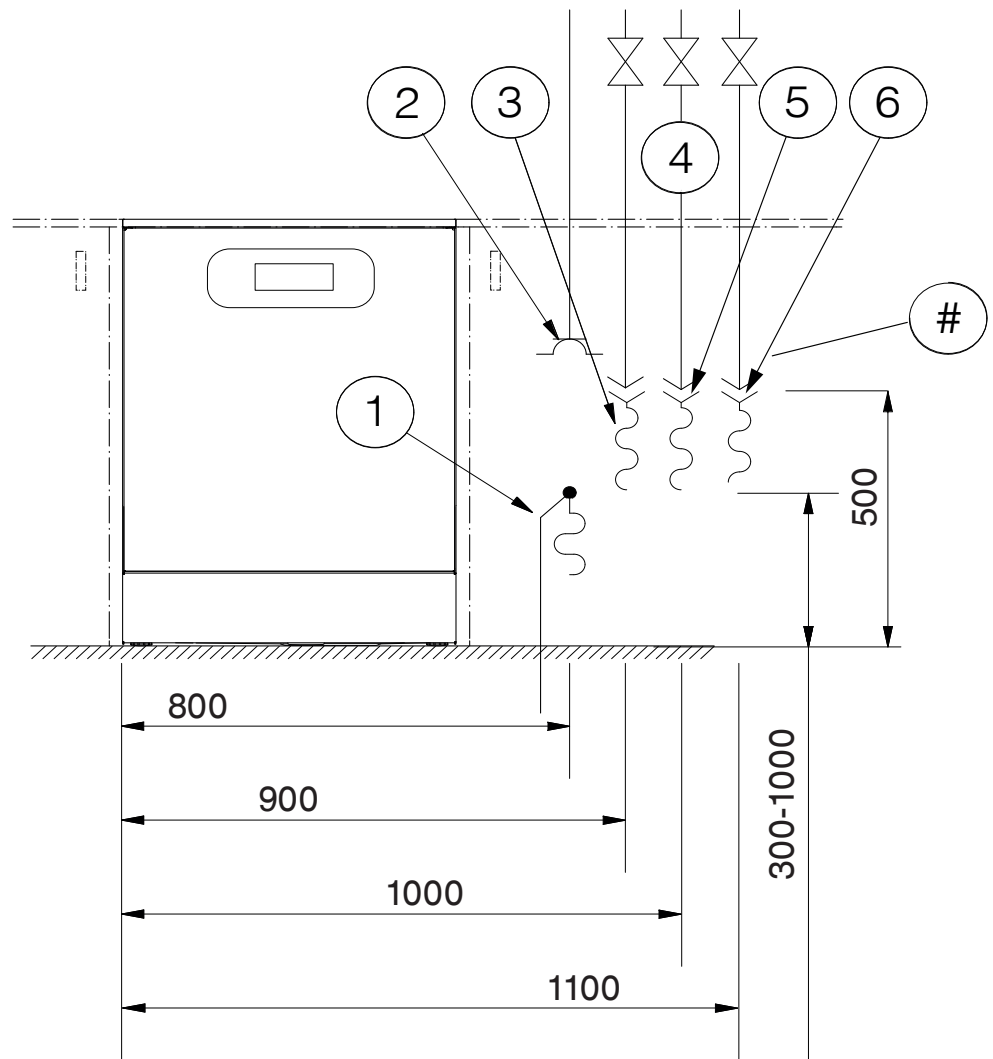- ① Waste water
- ② Mains connection cable
- ③ Demineralised water
- ④ Stopcocks
- ⑤ Cold water
- ⑥ Hot water

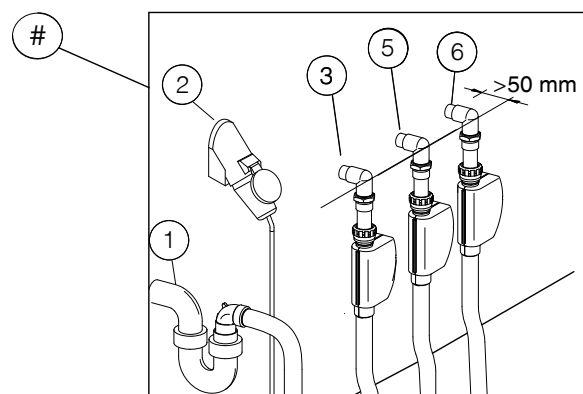
### External dispensing
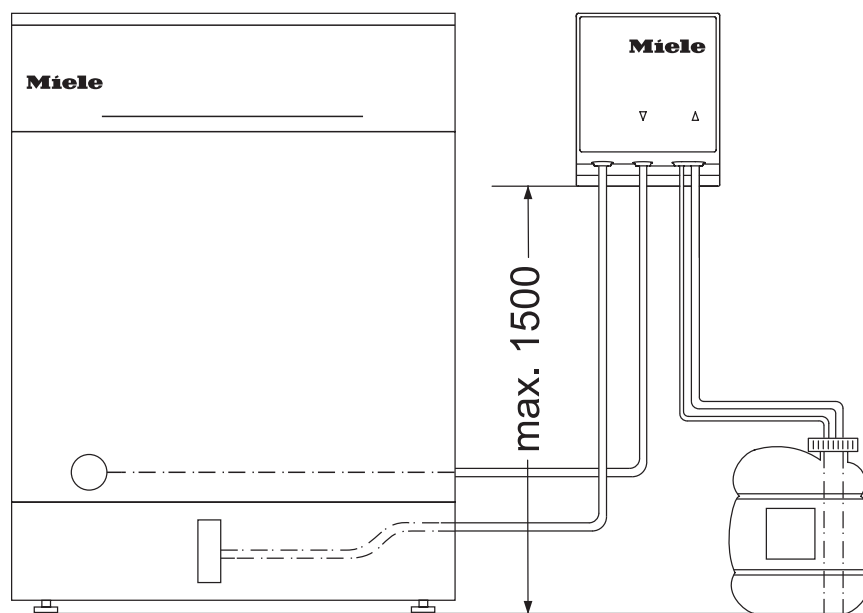
|                                                                 |       |
|-----------------------------------------------------------------|-------|
| Max. delivery head                                              | 1.5 m |
| Length of dispensing hose, DOS module to suction lance          | 1.8 m |
| Length of dispensing hose, back of machine to DOS module        | 2.8 m |
| Length of mains connection cable, back of machine to DOS module | 2.8 m |

Place the container next to the machine on the floor or in an adjacent cabinet.  
The container must not be placed on the machine or above the machine.

External feed  
pump (depending  
on country)

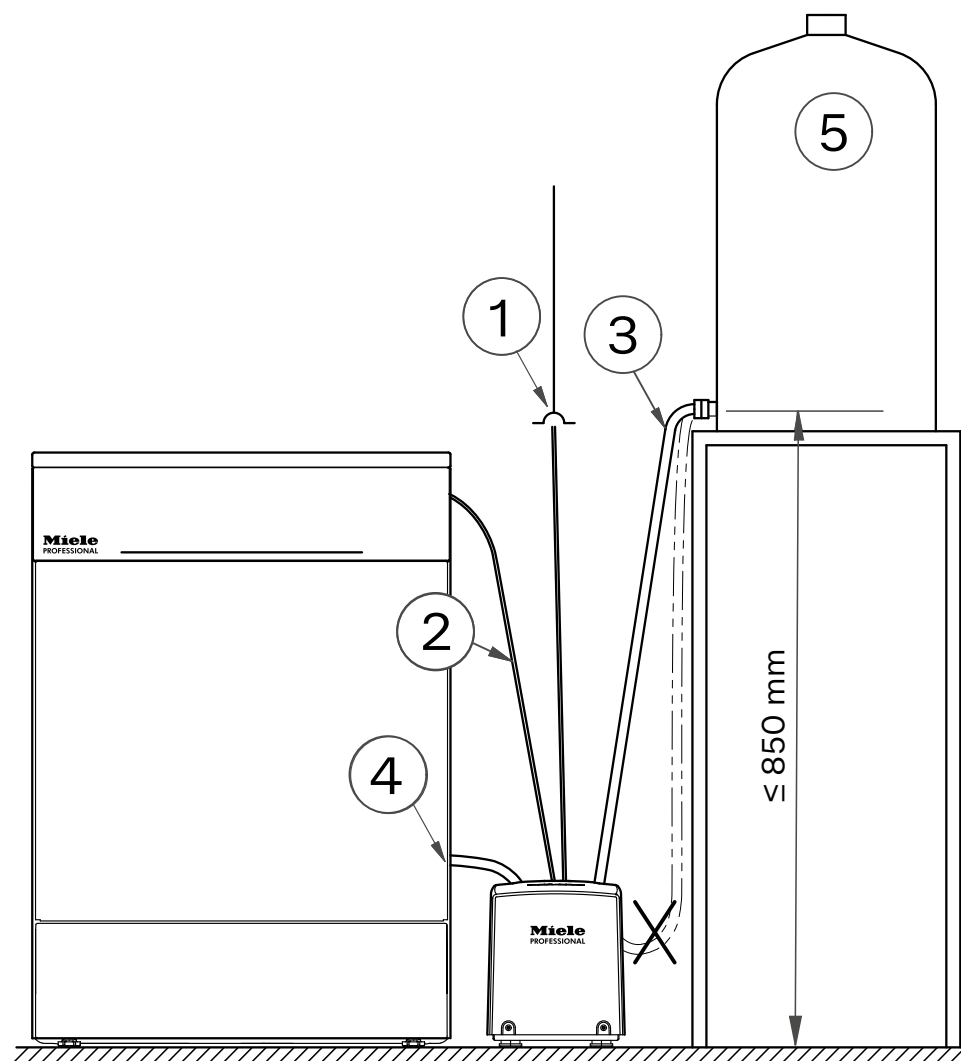|   |                                                            |       |
|---|------------------------------------------------------------|-------|
| ① | Mains connection cable length                              | 2.5 m |
| ② | Length of mains connection cable to machine controls       | 2.5 m |
| ③ | Length of water inlet hose from water container to module  | 1.5 m |
| ④ | Length of water inlet hose from module to cleaning machine | 1.3 m |
| ⑤ | Water container                                            |       |

Place the water container next to the machine on a cabinet or pedestal.  
The water container must not be placed on the machine.

### Environmental requirements

Condensation can build up in the area surrounding the machine. Any furniture and fittings in the room must therefore be suitable for purpose.

#### Vapour barrier for built-under machines

The vapour barrier supplied protects the worktop from damage caused by steam when the door is opened. Fit the vapour barrier above the door under the worktop.

### Electrical connection

All work on the electrical connection must be carried out by Miele Customer Service or a qualified electrician.

The customer's installation of the machine must comply with over-voltage category CAT II in accordance with IEC 61010-1.

The mains voltage must not fluctuate by more than  $\pm 10\%$  around the nominal voltage.

#### Main switch

Install a main switch on the customer's site that is capable of disconnecting the machine at all poles. The main switch must:

- Have a contact gap of at least 3 mm
- Be designed to operate at the rated current of the machine
- Be able to be locked in the zero position
- Be accessible after the machine has been installed

The mains connection cable must be laid protected from the risk of thermal damage.

#### Plug connection

Connect the cleaning device to the mains power supply using a suitable socket that complies with national regulations.

#### Hard-wired

Alternatively, the machine can be connected via a hard-wired connection.

#### Residual current device (RCD)

For added safety, every machine should be protected by an RCD type A with a trip current of 30 mA.

The RCD must be installed so that it is easily accessible.

#### Equipotential bonding

The washer is able to accommodate equipotential bonding. The connection screw (M 8) is located on the rear of the machine.

Equipotential bonding should be carried out if possible on the customer's site.

### **Electromagnetic compatibility (EMC)**

The machine has been tested for electromagnetic compatibility (EMC, class B) in accordance with EN 61326-1 and is suitable for operation in commercial environments, such as hospitals, medical practices and laboratories and other similar environments which are connected to the public supply network.

The machine's high frequency (HF) energy emissions are very low and are therefore unlikely to interfere with other electronic machines in the vicinity.

Flooring at the installation site must be wood, concrete or tiled. Synthetic flooring must be able withstand a relative humidity level of at least 30 % to minimise the risk of electrostatic discharges.

The quality of the supply voltage should comply with that found in a typical commercial or hospital environment. Check that the power supply voltage is within a range of +/-10 % of its nominal value.

### Network connection

The machine can be integrated into a customer's network via an Ethernet interface.

The operator is responsible for the secure configuration and operation of the network. Creating a secure configuration reduces the scope for the machine to be attacked. Non-secure configurations can make it easier for potential attackers to gain unauthorised access.

 **Unauthorised access poses a risk.**  
Settings in the machine may be changed as a result of unauthorised access via the network.  
Under no circumstances should it be possible to access the machine via public or unsecured networks, either directly or indirectly (e.g. using port forwarding).  
The IP addresses from 192.168.10.1 to 192.168.10.255 are reserved for Customer Service. They must never be set in the internal network.

Please also note the following recommendations for the network configuration of the machine:

1. Configure the network in which the machine is located with as many restrictions as possible:
  - Only allow machines or persons access to the network if absolutely necessary
  - Use a VLAN for network segmentation, for example
2. Make sure to create a secure configuration for all machines connected to the network, e.g. with the help of:
  - The information on network security in the operating instructions for the connected machines
  - The recommendations of the German Federal Office for Information Security (<https://www.bsi.bund.de>) for Europe and relevant cyber security centres in your country.
3. Pay particular attention to the following areas:
  - User management
  - Staff entering and exiting
  - Authorisation
  - Authentication
  - Update management, etc.

## Water connection

### Water inlet

The machine must only be connected to fully vented pipework. A brief increase in the water pressure can damage components of the machine.

The quality of the incoming water must at least correspond to the drinking water specification of the country in which the machine is being operated.

The machine must be connected to cold and hot water. If hot water is not available, the inlet hose for hot water must be connected to the cold water supply. For this purpose, a Y-piece can be used on the cold water inlet on the customer side.

To be able to use the programmes supplied by the factory, a DI water connection is also required.

The connections with stopcocks at the customer's site must have 3/4" flat-sealing external threads.

The stopcocks must be accessible after the machine has been installed.

### DI water ring line

The machine can be connected to a ring line system for DI water. For this purpose, a retrofitting kit must be installed in the machine by Miele Customer Service or an authorised technician and the control system must be set accordingly.

### DI water inlet for 8,5-60 kPa (optional)

For the connection to 8,5-60 kPa overpressure, the machine must be retrofitted with a feed pump by Miele Customer Service or an authorised technician.

The outlet nozzle of the unpressurised demineralised water container must be at least level with the top edge of the machine.

### Drainage

The machine should be connected to a separate on-site drainage system for the washer only. If a separate connection is not available, we recommend connecting the hose to a dual-chamber siphon. The waste water pipes must be temperature-resistant to at least 93 °C. If the hose is to be fitted directly to the drainage system on site, use the supplied hose clip.

The on-site connector for the drain hose can be adapted to different hose diameters. If the connector extends more than 30 mm into the drain hose, it must be shortened. Otherwise, the drain hose can become blocked.

Lay the drain hose so that it does not kink and is not being subjected to pressure or tension.

If the on-site drain connection is situated lower than the guide path for the lower basket rollers in the open door, a siphoning effect during a programme can cause the wash cabinet to empty itself of water. In this case, lay the drain hose with a bend in it so that its highest point is at least level with the guide path for the lower basket rollers.

## Electrical connection

|                                      | Default                 | Voltage variant *       |
|--------------------------------------|-------------------------|-------------------------|
| Voltage                              | 3N AC 400 V             | AC 230 V                |
| Frequency                            | 50 Hz                   | 50 Hz                   |
| Fuse rating                          | 3 x 16 A                | 16 A                    |
| Plug                                 | CEE **                  | Typ F                   |
| Mains connection cable length        | 1.9 m                   | 1.9 m                   |
| Mains connection cable cross-section | 5 x 2.5 mm <sup>2</sup> | 3 x 1.5 mm <sup>2</sup> |
| Heater rating                        | 8.5 kW                  | 2.5 kW                  |
| Total rated load                     | 9.3 kW                  | 3.3 kW                  |

## Versions for BE, NO

|                                      | Default                 | Voltage variant *       |
|--------------------------------------|-------------------------|-------------------------|
| Voltage                              | 3 AC 230 V              | AC 230 V                |
| Frequency                            | 50 Hz                   | 50 Hz                   |
| Fuse rating                          | 3 x 20 A                | 16 A                    |
| Plug                                 | -                       | Typ F                   |
| Mains connection cable length        | 1.9 m                   | 1.9 m                   |
| Mains connection cable cross-section | 4 x 2.5 mm <sup>2</sup> | 3 x 1.5 mm <sup>2</sup> |
| Heater rating                        | 6.7 kW                  | 2.5 kW                  |
| Total rated load                     | 7.5 kW                  | 3.3 kW                  |

\* Conversion work done by the Miele Customer Service Department

\*\* CEE plug only for AT, BE, CH, CZ, DE, DK, ES, FI, FR, GR, IE, IT, LU, NL, NO, PL, PT, SE, SK, TR

## en - Technical data

---

### Water inlet

|                                                                                  |               |
|----------------------------------------------------------------------------------|---------------|
| Water connections                                                                |               |
| - Cold water                                                                     | 1             |
| - Hot water                                                                      | 1             |
| - Demineralised water                                                            | 1             |
| Water temperature                                                                |               |
| - Cold water                                                                     | 5–20 °C       |
| - Hot water                                                                      | 5–65 °C       |
| - Demineralised water                                                            | 5–20 °C       |
| Connection hose length                                                           | 1.7 m         |
| Connection at the customer's site with stopcock,<br>flat-sealing external thread | 3/4"          |
| Water connection pressure                                                        | 200–1000 kPa  |
| Flow rate                                                                        | 7.5 l/min     |
| Water hardness                                                                   | 0–10.7 mmol/l |
| Chloride content                                                                 | <100 mg/l     |
| Demineralised water                                                              |               |
| - Recommended conductivity                                                       | <15 µS/cm     |
| - pH value                                                                       | 5–8           |

### Drainage

|                               |          |
|-------------------------------|----------|
| Drain hose length             | 1.4 m    |
| Max. drain hose length        | 4 m      |
| Max. delivery head from floor | 1 m      |
| Max. water temperature        | 93 °C    |
| Max. transient flow rate      | 16 l/min |
| Hose inner diameter (Ø)       | 22 mm    |
| Waste water pipe              | ≥ DN 50  |

## Dimensions

|                                |         |
|--------------------------------|---------|
| Height of freestanding machine | 835 mm  |
| Height of built-under machine  | 820 mm  |
| Height adjustment              | +60 mm  |
| Width                          | 598 mm  |
| Depth                          | 598 mm  |
| Depth including control panel  | 639 mm  |
| Depth with door open           | 1193 mm |

## Weights

|                                                    |                        |
|----------------------------------------------------|------------------------|
| <b>Freestanding machine</b>                        |                        |
| Net weight                                         | 74 kg                  |
| Max. operating weight                              | 130,5 kg               |
| Maximum load capacity of open door                 | 37 kg                  |
| Maximum load capacity of machine lid               | 69 kg                  |
| Max. individual load (Ø 38 mm)                     | 330 N                  |
| Max. surface load (load-bearing capacity of floor) | 3,65 kN/m <sup>2</sup> |
| <b>Built-under machine</b>                         |                        |
| Net weight                                         | 72 kg                  |
| Max. operating weight                              | 128,5 kg               |
| Maximum load capacity of open door                 | 37 kg                  |
| Max. individual load (Ø 38 mm)                     | 320 N                  |
| Max. surface load (load-bearing capacity of floor) | 3,59 kN/m <sup>2</sup> |

## en - Technical data

---

### Emission values

#### Heat dissipation

|                           |          |
|---------------------------|----------|
| Heat output into the room | 5.4 MJ/h |
|---------------------------|----------|

#### Sound pressure level

|                                                     |                     |
|-----------------------------------------------------|---------------------|
| Sound pressure level LpA during cleaning and drying | <65 dB(A) re 20 µPa |
|-----------------------------------------------------|---------------------|

### Operating conditions

|                     |         |
|---------------------|---------|
| Ambient temperature | 5–40 °C |
|---------------------|---------|

Relative humidity:

|         |      |
|---------|------|
| Minimum | 10 % |
|---------|------|

|                                       |      |
|---------------------------------------|------|
| Maximum, for temperatures up to 31 °C | 80 % |
|---------------------------------------|------|

|                                                |      |
|------------------------------------------------|------|
| Linear decreasing for temperatures up to 40 °C | 50 % |
|------------------------------------------------|------|

|                               |          |
|-------------------------------|----------|
| Max. altitude above sea level | 2000 m * |
|-------------------------------|----------|

\* If installed above 1500 m, the boiling point of the wash water will be lower. In this case, Customer Service must adjust the disinfection temperature and the holding time.

### Storage and transportation conditions

|                     |               |
|---------------------|---------------|
| Ambient temperature | -20 to +60 °C |
|---------------------|---------------|

|                   |         |
|-------------------|---------|
| Relative humidity | 10–85 % |
|-------------------|---------|

|              |              |
|--------------|--------------|
| Air pressure | 500–1060 hPa |
|--------------|--------------|

|                            |       |
|----------------------------|-------|
| Weight including packaging | 79 kg |
|----------------------------|-------|

|                                       |        |
|---------------------------------------|--------|
| Maximum floor load on transport route | 790 kN |
|---------------------------------------|--------|

### Site access dimensions including transport pallet

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