



NeuroNexus IST-nDrive User Manual

[Home](#) » [NeuroNexus](#) » NeuroNexus IST-nDrive User Manual 




IST-nDrive User Manual

IST-nDrive

This user manual will show you how to prepare your probe and nDrive for implantation.


Step 1

Prepare the nDrive, nDrive base, and IST-nDrive. Do not attach the probe or probe mount yet.

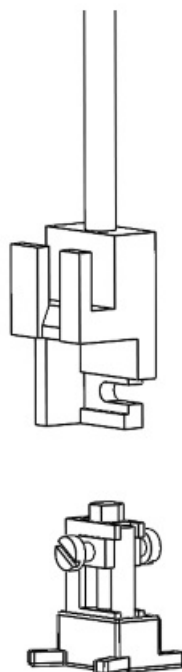


Step 2

Loosen the back screw on the nDrive. Slide the nDrive assembly onto the IST-nDrive.

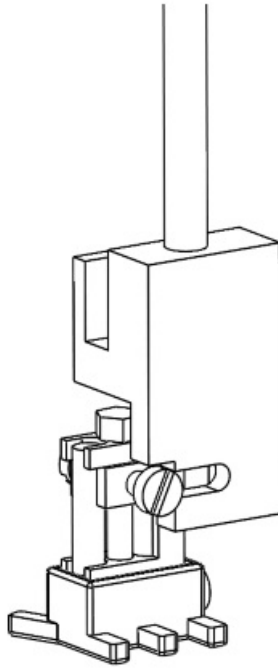


Step 3

Tighten the screw to lock the nDrive onto the IST-nDrive.

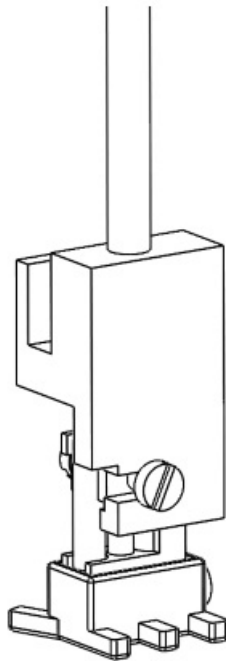


Step 4

Align the probe mount with the front screw.

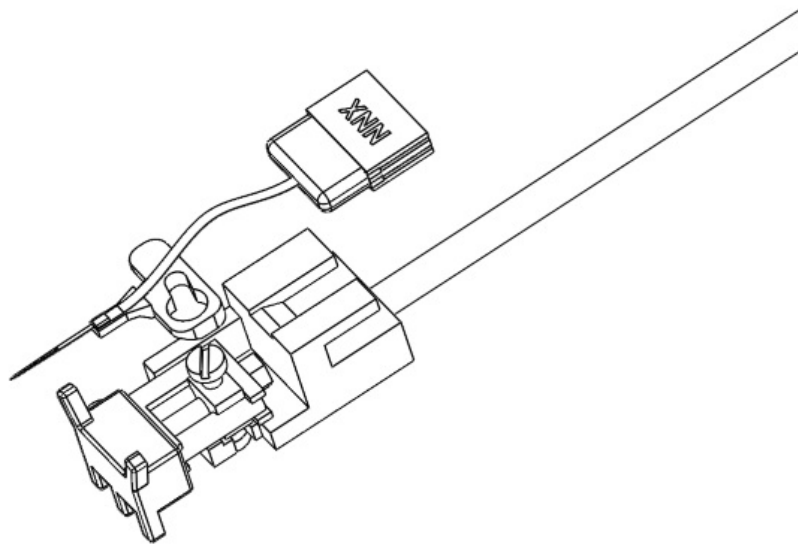


Step 5

Push the front screw through the larger probe mount screwhead opening.

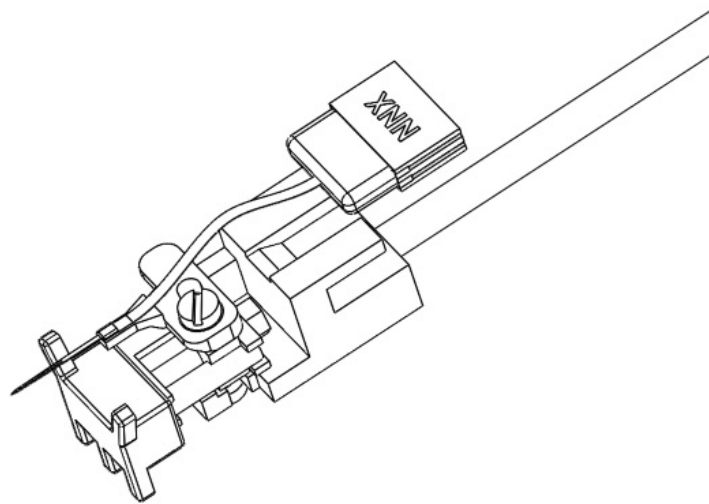


Step 6

Slide the probe mount to the right and tighten the screw to lock it in place.

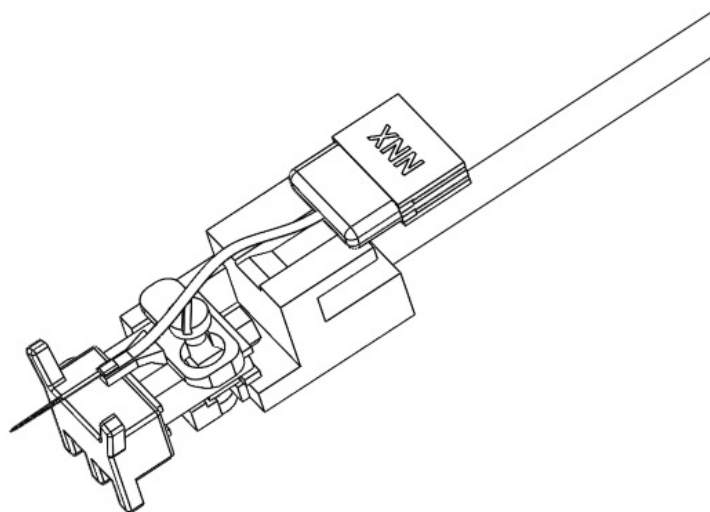

Step 7

Place the probe connector into the connector bracket.

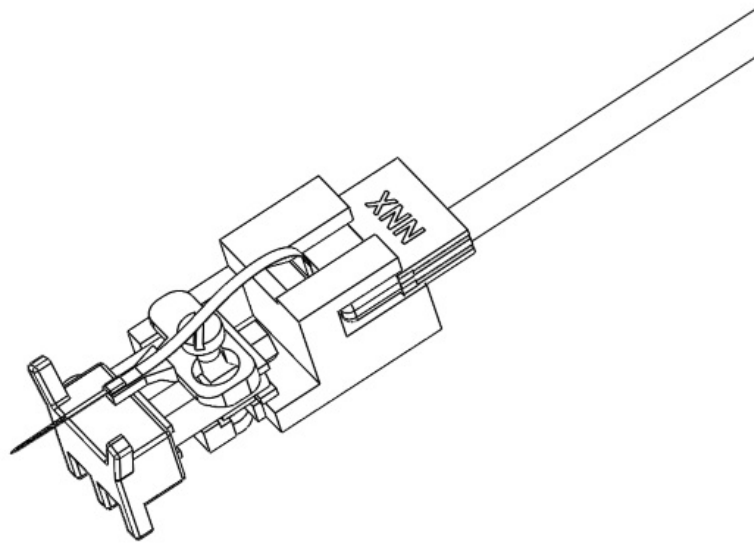


Step 8

Mount the IST drive on a stereotaxic frame.


Perform insertion at target location. Cement the drive base once initial insertion has been performed.

NOTE: Apply cement to only the nDrive base, leaving room to access the screw to release the nDrive unit after the experiment.

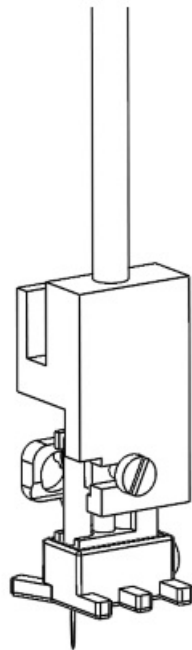

Step 9

Once the nDrive base has been cemented, loosen the back screw and carefully slide the IST-nDrive away.

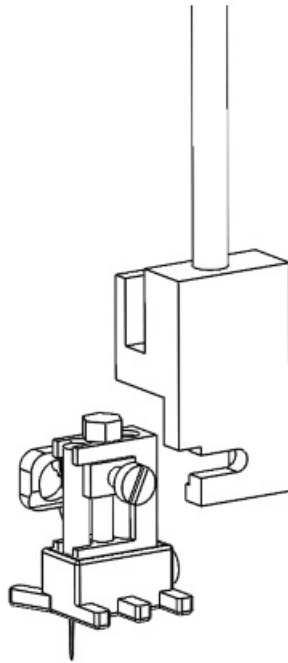


NeuroNexus Technologies, Inc. ©2021 | 5200 S. State Street, Suite 200, Ann Arbor, Michigan USA
Telephone: +1.734.913.8858 | Fax: +1.734.786.0069 | support@neuronexus.com

Contents

- [1 Documents / Resources](#)
- [2 Related Posts](#)

Documents / Resources

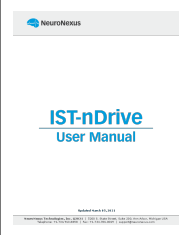	NeuroNexus IST-nDrive [pdf] User Manual IST-nDrive
---	---