



nuwave LumAir LA100 Pathogen Detection System User Guide

[Home](#) » [nuwave](#) » nuwave LumAir LA100 Pathogen Detection System User Guide 



LumAir LA100 Pathogen Detection System Cartridge Loading/Unloading Guide

This guide is intended to provide instructions on how to insert or remove a cartridge from an installed LA100 Pathogen Detection System.



Note: Please see the cartridge data sheet for detailed information on the application, use, and disposal of

Contents

- [1 Introduction](#)
- [2 Package Contents](#)
- [3 Step 1: Inserting a New Cartridge \(visual steps\)](#)
- [4 Step 2 – Removing the Cartridge](#)
- [5 Documents / Resources](#)
- [6 Related Posts](#)

Introduction


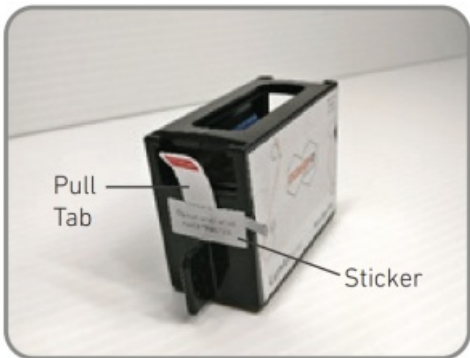

This guide is intended to provide instructions on how to insert or remove a cartridge from an installed LumAir LA100 Pathogen Detection System.



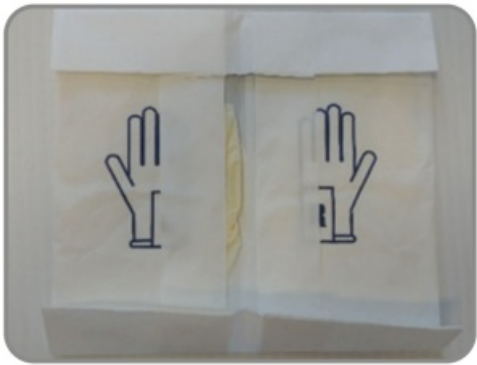




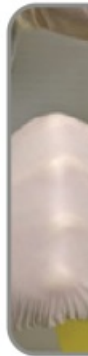
Package Contents

	
Cartridge (in Foil Bag)	
	
Bio-Hazard Safety Disposal Bag	

Step 1: Inserting a New Cartridge (visual steps)

1. Put on sterile gloves using sterile technique	2. Remove Cartridge from f
	
3. Peel back sticker to remove cartridge pull tab	4. Once the pull tab is remo
	
5. Insert cartridge fully into the unit	6. LED status changes to g
	

Step 2 – Removing the Cartridge


1. Put on sterile gloves using sterile technique	2. Push button, turning the LED status li
	
3. Peel back the sticker fully	4. Insert the locking slide (included in ca
	
5. Ensure locking slide is completely inserted	6. Remove cartridge and place in yellow
	

For questions not covered in this guide:

Email: support@nuwavesensors.com

Tel: (+353) 1 254 4188 (International)

Documents / Resources

	nuwave LumAir LA100 Pathogen Detection System [pdf] User Guide nuwave, LumAir, LA100, Pathogen, Detection, System, Cartridge Loading, Unloading Guide
---	--