



LabCon Map PC Recording 4 Channel Mapping PC Software User Manual

[Home](#) » [PPI](#) » LabCon Map PC Recording 4 Channel Mapping PC Software User Manual 

Contents

- [1 LabCon Map PC Recording 4 Channel Mapping PC Software](#)
- [2 OPERATOR PAGE PARAMETERS](#)
- [3 SUPERVISORY](#)
- [4 FACTORY](#)
- [5 FRONT PANEL KEYS](#)
- [6 PV ERROR INDICATIONS](#)
- [7 ELECTRICAL CONNECTIONS](#)
- [8 Documents / Resources](#)
 - [8.1 References](#)
- [9 Related Posts](#)



LabCon Map PC Recording 4 Channel Mapping PC Software



This brief manual is primarily meant for quick reference to wiring connections and parameter searching. For more details on operation and application; please log on to www.ppiindia.net

OPERATOR PAGE PARAMETERS

Parameters	Settings (Default Value)
Time Start Command >>	Yes No
Time Abort Command >>	(Default : No)
Time Interval (H:M) >>	0.00 to 500.00 (HH:MM) (Default : 0.10)
Ctrl Set Value >>	Setpoint LO limit to Setpoint HI limit (<i>Resolution 0.1°C for RTD / DC Linear & 1°C for Thermocouple</i>) (Default : 25.0)
Ctrl Lo Deviation >>	For RTD & DC Linear : 0.2 to 99.9 For Thermocouple : 2 to 99 (Default : 2.0)
Ctrl Hi Deviation >>	For RTD & DC Linear : 0.2 to 99.9 For Thermocouple : 2 to 99 (Default : 2.0)
Change Password >>	1 to 100 (Default : 0)

SUPERVISORY

SUPERVISORY > SENSOR INPUT

Parameters	Settings (Default Value)
Ctrl Zero Offset >>	-50 to 50 <i>(Resolution 0.1°C for RTD / DC Linear & 1°C for Thermocouple)</i> (Default : 0.0)
Map-1 Zero Offset >> Map-2 Zero Offset >> Map-3 Zero Offset >> Map-4 Zero Offset >>	-50 to 50 <i>(Resolution 0.1°C for RTD / DC Linear & 1°C for Thermocouple)</i> (Default : 0.0)

SUPERVISORY > RECORDING

Parameters	Settings (Default Value)
Rec. Interval (min) >>	0 to 250 Min (Default : 5 Min)
Alarm Toggle Record >>	Enable Disable (Default : Enable)
Del PC Records >>	Yes No (Default : No)
Del Display Record >>	Yes No (Default : No)
GSM Machine ID >> (refer note)	1 to 128 (Default : 1)
RTC Time >>	0 to 23 for Hour 0 to 59 for Min (Default : NA)
RTC Date >>	1 to 31 (Default : NA)
RTC Month >>	1 to 12 (Default : NA)
RTC Year >>	2000 to 2099 (Default : NA)
Device Id >>	1 to 127 (Default : 1)

SUPERVISORY > CONTROL

Parameters	Settings (Default Value)
Tune >>	Yes No (Default : No)
Setpoint LO Limit >>	Min Range for the Selected Input Type to Setpoint HI Limit (<i>Resolution 0.1°C for RTD / DC Linear & 1°C for Thermocouple</i>) (Default : 0.0)
Setpoint	Setpoint LO Limit to
HI Limit >>	Max Range for the Selected Input Type
	(<i>Resolution 0.1°C for RTD /</i>
	<i>DC Linear & 1°C for Thermocouple</i>)
	(Default : 600.0)
Compressor	0 to 100
Setpoint	(<i>Resolution 0.1°C for RTD / DC Linear & 1°C for Thermocouple</i>)
	(Default : 45.0)
Compressor Hyst >>	0.1 to 99.9 (Default : 2.0)
Heat Ctrl Action >>	ON-OFF PID (Default : PID)
Heat Hyst >>	0.1 to 99.9 (Default : 0.2)

Heat Only Control	Heat + Cool Control Zone : Single	Heat + Cool Control Zone : Dual
Proportional Band >> 0.1 to 999.9 (Default : 50.0)	Proportional Band >> 0.1 to 999.9 (Default : 50.0)	Cz Prop Band >> <i>Proportional Band for Cool Pre-dominant zone</i> 0.1 to 999.9 (Default : 50.0)

Integral Time >> 0 to 3600 sec (Default : 100 sec)	Integral Time >> 0 to 3600 sec (Default : 100 sec)	Cz Integral Time >> <i>Integral Time for Cool Pre-dominant zone</i> 0 to 3600 sec (Default : 100 sec)
Derivative Time >> 0 to 600 sec (Default : 16 sec)	Derivative Time >> 0 to 600 sec (Default : 16 sec)	Cz Derivative Time >> <i>Derivative Time for Cool Pre-dominant zone</i> 0 to 600 sec (Default : 16 sec)
Cycle Time >> 0.5 to 100.0 sec (Default : 10.0 sec)	Cycle Time >> 0.5 to 100.0 sec (Default : 10.0 sec)	H_z Prop Band >> <i>Proportional Band for Heat Pre-dominant zone</i> 0.1 to 999.9 (Default : 50.0)
Overshoot Inhibit >> Enable Disable (Default : Disable)	Overshoot Inhibit >> Enable Disable (Default : Disable)	H_z Integral Time >> <i>Integral Time for Heat Pre-dominant zone</i> 0 to 3600 sec (Default : 100 sec)
Cutoff Factor >> 1.0 to 2.0 sec (Default : 1.2 sec)	Cutoff Factor >> 1.0 to 2.0 sec (Default : 1.2 sec)	H_z Derivative Time >> <i>Derivative Time for Heat Pre-dominant zone</i> 0 to 600 sec (Default : 16 sec)
		Cycle Time >> 0.5 to 100.0 sec (Default : 10.0 sec)
		Overshoot Inhibit >> Enable Disable (Default : Disable)
		Cutoff Factor >> 1.0 to 2.0 sec (Default : 1.2 sec)

SUPERVISORY > PASSWORD

Parameters	Settings (Default Value)
Change Password >>	1000 to 1999 (Default : 123)

SUPERVISORY > EXI

Parameters	Settings (Default Value)
Exit Setup Mode >>	Yes No (Default : No)

FACTORY

FACTORY > CONTROL SENSOR INPUT

Parameters	Settings (Default Value)
Input Type >>	Refer Table 1 (Default : RTD Pt100)
Signal LO >>	
Signal HI >>	
Range LO >>	-199.9 to RANGE HI (Default : 0.0)
Range HI >>	RANGE LO to 999.9 (Default : 100.0)

FACTORY > MAPPING PARAMETERS

Parameters	Settings (Default Value)		
Mapping Points >>	0 to 4 (Default : 0)		
Input Type >>	Refer Table 1 (Default : RTD Pt100)		
Signal LO >>	Input Type	Settings	Default
	0 to 20mA	0.00 to Signal High	0.00
	4 to 20mA	4.00 to Signal High	4.00
	0 to 5V	0.000 to Signal High	0.000
	0 to 10V	0.00 to Signal High	0.00
	1 to 5V	1.000 to Signal High	1.000
Signal HI >>	Input Type	Settings	Default
	0 to 20mA	Signal Low to 20.00	20.00
	4 to 20mA	Signal Low to 20.00	20.00
	0 to 5V	Signal Low to 5.000	5.000
	0 to 10V	Signal Low to 10.00	10.00
	1 to 5V	Signal Low to 5.000	5.000
Range LO >>	-199.9 to RANGE HI (Default : 0.0)		
Range HI >>	RANGE LO to 999.9 (Default : 100.0)		

FACTORY > ALARM PARAMETERS

Parameters	Settings (Default Value)
Hysteresis >>	0.1 to 99.9 (Default : 0.2)
Inhibit >>	Yes No (Default : Yes)

FACTORY > HEAT COOL SELECT

Parameters	Settings (Default Value)
Control Strategy >>	Heat Only Cool Only Heat + Cool (Default : Heat + Cool)
Control Strategy : Cool Only	
Time Delay (sec) >>	0 to 1000 Sec (Default : 200 Sec)
Control Strategy : Heat + Cool	
Compressor Strategy >>	CONT. OFF CONT. ON SP BASED PV BASED (Default : CONT. ON)

CONT. ON	SP BASED	PV BASED
Time Delay (sec) >> 0 to 1000 Sec (Default : 200 Sec)	Boundary Set Value >> 0 to 100 (<i>Resolution 0.1°C for RTD / DC Linear</i> & 1°C for Thermocouple)	Time Delay (sec) >> 0 to 1000 Sec (Default : 200 Sec)
	Control Zones >> Single Dual (Default : Single)	
	Time Delay (sec) >> 0 to 1000 Sec (Default : 200 Sec)	

FACTORY > SOAK TIMER PARAMETERS

Parameters	Settings (Default Value)
Enable >>	No Yes (Default : No)
Start Band >>	0 to 999.9 (Default : 0.5)
Holdback Strategy >>	None Up Down Both (Default : None)
Hold Band >>	0.1 to 999.9 (Default : 0.5)
Heat Off >>	No Yes (Default : No)
Cool Off >>	No Yes (Default : No)
Power Recovery >>	Abort Restart Continuous (Default : Restart)

FACTORY > DOOR OPEN

Parameters	Settings (Default Value)
Enable >>	No Yes (Default : No)
Start Band >>	0 to 999.9 (Default : 0.5)
Holdback Strategy >>	None Up Down Both (Default : None)
Hold Band >>	0.1 to 999.9 (Default : 0.5)
Heat Off >>	No Yes (Default : No)
Cool Off >>	No Yes (Default : No)
Power Recovery >>	Abort Restart Continuous (Default : Restart)

FACTORY > MAINS FAILURE

Parameters	Settings (Default Value)
Enable >>	No Yes (Default : No)
Start Band >>	0 to 999.9 (Default : 0.5)
Holdback Strategy >>	None Up Down Both (Default : None)
Hold Band >>	0.1 to 999.9 (Default : 0.5)
Heat Off >>	No Yes (Default : No)
Cool Off >>	No Yes (Default : No)
Power Recovery >>	Abort Restart Continuous (Default : Restart)

FACTORY > SMS ALERT

Parameters	Settings (Default Value)
Enable >>	No Yes (Default : No)
Start Band >>	0 to 999.9 (Default : 0.5)
Holdback Strategy >>	None Up Down Both (Default : None)
Hold Band >>	0.1 to 999.9 (Default : 0.5)
Heat Off >>	No Yes (Default : No)
Cool Off >>	No Yes (Default : No)
Power Recovery >>	Abort Restart Continuous (Default : Restart)

FACTORY > FACTORY DEFAULT

Parameters	Settings (Default Value)
Set To Default >>	Yes No (Default : No)

FACTORY > PASSWORD

Parameters	Settings (Default Value)
Set To Default >>	Yes No (Default : No)

FACTORY > EXIT

Parameters	Settings (Default Value)
Exit Setup Mode >>	Yes No (Default : No)

TABLE 1

TABLE 1		
What it means	Range (Min. to Max.)	Resolution
Type J Thermocouple	0 to +960°C	Fixed 1°C
Type K Thermocouple	-200 to +1376°C	
Type T Thermocouple	-200 to +385°C	
Type R Thermocouple	0 to +1770°C	
Type S Thermocouple	0 to +1765°C	
Type B Thermocouple	0 to +1825°C	
Type N Thermocouple	0 to +1300°C	
Reserve	Reserved for customer specific Thermocouple type not listed above. The type shall be specified in accordance with the ordered (optional on request) Thermocouple type.	
3-wire, RTD Pt100	-199.9 to 600.0°C	Fixed 0.1°C
0 to 20mA DC current	-199.9 to 999.9 units	Fixed 0.1 unit
4 to 20mA DC current		
0 to 5.0V DC voltage		
0 to 10.0V DC voltage		
1 to 5.0V DC voltage		

FRONT PANEL KEYS


Symbol	Key	Function
	Scroll	Press to scroll through various Process Information Screens in Normal Operation Mode.
	Alarm Acknowledge	Press to acknowledge and mute (if active) alarm output.
	DOWN	Press to decrease the parameter value. Pressing once decreases the value by one count; keeping pressed speeds up the change.
	UP	Press to increase the parameter value. Pressing once increases the value by one count; keeping pressed speeds up the change.
	SET-UP	Press to enter or exit set-up mode.
	ENTER	Press to store the set parameter value and to scroll to the next parameter.

PV ERROR INDICATIONS

Message	Error Type	Cause
OPEN	Sensor Open	Sensor (RTD Pt100) Broken / Open
OVR	Over-range	Temperature above Max. Specified Range
UNR	Under-range	Temperature below Min. Specified Range

ELECTRICAL CONNECTIONS

4 Mapping Channels

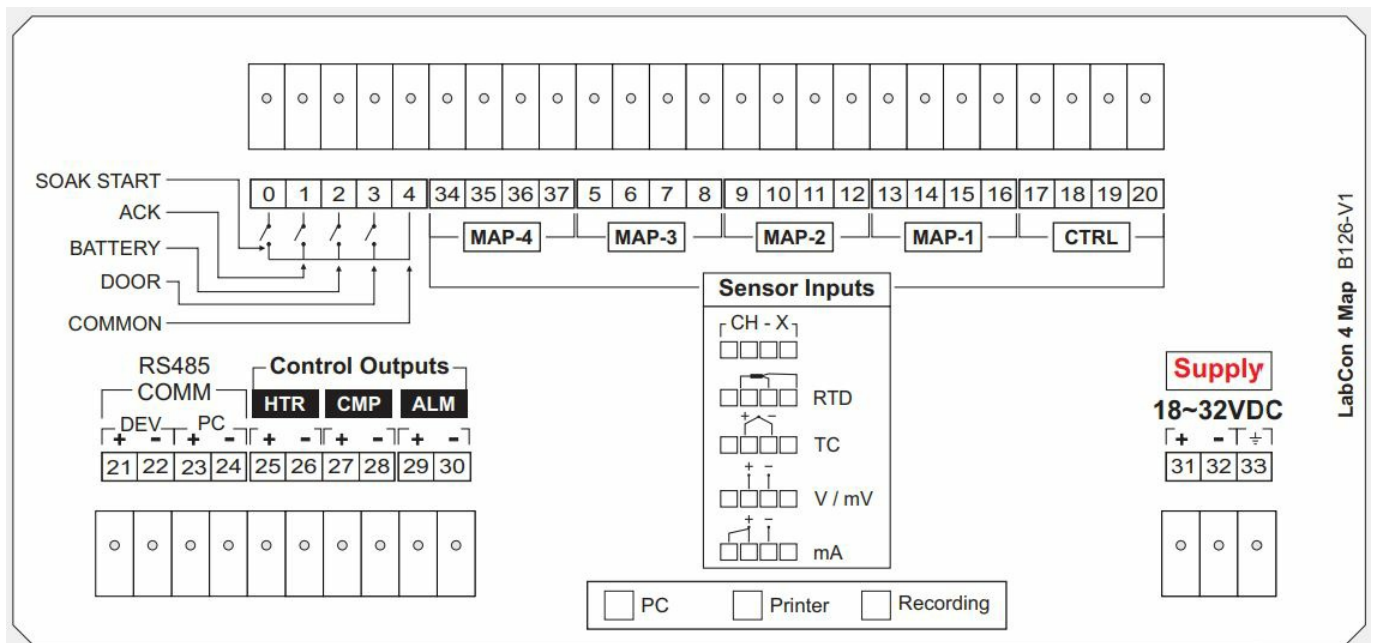

CONTACT

Address: 101, Diamond Industrial Estate, Navghar, Vasai Road (E), Dist. Palghar – 401 210.

Sales: 8208199048 / 8208141446

Support: 07498799226 / 08767395333

E: sales@ppiindia.net,

support@ppiindia.net

Documents / Resources

	PPI LabCon Map PC Recording 4 Channel Mapping PC Software [pdf] User Manual LabCon Map PC Recording 4 Channel Mapping PC Software, LabCon Map PC, Recording 4 C hannel Mapping PC Software, 4 Channel Mapping PC Software, Mapping PC Software, PC Sof tware
--	---

References

- [Factory Automation Products India | Automation Solutions in India](#)