



QAREQU C60 Microscope Camera User Manual

[Home](#) » [QAREQU](#) » QAREQU C60 Microscope Camera User Manual 

QAREQU C60 Microscope Camera



Contents

- [1 Connection](#)
- [2 Buttons](#)
- [3 Specifications](#)
- [4 Appendix: About Micro SD Card](#)
- [5 Appendix: Using On Computer](#)
- [6 Customer Support](#)
- [7 Documents / Resources](#)
 - [7.1 References](#)
- [8 Related Posts](#)

Connection






HDMI

Connect camera to monitor directly with HDMI cable, monitor must support 1080P resolution, otherwise you cannot see full size video.

USB

Connect camera to computer with USB cable, it will auto turn on and does not require DC 12 power. Refer to Appendix: Using on computer for more information

Buttons

-  **MENU** When power supply is connected it will turn on after 3-second. Indicator light turns to green. Press and hold the button for 3-second to power off.
- After power on, short-press to enter Menu, press again back to previous menu level.
-  Digital zoom in (max 7X) under camera mode.
- Move cursor up in Menu.
-  Digital zoom out under camera mode.
- Move cursor down in Menu.
-  Enter video record mode, press Camera icon to start/ stop record.
- Back to previous menu level in Menu.
-  Take photo, indicator light will turn to red.
- Select/ Confirm in Menu.

Specifications

Photo Pixel (max)	60MP (4:3) 40MP (16:9)	Lens Mount	C-mount
Video Resolution	2K 30FPS 1080P 60FPS	Micro SD Storage	Min 4GB Max 64GB
HDMI Output	1080P 60FPS	DC Power	12V 1A
CMOS Sensor	1/2.33 inch	Size	63 x 73 x 37mm (2.5 x 2.9 x 1.5")
CMOS Pixel	1.33 x 1.33um	Weight	150g (0.33 lb)

Appendix: About Micro SD Card

- Camera must under power off mode when insert/withdraw card, otherwise it will turn off and need to power on again.
- Card must be formatted inside camera before record or taking photo (Menu > Format > OK)

Appendix: Using On Computer

- It is driver- free on Windows/ Mac computer, not for mobile devices.
- Micro SD card cannot be inserted, other wise it will work as removable drive for files preview/ playback/ transfer.
- Download freeware like OBS or AMCap, or other compatible software for operation.
- Connect USB cable from camera to computer, it will auto turn on, DC power supply is not required.
- Device must be correctly selected from software menu if you have webcam or other camera connected.

6. All buttons and menu of camera will be disabled, operation is done on software only.

Customer Support

 facebook.com/qarequ

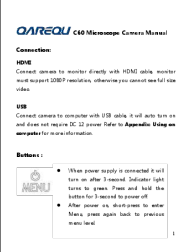
 qarequ@outlook.com

 www.qarequ.com

Copyright @ QAREQU All Rights Reserved

QAREQU

Documents / Resources

	QAREQU C60 Microscope Camera [pdf] User Manual C60 Microscope Camera, C60, Microscope Camera, Camera
---	---

References

- [User Manual](#)

[Manuals+](#), [Privacy Policy](#)

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.