

[Skip to content](#)

## Manuals+

User Manuals Simplified.



# TraceableLIVE Thermometer App Instruction Manual

[Home](#) » [TraceableLive](#) » TraceableLIVE Thermometer  
App Instruction Manual



## TraceableLIVE™ QUICK START GUIDE

Contents [hide](#)

[1 Thermometer App](#)

[2 Documents / Resources](#)

[2.1 References](#)

[3 Related Posts](#)

## Thermometer App

### 1. Create Account

Visit <https://www.traceablelive.com>

A screenshot of the TraceableLIVE web interface. It features a light blue background with a large blue 'T' logo at the top. Below the logo are two white input fields for 'Username' and 'Password'. A link for 'Forgot password?' is positioned below the password field. At the bottom, there are two buttons: 'New Account' and 'Login'.

Click on New Account.

### 2. Enter Account/User

Information

Fill in all required fields.

**Note:** Password must be a minimum of 8 characters including a number and a special character (!@#\$\$%^&\*).

Register a new account

\*required fields

<b>COMPANY*</b> Company	<b>ADMINISTRATOR FIRST NAME*</b> Administrator First Name
<b>ADDRESS*</b> Address Line 1: Address Line 2: City: State: ZIP Code: Country:	<b>ADMINISTRATOR LAST NAME*</b> Administrator Last Name  <b>ADMINISTRATOR EMAIL*</b> Administrator Email  <b>ADMINISTRATOR PHONE NUMBER*</b> Administrator Phone Number
<b>SUBSCRIPTION*</b> <input checked="" type="radio"/> Monthly <input type="radio"/> Annual	<b>PASSWORD*</b> Password
<b>ADD-ONS</b> <input type="radio"/> SMS	<b>PASSWORD CONFIRMATION*</b> Password Confirmation

Click Register

Register

Payment Information

Congrats! You're almost there - Your TraceableLive account has been created but we need you to fill out your payment preferences before you can add a new device. Your account will only be charged when you add a new device.

**PAYMENT INFORMATION\***

Card Number: \_\_\_\_\_

Expiration: - Select - / - Select -

CVV: \_\_\_\_\_

Cardholder: \_\_\_\_\_

Your account will only be charged after devices or services have been added.

**BILLING ADDRESS\***

Address Line 1: \_\_\_\_\_

Address Line 2: \_\_\_\_\_

City: \_\_\_\_\_

State: \_\_\_\_\_

ZIP Code: \_\_\_\_\_

Country: - Select One -

Save Payment Info

- Enter Payment Info./ Confirm Charge screen does not appear, go to log in screen, account settings and add payment information.

If this

**Note:** Payment must be set up prior to adding a device to your service.

**Click Save Payment Info**



4. Begin WiFi Setup

“AP” will appear on display.

On device Press/HOLD WiFi button for 3 secs.

Wait for a few seconds then press (no need to hold) WiFi button, “AP rEAdY” will appear on display.



800-517-8431

5 Commonwealth Ave Woburn, MA 01801 Phone 781-665-1400 Toll Free 1-




5. Connect Unit to WiFi network via Mobile device

device (smart phone, tablet, etc.), find CC6500-XXXX in available networks.

XXXX will be last 4 digits of unit's S/N# (on back of unit).

Using a WiFi enabled

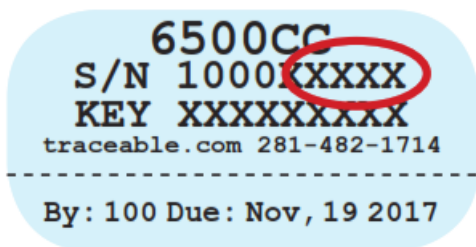


6. Open Web Browser on your mobile device  
192.168.1.1

Enter into web browser:

Select desired local network to which the device will be connected, and enter password for network. The most common network is WPA2.

7. Press Add/Confirm device restart

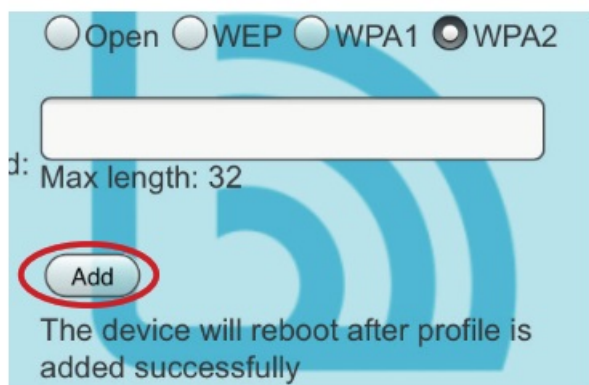

Click Add  
posting data.

Unit will reboot, connect to WiFi, and begin

**Note:** Upon successful connection, you will receive an error message in your browser (Example: Safari cannot open the page because the network connection was lost.), and device reboot may take up to 1 minute, mobile will exit WiFi.

**Note:** The WiFi symbol will flash on device display until the device posts data successfully.

**Note:** ALLOW UP TO 30 MINS FOR DEVICE DATA TO UPDATE AND START TRANSMITTING DATA.

If an ERR message occurs on device, press acknowledge button (✓) and repeat steps 4-7.

Log-in to TraceableLIVE account to manage device settings, alarms, users, and payment options.


8. **Enter Device Information (located on back of unit)**

Go to [www.traceablelive.com](http://www.traceablelive.com) Click on Add New Device

Complete all required

fields.

Enter S/N & Device Key traceable.com 281-482-1714 By: 100 Due: Nov, 19 2017 Device and account setup is

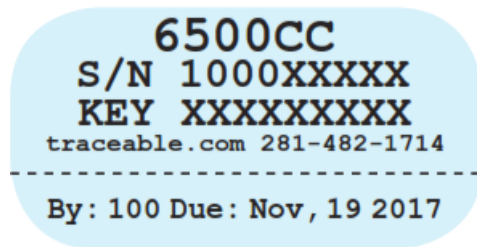

now complete.

## Documents / Resources



[TraceableLIVE Thermometer App](#) [pdf] Instruction Manual  
WD-99460-10, App

## References

- [TP Traceable Products](#)
- [traceablelive.com](#)
- [traceablelive.com](#)

## Manuals+,

- [home](#)
- [privacy](#)