



Link Report Creator Layout



TSI Link Report Creator Layout User Guide

[Home](#) » [tsi](#) » TSI Link Report Creator Layout User Guide 

Contents

- [1 TSI Link Report Creator Layout](#)
- [2 Product Information](#)
- [3 Product Usage Instructions](#)
- [4 Overview](#)
- [5 Prerequisites](#)
- [6 How to Use the Layout View](#)
- [7 Pictures in the Layout View](#)
- [8 Documents / Resources](#)
 - [8.1 References](#)
- [9 Related Posts](#)



TSI Link Report Creator Layout



Product Information

Specifications

- **Product Name:** TSI LinkTM Report Creator
- **Feature:** Layout View
- **Manufacturer:** TSI Incorporated
- **Website:** www.tsi.com

Product Usage Instructions

Opening the Layout View

To open the Layout View, select the Layout View option and click on the Add (+) button in the worksheet.

Selecting an Image

Update the Name of the Layout View and click on the IMPORT IMAGE button to select a file for visualization.

Placing Markers and Data on the Layout View

Choose the metric to display on the layout view from the drop-down menu. Place individual tests accordingly.

Inserting the Layout View into the Report Creator

After placing markers and data, insert the image with marker locations into the worksheet template. The image can be edited or deleted as needed.

Moving Images in the Layout View

To move the image within the worksheet, click on it, hold down the mouse button, and drag it to the desired location.

FAQ

- **Q: How can I edit marker locations on the Layout View?**

A: To edit marker locations, select the marker and drag it to the new desired location within the Layout View.

• Q: Can I add multiple images to the Layout View?

A: Yes, you can add multiple images to the Layout View by repeating the process of selecting an image and placing markers on each image.

Overview

The Layout View is a feature within many different TSI Link™ Report Creator Worksheets. It provides a spatial representation of your study, showing how your readings vary across the floor plan, around a specific work area, or along your fence line.

Prerequisites


Using the Layout View requires:

- TSI Link™ user account, i.e. a login and password
- TSI Link™ Report Creator subscription
- TSI Link™ Report Creator Microsoft® Excel® Add-in
- Data from a compatible TSI® device Loaded into a Report Creator Worksheet.

How to Use the Layout View


Opening the Layout View

By selecting the Layout View and using the Add (+) button a visualization of the data can be included in the worksheet.

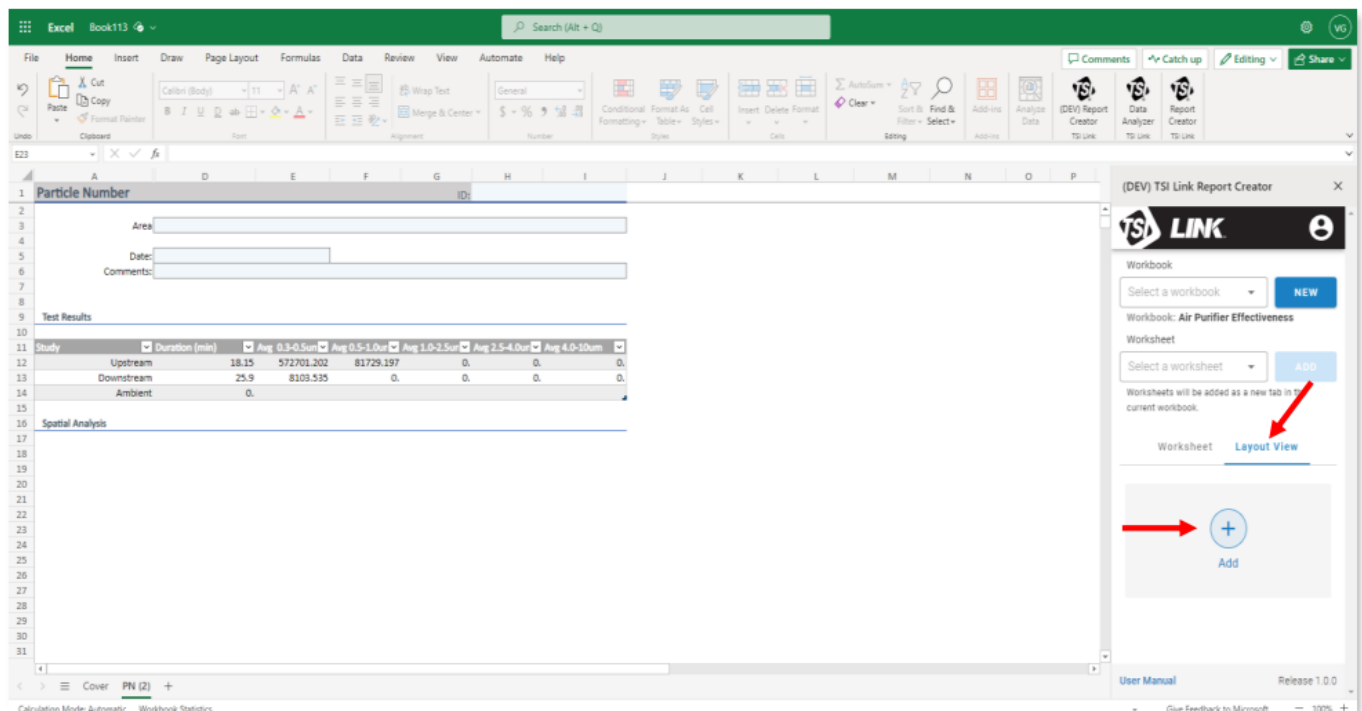


Selecting an Image

A new dialog is opened. Update the Name of the Layout View and use the IMPORT IMAGE button to select a file.

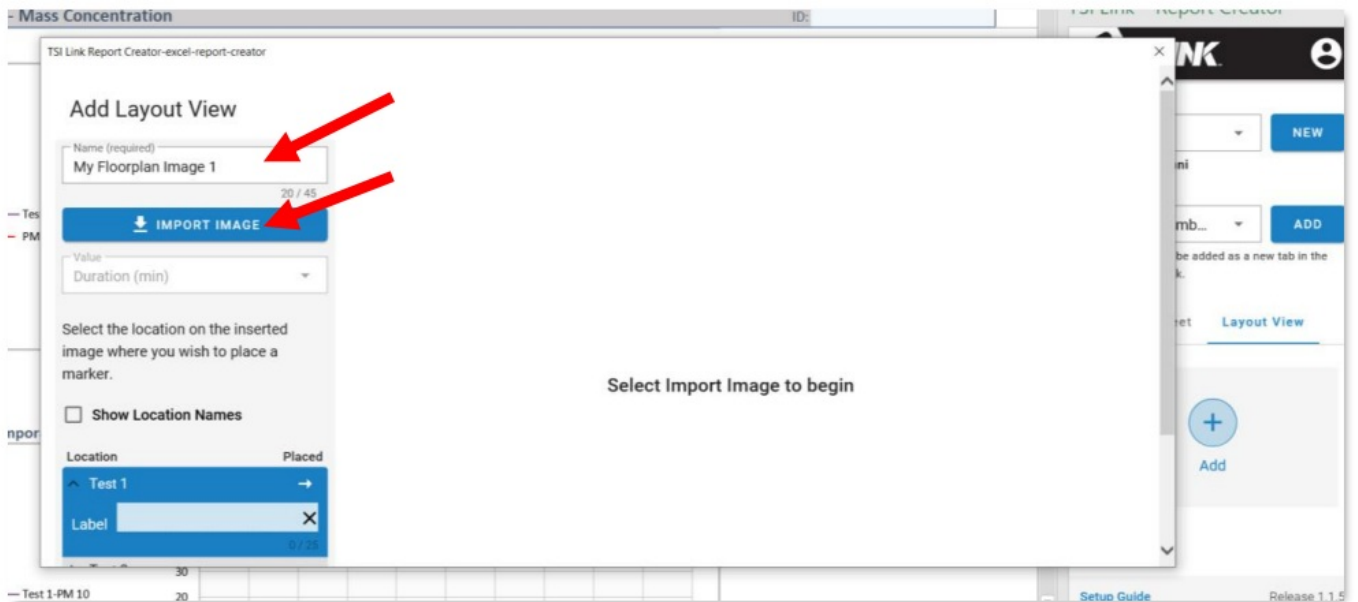

After using the IMPORT IMAGE button, a file picker will appear to help select an image file

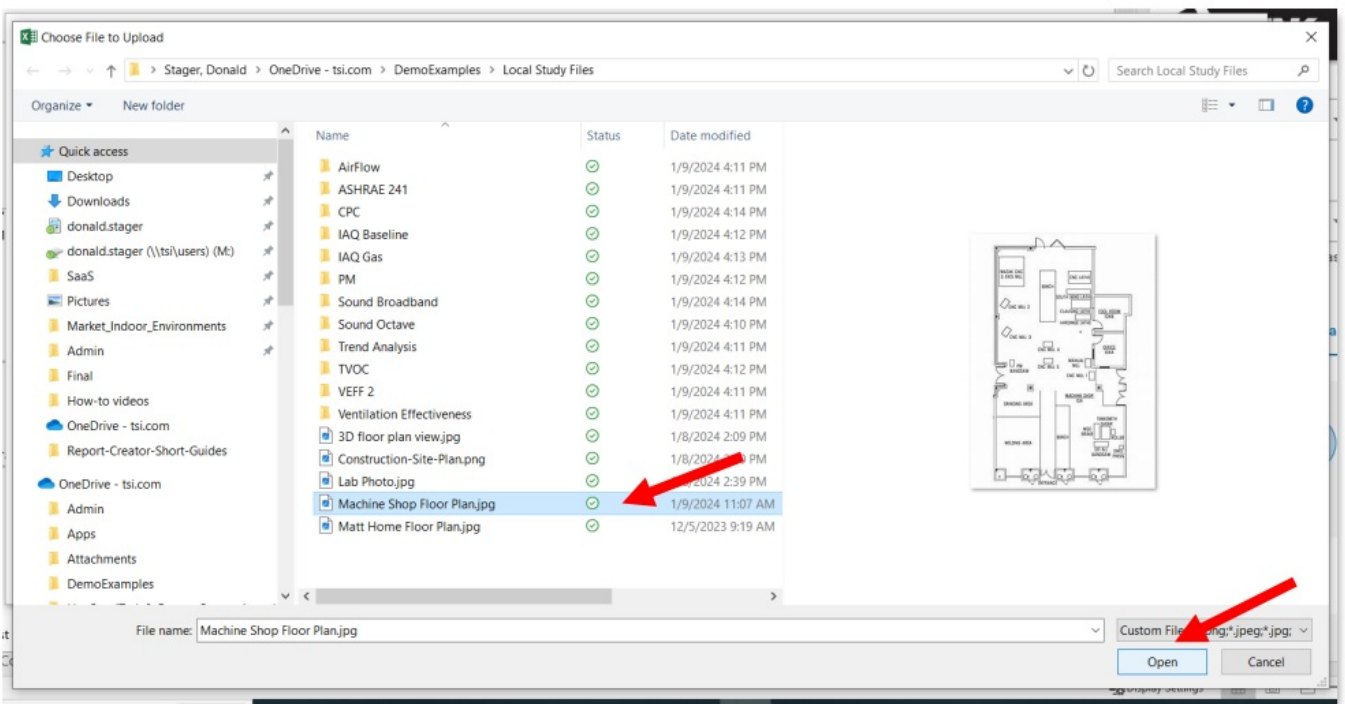

Placing the Markers and Data on the Layout View

Choose the metric you want to display on the layout view via:

- A – Drop down Menu

TSI Link Report Creator-excel-report-creator

Add Layout View

Name (required)
My Floorplan Image 1 20 / 45

IMPORT IMAGE

Value
Duration (min) **A**


Select the location on the inserted image where you wish to place a marker.

☐ Show Location Names

Location	Placed
Test 1	→
Label	✕
Test 2	

Expand All | Collapse All

CANCEL



- **B** – Scrolling to choose the desired metric

TSI Link Report Creator-excel-report-creator

Add Layout View

Name (required)
My Floorplan Image 1 20 / 45

IMPORT IMAGE

Value
Duration (min) **B**

Avg PM1.0

Avg PM2.5

Avg PM4.0

Avg PM10

Max PM1.0

Max PM2.5

Test 2

Expand All | Collapse All

CANCEL




Place the individual Tests:

- **C** – Choose Test
- **D** – Click on where image where you want the Marker

TSI Link Report Creator-excel-report-creator

Add Layout View

Name (required)
My Floorplan Image 1 20 / 45

 IMPORT IMAGE

Value
Avg PM2.5


Select the location on the inserted image where you wish to place a marker.

☒ Show Location Names

Location	Placed
> Test 1	✓
^ Test 2	→
Label	✗

Expand All | Collapse All

0 / 25




After the marker has been place it can be:

- **E** – Labeled
- **F** – Deleted, for example if the location is not quite right

TSI Link Report Creator-excel-report-creator

Add Layout View

Name (required)
My Floorplan Image 1 20 / 45

 IMPORT IMAGE

Value
Avg PM2.5

Select the location on the inserted image where you wish to place a marker.


☒ Show Location Names

Location	Placed
^ Test 1	✓
Label Bench	✗
^ Test 2	✓
Label	✗

Expand All | Collapse All


5 / 25

0 / 25

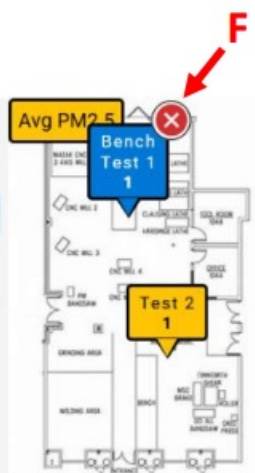


Inserting the Layout View into the Report Creator

The image with the marker locations can then be inserted into the worksheet template. It can also be edited or deleted.

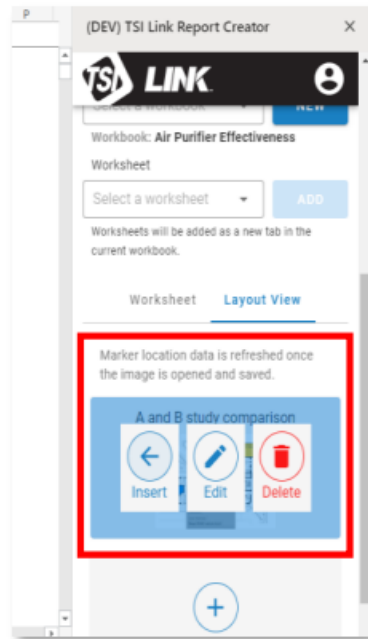

The image can be moved within the worksheet by clicking on it, holding down the mouse button and dragging it to the desired location.

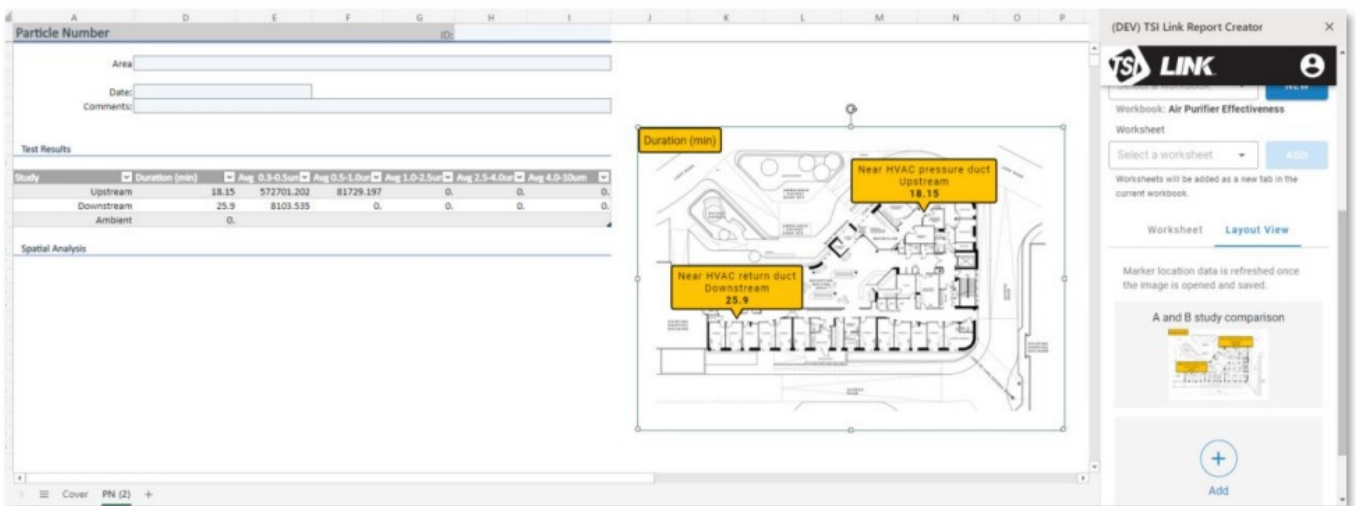

Pictures in the Layout View

Provide another strong representation.

TSI and the TSI logo are registered trademarks of TSI Incorporated in the United States and may be protected under other country's trademark registrations. Microsoft, Windows and Excel are registered trademarks of Microsoft Corporation

TSI Incorporated – Visit our website www.tsi.com for more information.

USA Tel: +1 800 680 1220

UK Tel: +44 149 4 459200

France Tel: +33 1 41 19 21 99

Germany Tel: +49 241 523030

India Tel: +91 80 67877200

China Tel: +86 10 8219 7688

Singapore Tel: +65 6595 6388

Documents / Resources

TSI Incorporated

TSI Link™ Report Creator

Layout View Guide



Contents

Introduction

Getting Started

Using the Report Creator

Customizing the Report

Printing the Report

Appendix

Overview

Getting Started

Using the Report Creator

Customizing the Report

Printing the Report

Appendix

Navigation

Getting Started

Using the Report Creator

Customizing the Report



Printing the Report

Appendix

[TSI Link Report Creator Layout](#) [pdf] User Guide

Standard Connected, Layout View Guide, Link Report Creator Layout, Link, Report Creator Lay out, Creator Layout, Layout

References

-  [Home | TSI](#)
-  [Home | TSI](#)
- [User Manual](#)

[Manuals+](#), [Privacy Policy](#)

This website is an independent publication and is neither affiliated with nor endorsed by any of the trademark owners. The "Bluetooth®" word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. The "Wi-Fi®" word mark and logos are registered trademarks owned by the Wi-Fi Alliance. Any use of these marks on this website does not imply any affiliation with or endorsement.