



Unity Lab Services Controlled Rate Freezer TSCM Battery Replacement Instructions

[Home](#) » [Unity Lab Services](#) » Unity Lab Services Controlled Rate Freezer TSCM Battery Replacement Instructions 

Contents

- [1 Unity Lab Services Controlled Rate Freezer TSCM Battery Replacement](#)
- [2 Installation Instructions](#)
- [3 Controlled Rate Freezer TSCM Battery Replacement](#)
- [4 Documents / Resources](#)
 - [4.1 References](#)
- [5 Related Posts](#)



Unity Lab Services Controlled Rate Freezer TSCM Battery Replacement



Installation Instructions

Revision Date: August 5, 2021


Controlled Rate Freezer TSCM Battery Replacement

1. Login to touchscreen. Turn unit off by pressing power button & message will pop up confirming shut down command. After screen processes power off command flip rocker switch on right side of unit and unplug from wall.
2. Remove screws on right side wall panel to gain access to internal deck
3. Remove side panel and battery is located on right side. Typically, the battery color is grey & blue or black.
4. Disconnect red and black wires from battery terminals
5. Remove battery bracket nut. Slide battery bracket towards the front of the unit and lift to remove. Remove old battery and replace with new one.
6. Complete Reverse order after new battery is installed.

Note: Ensure to connect wires correctly to battery terminals: Red (+) wire to red positive terminal & black (-) wire to negative black terminal

www.unitylabservices.com/contactus

Documents / Resources

	<p>Unity Lab Services Controlled Rate Freezer TSCM Battery Replacement [pdf] Instructions Controlled Rate Freezer TSCM Battery Replacement, Controlled Rate Freezer TSCM, TSCM B attery Replacement, Freezer TSCM</p>
---	---

References

- [U Contact Us](#)