



# Unity Lab Services Heraguard ECO Clean Bench User Guide

[Home](#) » [Unity Lab Services](#) » Unity Lab Services Heraguard ECO Clean Bench User Guide 



## Contents

- [1 Heraguard ECO Clean Bench](#)
- [2 Documents / Resources](#)
  - [2.1 References](#)
- [3 Related Posts](#)

## Heraguard ECO Clean Bench

### Troubleshooting guide

How do I activate the UV light on my Heraguard ECO?

#### Heraguard ECO

Revision Date: November 10, 2014


#### How do I activate the UV light on my Heraguard ECO?

The Heraguard Eco must be turned off for the UV light to function.


1. Turn the cabinet blowers off ( hold the “ON” key for 10 seconds)
2. Install the UV light resistant cover.



3. Press and hold the UV light button for 10 seconds

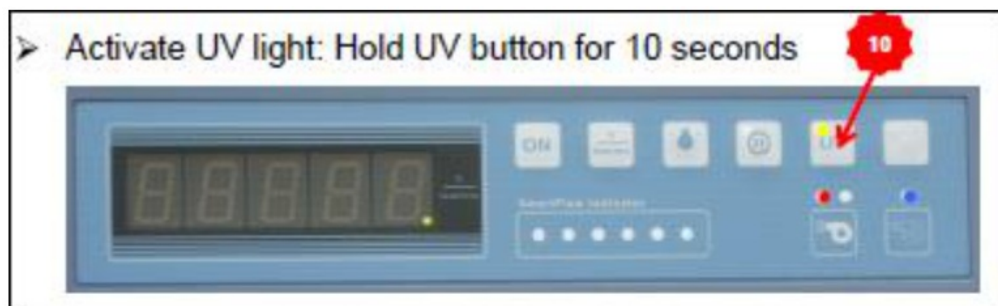

4. If the UV light still isn't functioning verify the UV bulb cover has been removed.

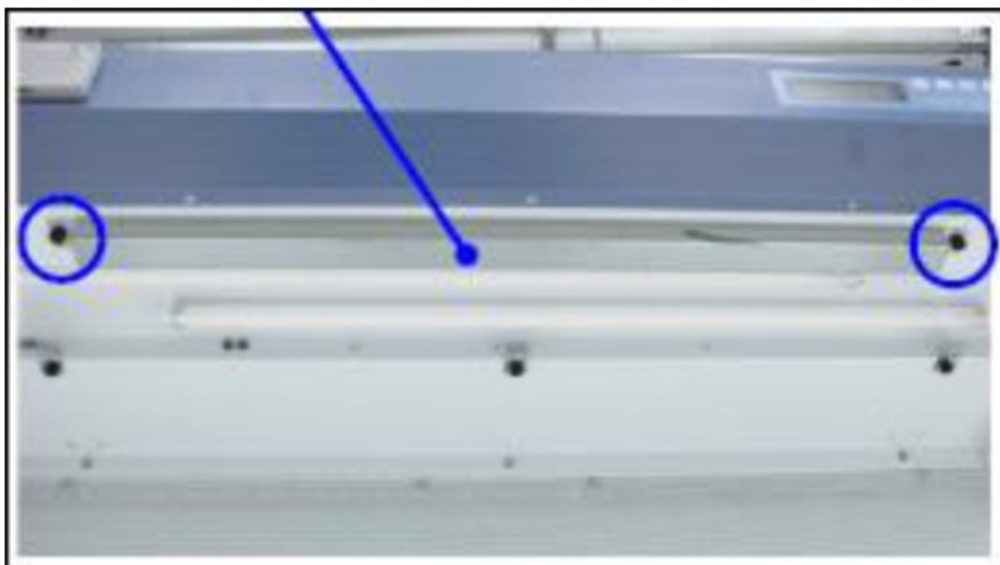

5. Please note the UV light resistant covers along with the UV lamp are recommended to be replaced every 1500 hours. See page 21, 37 in the operator's manual.

#### **UV bulb replacement**

- ▶ Align the pins of bulb with the sockets
- ▶ Insert the bulb into sockets
- ▶ Twist' turn until the bulb locks

- ▶ Test Install light cover



#### UV light resistant cover

- ▶ No interlock as with BSC
- ▶ Held in place with elastic cords



**Unity** Lab Services

A Thermo Fisher Scientific Brand

[www.unitylabservices.com/contactus](http://www.unitylabservices.com/contactus)

#### Documents / Resources

	<p><a href="#">Unity Lab Services Heraguard ECO Clean Bench</a> [pdf] User Guide Heraguard ECO Clean Bench, Heraguard ECO, Clean Bench, Bench</p>
--	---

#### References

- [U Contact Us](#)

